

# Oxis Research<sup>®</sup>

...ENHANCING OXIDATIVE STRESS TECHNOLOGY



Catalog and Technical Resource

# OxisResearch®

*A DIVISION OF OXIS INTERNATIONAL, INC.*

## Catalog and Technical Resource

Copyright © 2003 All rights reserved. OXIS International, Inc.

6040 North Cutter Circle, Suite 317  
Portland, OR 97217-3935 • USA

800.547.3686 • +1.503.283.3911  
fax: +1.503.283.4058

[www.oxisresearch.com](http://www.oxisresearch.com)  
[information@oxisresearch.com](mailto:information@oxisresearch.com)



*"Dedicated to helping researchers understand the relevance of Oxidative and Nitrosative Stress on biological systems"*

### *A Letter To Our Patrons...*

This catalog introduces fourteen new assay kits to compliment our already broad range of oxidative stress related testing products. In the past two years we have identified new biomarkers of relevance to oxidative and nitrosative stress while creating simpler, more robust tests for existing biomarkers where possible.

With the knowledge that commercially available tests related to oxidative protein modifications were lacking in the marketplace we tasked our researchers with identifying biomarkers that may be useful in this area. The result of this program was the development of two new assays for detection of enzyme activities subject to oxidative modification (i.e. **Aconitase-340** for mitochondrial and tissue related studies and  **$\alpha_1$ -Antiproteinase-410** for plasma relevant studies). The activity of both of these enzymes has been shown to be mediated by oxidative stressors and in the case of the  $\alpha_1$ -antiproteinase, the deactivation is thought to be reversible via the methionine sulfoxide/methionine sulfoxide reductase pathway. We also brought to market a new **Nitrotyrosine ELISA** kit based on its relevance to peroxynitrite mediated damage to proteins.

In the area of lipid peroxidation we sought to improve upon existing methods such as our popular **LPO-586** assay. With the release of our **MDA-586** and **HAE-586** assays we made it possible for the first time to measure highly reactive hydroxyalkenal by-products without the contribution of the more readily detectable malondialdehyde (MDA). At the same time we improved upon the MDA side of the assay by providing validated methods for detection of "free" and "total" MDA in biological samples. We also provide third derivative spectroscopy procedures for those interested in a more definitive test for plasma MDA using our **MDA-586** product. We recommend our **MDA-586** for plasma and tissue samples and our **HAE-586** for detection of HAE formation in low protein samples such as oils and other lipid extracts.

We now offer improved methods for determining 8-epi-prostaglandin  $F_{2\alpha}$  levels in various biological samples. Two new assays we introduced, one for determination of **urinary 8-isoprostane** that does not require solid phase extraction and an entirely new assay for the **metabolite of 8-isoprostane** in urine. We believe the metabolite assay may offer significant advantages over traditional measurements of the parent compound.

Other additions to our assay portfolio include a **glutathione-S-transferase** assay, **glucose-6-phosphate dehydrogenase** assay, and a **nitric oxide** assay.

We expect that the next two years will yield more and better technologies for the study of oxidative and nitrosative stress. We are committed to quickly identifying and facilitating their commercialization in order that we might remain your preferred provider in this area.

Thank you for your patronage, we look forward to being of continued service to you and hope we can be an integral part of your successful investigation.

Our very best regards,

The OXIS Team

*On the cover:*

Cytosolic, suborganelle, systemic and interstitial processes illustrating the potential biological roles of free radical reactions.

*On the back cover:*

Cytosolic, suborganelle, systemic and interstitial processes including references to applicable OxisResearch® products.

## OXIS International Inc.

OXIS International (OTCBB: OXIS) (Nouveau Marché: OXIS), headquartered in Portland, Oregon, focuses on developing technologies and products for research and commercialization related to oxidative stress. OXIS has been awarded more than 45 patents in the United States and internationally, sells 31 oxidative stress assays of which 19 were internally developed, and offers an extensive catalog of products focused on the study of oxidative stress. OXIS also holds the rights to three therapeutic classes of antioxidants for the treatment of diseases of oxidative stress.



### What is Oxidative Stress?

Oxidative stress is caused by the unchecked activities of free radicals and reactive oxygen species (ROS). These are highly reactive, unstable molecules that, in healthy cells, are created and neutralized as part of normal cell life. However, if the production of free radicals and ROS exceeds the body's natural defense mechanisms, oxidative damage to vital molecules in the body can occur, leading to cell death.

### OXIS is an expert in oxidative stress

OXIS develops, manufactures, and markets products and technologies related to oxidative stress. OXIS currently offers more than 80 research products for sale. At the heart of the *OxisResearch*<sup>®</sup> product line are research kits that simplify the testing of Oxidative, Antioxidant, Nitrosative and Inflammatory Biomarkers thought to play a central role in many human degenerative diseases including cancer, diabetes, atherosclerosis, stroke, Alzheimer disease, Parkinson disease and amyotrophic lateral sclerosis (ALS). Researchers at major universities throughout the world regularly use OXIS products for their oxidative-stress related testing needs. OXIS continues to develop new oxidative-stress products for the research market, releasing seven new assays in 2002. References to OXIS' research products can be found at the OXIS Citation Library ([www.oxisresearch.com](http://www.oxisresearch.com)), which also provides a link to Medline, where researchers can order full text citations and explore additional literature on related subjects.

OXIS is moving its research products into commercial applications. In 2002, OXIS announced plans to launch its Animal Health Profiling Program, which was created to bring to market screening tests to measure oxidative stress in cattle blood. A recent university study showed a correlation between an OXIS biomarker and the incidence of bovine respiratory disease. It is estimated that approximately 65% of the 38 million calves born in the U.S. will be affected with this disease, costing cattle producers \$1 billion annually. Early identification of susceptible cattle is essential to decrease these losses. OXIS believes that it will be the first company to offer a biomarker that provides a valuable analytical tool for managing the health of cattle.

OXIS is preparing to bring a unique compound called Ergothioneine into the highly lucrative nutraceuticals market. Ergothioneine is one of the more powerful antioxidants available. With its proprietary method for commercially producing Ergothioneine, OXIS plans to become the primary worldwide source for this molecule. OXIS is testing and plans to commercialize Ergothioneine as a nutraceuticals dietary supplement. Additionally, several companies, including two major international cosmetics firms, currently use Ergothioneine as an anti-aging agent in skin-care products.

The OXIS leading therapeutic drug candidate, BXT-51072, completed a Phase IIA clinical trial in inflammatory bowel disease (IBD) in 1999 and further clinical studies continue to be reviewed based on the Company's ability to fund these activities. Two other therapeutic programs are in the pre-clinical stage of development.

A Letter To Our Patrons ..... 2

About OXIS International, Inc. .... 3

Ordering Instructions ..... 6

International Distributors ..... 8

Oxisresearch.com Online ..... 10

Selecting OxisResearch® Products ..... 12

OxisResearch® Products ..... 16

Bioxytech® Test Kits

*Antioxidant Biomarkers*

21010 SOD-525 Assay ..... 16

21011 GSH-400 Assay ..... 18

21014 Plasma GPx-EIA Assay ..... 20

21017 cGPx-340 Assay ..... 22

21018 GR-340 Assay ..... 24

21023 GSH-420 Assay ..... 26

21040 GSH/GSSG-412 Assay ..... 28

21042 Catalase-520 Assay ..... 30

21045 G6PD ; 6PGD-340 Assay ..... 32

21046 GST-340 Assay ..... 34

21052 AOP-490 Assay ..... 36

*Oxidative Biomarkers*

21012 LPO-586 Assay ..... 38

21019 8-Isoprostane Assay ..... 40

21024 H<sub>2</sub>O<sub>2</sub>-560 Assay ..... 42

21025 LPO-560 Assay ..... 44

21026 8-OHdG Assay ..... 46

21041 Aconitase-340 Assay ..... 48

21043 HAE-586 Assay ..... 50

21044 MDA-586 Assay ..... 52

21047 α<sub>1</sub>-AP-410 Assay ..... 54

21048 Urinary 8-EPI-PGF<sub>2α</sub> Assay ..... 56

21049 Urinary F<sub>2</sub>-Isoprostane Metabolite Assay ..... 58

*Inflammatory Biomarkers*

21013 MPO-EIA Assay ..... 60

21015 Lactof-EIA Assay ..... 62

*Nitrosative Biomarkers*

21055 Nitrotyrosine-EIA Assay ..... 64

22110 Nitric Oxide Assay (enzymatic) ..... 66

22111 Nitric Oxide Assay (non-enzymatic) ..... 68

22112 Nitric Oxide Synthase Assay (radioactive arginine) ..... 70

22113 Nitric Oxide Synthase Assay (colorimetric) ..... 72

Antioxidants

26511 L-Ergothioneine (100 mg) ..... 74

26524 L-Ergothioneine (250 mg) ..... 74

## Controls

27617	Cellular Glutathione Peroxidase Control .....	76
27618	Glutathione Reductase Control .....	76
27619	Cu/Zn SOD Control .....	76
27622	Catalase Control ( <i>Aspergillus Niger</i> ) .....	77

## Antibodies

24310	Anti-Myeloperoxidase, (Monoclonal) .....	78
24311	Anti-Lactoferrin, (Monoclonal) .....	78
24312	Anti-Nitrotyrosine .....	78
24316	Anti-Catalase .....	78
24317	Anti-Cytochrome P450 2E1 .....	79
24318	Anti-Cytochrome P450 3A4 .....	79
24319	Anti-Glutathione S-Transferase A1-1 .....	80
24320	Anti-Glutathione S-Transferase M1-1 .....	80
24321	Anti-Glutathione S-Transferase P1-1 .....	80
24322	Anti-Cu/Zn - Superoxide Dismutase .....	81
24323	Anti-Mn Superoxide Dismutase .....	81
24325	Anti-4-Hydroxy-2-nonenal (100mg) .....	82
24326	Anti-8-Hydroxy-2'-deoxyguanosine (100mg) .....	82
24327	Anti-4-Hydroxy-2-nonenal (20mg) .....	82
24328	Anti-8-Hydroxy-2'-deoxyguanosine (20mg) .....	82
25310	Anti-Endothelial Nitric Oxide Synthase .....	82
25311	Anti-Inducible Nitric Oxide Synthase .....	83
25312	Anti-Brain Nitric Oxide Synthase .....	83

## Enzymes

25447	Bovine Cu/Zn Superoxide Dismutase (bSOD) (25mg) .....	84
25414	Bovine Cu/Zn Superoxide Dismutase (bSOD) (250mg) .....	84
25415	Bovine Cu/Zn Superoxide Dismutase (bSOD) (500mg) .....	84
25416	Bovine Cu/Zn Superoxide Dismutase (bSOD) (1g) .....	84
25430	Catalase, <i>Aspergillus niger</i> .....	84
25431	Catalase, Bovine Liver .....	85
25432	Catalase, Human Erythrocyte .....	85
25433	Glutathione Reductase, Yeast .....	86
25434	Glutathione-S-Transferase A1-1 Recombinant Human .....	86
25435	Glutathione-S-Transferase M1-1 Recombinant Human .....	87
25436	Glutathione-S-Transferase P1-1 Recombinant Human .....	87
25437	Myeloperoxidase, Human Polymorphonuclear Leukocytes .....	88
27510	Brain Nitric Oxide Synthase, Recombinant Rat .....	88
27511	Constitutive Nitric Oxide Synthase Recombinant Rat Neuronal .....	89
27512	Constitutive Nitric Oxide Synthase Recombinant Human, High Purity .....	89
27513	Inducible Nitric Oxide Synthase, Mouse Macrophage .....	90

## Markers

26595	4-Hydroxy-non-enal-diethylacetal (50mg) .....	92
26515	4-Hydroxy-non-enal-diethylacetal (100mg) .....	92
26516	4-Hydroxy-non-enal-diethylacetal (1g) .....	92
26594	4-Hydroxy-hex-2-enal-diethylacetal (50mg) .....	92
26518	4-Hydroxy-hex-2-enal-diethylacetal (100mg) .....	92
26519	4-Hydroxy-hex-2-enal-diethylacetal (1g) .....	92
26525	8-Hydroxyguanosine .....	93
26526	4-Hydroxy-2-nonenal .....	93
26527	3-Nitro-L-Tyrosine .....	94
26571	8-Hydroxy-2'-deoxyguanosine .....	94

## Specialty Chemicals

26510	2,3-dimethoxy-1-naphthoquinone .....	95
26513	1-Methyl-4-vinyl-pyridium Triflouromethane Sulfonate .....	95

Product Index ..... 96

Product Numeric Index ..... 98

BIOXYTECH® is a registered trademark of OXIS International, Inc.

## Ordering Instructions

---

### Pricing and Terms of Sale

U.S. payment terms are net 30 days with approved credit. Bank and wire fees are not to be deducted from the invoice amount. Outside the U.S., payment terms may vary by country and distributor. We reserve the right to change prices without prior notification.

### Ordering Information

We accept orders via mail, phone, fax, e-mail and online. Our business hours are Monday thru Friday, 8:00 am to 5:00 pm PST. No orders are shipped over the weekend due to the temperature requirements for most of our products. Although written confirmation of an order is not required, please mark it "confirming order, do not duplicate" if one is sent.

When ordering, please provide the following information to ensure prompt and accurate service:

- Name of Institution
- Complete Billing and Shipping Address
- Purchase Order Number
- Catalog Number and Quantity Desired
- Contact Information

Orders must be backed by a valid Purchase Order or a Visa/Mastercard number. (Please have the account number and expiration date available)

Phone: 800.547.3686 USA  
503.283.3911  
Fax: 503.283.4058  
Internet: [www.oxisresearch.com](http://www.oxisresearch.com)  
Email: [information@oxisresearch.com](mailto:information@oxisresearch.com)

### International Orders

For all of our customers outside the U.S., please contact the OXIS distributor nearest you. (International Distributor list is on page 8) If your country isn't listed, feel free to contact us directly.

### Shipping

All shipments are FOB Portland, Oregon, USA. Unless otherwise specified, OXIS will select the freight carrier to be used. Shipping and handling charges are prepaid and added to the invoice utilizing a standard two-day delivery service. Overnight service is available upon request and is used when transporting frozen items.

The shipment of hazardous items is governed by the Department of Transportation (DOT) and the International Air Transport Association (IATA). These products require special packaging and will incur additional shipping charges as defined by freight carriers. Items shipped under refrigeration or frozen will incur additional charges as well.

### Material Safety and Data Sheets (MSDS)

Material Safety and Data Sheets will be provided as required by OSHA's Hazard Communication Law for products that are hazardous. Many OXIS products are new and experimental and have not been tested for toxicity. THE ABSENCE OF A WARNING STATEMENT DOES NOT IMPLY THAT THE PRODUCT IS NOT HAZARDOUS. Research products should only be utilized by qualified investigators or technically trained personnel working under the direct supervision of qualified investigators. It is the investigator's responsibility to ensure the safe handling of all products.

**Damaged or Lost Goods**

If a damaged product is found, please follow these steps:

1. Immediately report the damage or loss to the carrier.
2. Notify your distributor or OXIS customer service to process a replacement order.

**Return Policy**

A return materials authorization number must be obtained from an OXIS Customer Service Representative prior to returning any item. Return authorization requests must be made within ten working days of order receipt. Because of the perishable nature of all OXIS products, returns of incorrectly ordered items will not be accepted. Claims regarding suspected deficient product may be made at any time to OXIS International, Inc. Technical Services Department.

Outside the U.S., return policies may vary by country. Contact your distributor for assistance.

**Definitions and Abbreviations**

Room temperature:	20° C
Refrigeration:	2°– 8° C
Frozen:	–20° C
Deep Freeze:	–70° C

**Warranty**

All products offered by OxisResearch®, Inc. are guaranteed to meet label claims during the period covered by the expiration date on the product, provided storage and use instructions are followed. Items that do not require an expiration date are guaranteed for the stability period as stated on the package insert provided the storage and use instructions are followed. Products shipped in accordance with purchase orders placed from this catalog conform to the composition and purity as described on product labelling at the time of shipment. OXIS reserves the right to modify composition and purity as listed in this catalog due to changes in market conditions, raw materials or other factors. ALL OTHER GUARANTEES, WARRANTIES, CONDITIONS AND REPRESENTATIONS, EITHER EXPRESSED OR IMPLIED, WHEREBY ARISING UNDER ANY STATUTE, LAW, COMMERCIAL USAGE OR OTHERWISE, INCLUDING IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE, ARE HEREBY EXCLUDED.

# OxisResearch®

Division of OXIS International, Inc.  
6040 North Cutter Circle, Suite 317  
Portland, OR 97217-3935 • USA

800.547.3686 • +1.503.283.3911  
fax: +1.503.283.4058

[www.oxisresearch.com](http://www.oxisresearch.com)  
[information@oxisresearch.com](mailto:information@oxisresearch.com)



## OxisResearch® International Distributors

**Australia**

Sapphire Bioscience Pty. Ltd.  
 phone: +61 2 9966 9662  
 fax: +61 2 9966 9661  
 e-mail: info@sapphirebio.com.au  
 website: www.sapphirebio.com.au

**Denmark**

Electra-Box Diagnostica ApS  
 phone: +45 4453 6211  
 fax: +45 4453 6212  
 e-mail: henning.andersen@electrabox.com  
 website: www.electrabox.se

**Austria**

Sanova Pharma GesmbH Medical Systems  
 phone: +43 1 80104 256  
 fax: +43 1 80104 204  
 e-mail: sanova.diagnostik@sanova.at  
 website: www.sanova.at

**Eastern Europe**

Intimex (Holdings) Ltd.  
 phone: +44 1256 766073  
 fax: +44 1256 766701  
 e-mail: sales@intimex.org

**Benelux**

ImmunoSource BV  
 phone: +32 3385 36 85  
 fax: +32 3384 38 18  
 e-mail: info@immunosource.com  
 website: www.immunosource.com

**Finland**

Electra-Box Diagnostica AB  
 phone: +46 8 712 3000  
 fax: +46 8 712 6509  
 e-mail: peter.blom@electrabox.com  
 website: www.electrabox.se

**Canada**

Medicorp, Inc.  
 phone: +1 514.733.1900  
 fax: +1 514.733.1212  
 e-mail: mktg@medicorp.com  
 website: www.medicorp.com

**France**

Tebu-Bio  
 phone: +33 0173 3042 1880  
 fax: +33 0173 3042 1882  
 e-mail: uk@tebu-bio.com  
 website: www.tebu-bio.com/contact/france.htm

**China**

Dakewe Biotech Co.  
 phone: +86 755 264 12280  
 fax: +86 755 264 12562  
 e-mail: info@dakewe.com  
 website: www.dakewe.com

**Germany**

NatuTec GmbH  
 phone: +49 069 67 73 78 0  
 fax: +40 069 67 73 78 79  
 e-mail: info@natutec.com  
 website: www.natutec.com

**Germany** (continued)

## Tebu-Bio

phone: +69 801013 0  
 fax: +69 801013 20  
 e-mail: germany@tebu-bio.com  
 website: www.tebu-bio.com/contact/  
 germany.htm

**Iceland**

## Electra-Box Diagnostica ApS

phone: +45 4453 6211  
 fax: +45 4453 6212  
 e-mail: peter.blom@electrabox.com  
 website: www.electrabox.se

**Italy**

## Prodotti Gianni SpA

phone: +39 2 5097 232  
 fax: +39 2 5097 276  
 e-mail: rravizza@prodottigianni.com  
 website: www.prodottigianni.com

**Japan**

## Funakoshi Co., Ltd.

phone: +81 3 5684 1620  
 fax: +81 3 5684 1775  
 e-mail: reagent@funakoshi.co.jp  
 website: www.funakoshi.co.jp

## WAKO Pure Chemical Industries, Ltd.

phone: +81 6 203 2371  
 fax: +81 6 222 1203  
 website: www.wako-chem.co.jp

**Korea**

## Woongbee MediTech Inc.

phone: +82 2 569 0781  
 fax: +82 2 569 0784  
 e-mail: woongbee@woongbee.com  
 website: www.woongbee.com

**New Zealand**

## Sapphire Bioscience Pty. Ltd.

phone: +61 2 9966 9662  
 fax: +61 2 9966 9661  
 e-mail: info@sapphirebio.com.au  
 website: www.sapphirebio.com.au

**Norway**

## Electra Box AS

phone: +47 32 23 7950  
 fax: +47 32 23 7949  
 e-mail: bente.skar@electrabox.com  
 website: www.electrabox.se

**Poland**

## Biokom

phone: +48 (10xx) 22 720 0145  
 fax: +48 (10xx) 22 720 5443  
 e-mail: Guest@biokom.com.pl  
 website: www.biokom.com.pl

**Spain**

## Deltaclon

phone: +34 91 764 53 64  
 fax: +34 91 381 42 75  
 e-mail: info@deltaclon.com  
 website: www.deltaclon.com

**Sweden**

## Electra-Box Diagnostica AB

phone: +46 8 712 3000  
 fax: +46 8 712 6509  
 e-mail: peter.blom@electrabox.com  
 website: www.electrabox.com

**Taiwan**

## Level Biotechnology

phone: +886 2269 59935  
 fax: +886 2269 50403  
 e-mail: info@level.com.tw  
 website: www.level.com.tw

**United Kingdom**

## Bio Stat Limited

phone: +44 161 483 5884  
 fax: +44 161 483 5778  
 e-mail: sales@bio-stat.com  
 website: www.bio-stat.com

## Tebu-Bio

phone: +44 01733 421880  
 fax: +44 01733 421882  
 e-mail: uk@tebu-bio.com  
 website: www.tebu-bio.com/contact/  
 great\_britain.htm

## Online with OxisResearch.com

*Your complete online source for oxidative and nitrosative stress research products, services, applications and assistance...*

### OxisResearch® Products

Offering full text descriptions, specifications, application information, insert bulletins, references and citations.

Oxisresearch.com is the leading online source for oxidative and nitrosative stress research products including assays, antioxidants, antibodies, enzymes, controls, and specialty chemicals.

### Application Resources

Review a variety of helpful OXIS citations, informative articles and product references on oxidative and nitrosative stress research.

### Product Support

Access frequently asked product questions, visit our literature library for ordering and downloading, and find the contact information for our expert application staff to help you solve your research problems.

### News Releases

Learn about our latest oxidative and nitrosative stress developments and products.

### Company Resources

Access our calendar of events, our corporate news releases and much more...



## OxisResearch® Products Online

*Your convenient online source for oxidative and nitrosative stress research products.*

### Sales Assistance

Click on "Order" on any page in our web site to use our OxisResearch® Products Online area to check pricing and place orders.

If you have any questions about ordering OxisResearch® products, we offer quick links to find OXIS distributors worldwide, or to access contact information for our OXIS sales team always ready to help you in your selection process.



**Visit [www.oxisresearch.com](http://www.oxisresearch.com)**

*"Dedicated to helping researchers understand the relevance of Oxidative and Nitrosative Stress in biological systems"*

## Interactive Product Citations Library

Quickly find informative summaries of oxidative and nitrosative stress related research citations using OxisResearch® test kits in a variety of applications...

You can easily sort our citation database online by choosing criteria to match your model system from simple pull-down menus. You can reference the biomarker category, animal species, sample type, and/or the disease condition parameters of interest to your research model.

**OxisResearch**  
A Division of Oxis International, Inc.

CONTACT US | ORDER | HOME

CORPORATE | OXISRESEARCH | PRODUCTS | WHAT'S NEW | APPLICATION

**Citation Library** - Select your desired application criteria and click on the "Search" button:

Biomarker Category: Any Biomarker  
Species: Any Species  
Sample Type: Any Sample  
Condition: Any Condition  
OXIS Product (Analyte): Any Product

**SEARCH**

*Easily access specific product citations using simple pull-down menus*

**Citations:**  
You found 140 citations that meet the criteria: Biomarker = "Any Biomarker", Species = "Any Species", Sample = "Any Sample", Product = "Any Product", and Condition = "Any Condition"...

Number: 3  
PMUI Link: [11127413](#)  
Product Info: [LPO-586 \(Malondialdehyde and/or 4-Hydroxyalkenals\)](#)  
Link:

Biomarker: Oxidative-Lipids  
Species: Mouse  
Samples: Kidney, Liver, Lung, Spleen, Thymus  
Condition: Toxicology  
Citation:

*Quick links are provided for easy access to Medline Citations*

Dermal and systemic toxicity after application of semisynthetic metal-working fluids (MWF) in B6C3F1.  
Al-Humad et al., 2000. J. Toxicicol. Env. Health 61: 579-89

Summary:  
Mice were fed with/without 5% ethanol in water and Metal Working Fluids (MFWs) were dispensed on unshaved back skin from cervical to sacral region.  
Lipid Peroxidation was assayed in liver, kidney, lung, spleen and thymus using the OXIS LPO-586 test for malondialdehyde (MDA).

Click on the search button and our interactive OxisResearch® Product Citations Library quickly matches your chosen project criteria against all of the summaries in our collection.

The resulting list of citations displays:

- OxisResearch® products used in each application with a link to their complete specifications,
- All biomarker categories, animal species, sample types and disease conditions described in the research project,
- A citation summary of the researcher's particular methods and results using OXIS products,
- Background information where helpful,
- Citation PMUI number linking you directly to the referenced full text Medline® citation.

**Visit [www.oxisresearch.com](http://www.oxisresearch.com)**

Valuable resources for your  
oxidative and nitrosative stress research...

## Exploring Oxidative And Nitrosative Stress With OxisResearch®

- I. Oxidative and Nitrosative Stress Overview
- II. Quantifying Oxidative and Nitrosative Stress
- III. Choosing OxisResearch® Products

I. Oxidative and Nitrosative Stress Overview

Every cell has chemical reactions involving the oxidation and reduction of molecules. These reactions or redox pathways can lead to the production of free radicals.

**Free Radicals**

A free radical is any chemical species capable of independent existence possessing one or more unpaired electrons. Biological free radicals are thus highly unstable molecules that have electrons available to react with various organic substrates.

Many free radicals are the result of naturally occurring processes such as oxygen metabolism and inflammatory processes. For example, when cells use oxygen to generate energy, free radicals are created as a consequence of ATP production by the mitochondria. Exercise can increase the levels of free radicals as can environmental stimuli such as ionizing radiation (from industry, sun exposure, cosmic rays, and medical X-rays), environmental toxins, altered atmospheric conditions (e.g. hypoxia and hyperoxia), ozone and nitrous oxide (primarily from automobile exhaust). Lifestyle stressors such as cigarette smoking and excessive alcohol consumption are also known to affect levels of free radicals. Radical species may combine to form other more damaging or toxic species such as peroxynitrite ( $O=NOO^-$ ), a product of superoxide and nitric oxide radical reaction.

Free radicals react with key organic substrates such as lipids, proteins, and DNA. Oxidation of these biomolecules can damage them, disturbing normal functions and may contribute to a variety of disease states.

It has been noted that certain organ systems are predisposed to greater levels of oxidative or nitrosative stress. Those organ systems most susceptible to damage are the pulmonary system (exposed to high levels of oxygen), the brain (exhibits intense metabolic activity yet has lower levels of endogenous antioxidants), the eye (constantly exposed to damaging UV light), circulatory system (victim to fluctuating oxygen and nitric oxide levels) and reproductive systems (at risk from the intense metabolic activity of sperm cells). Nearly every organ system can be found to have an Oxidative or Nitrosative "Achilles heel". With the current understanding that free radicals can act as cell signaling or "messenger" agents it is likely that they also play a role in normal cellular function as well as various disease etiologies.

Reactive Oxygen Species (ROS) is a term collectively describing radicals and other non-radical reactive oxygen derivatives. These intermediates may participate in reactions giving rise to free radicals or that are damaging to organic substrates. ROS in living organisms include the following:

<u>Radicals</u>		<u>Non-Radicals</u>	
Hydroxyl	$OH^\bullet$	Peroxynitrite	$ONOO^-$
Superoxide	$O_2^{\bullet-}$	Hypochloric acid	$HOCl$
Nitric oxide	$NO^\bullet$	Hydrogen Peroxide	$H_2O_2$
Thyl	$RS^\bullet$	Singlet Oxygen	$^1\Delta_g (^1O_2)$
Peroxyl	$RO_2^\bullet$	Ozone	$O_3$
Lipid peroxy	$LOO^\bullet$	Lipid peroxide	$LOOH$

Reactive Nitrogen Species (RNS) are radical nitrogen-based molecules that can act to facilitate nitrosylation reactions. Reactive Nitrogen Species (RNS) include:

Nitrous oxide	$NO^\bullet$	Nitrosyl cation	$NO^+$
Peroxynitrite	$OONO^-$	Nitrogen dioxide	$NO_2^\bullet$
Peroxynitrous acid	$ONOOH$	Dinitrogen trioxide	$N_2O_3$
Nitroxyl anion	$NO^-$	Nitrous acid	$HNO_2$
Nitryl chloride	$NO_2Cl$		

Many other radical species can be formed by biological reactions, for example: phenolic and other aromatic species are often formed during xenobiotic metabolism as part of natural detoxification mechanisms.

**Oxidative Stress**

Oxidative stress occurs when the generation of ROS in a system exceeds the system's ability to neutralize and eliminate them. The imbalance can result from a lack of antioxidant capacity caused by disturbance in production, distribution, or by an over-abundance of ROS from an environmental or behavioral stressor. If not regulated properly, the excess ROS can damage a cell's lipids, protein or DNA, inhibiting normal function. Because of this, oxidative stress has been implicated in a growing list of human diseases as well as in the aging process.

**Nitrosative Stress**

Nitrosative stress occurs when the generation of RNS in a system exceeds the system's ability to neutralize and eliminate them. Nitrosative stress may lead to nitrosylation reactions that can alter protein structure thus inhibiting normal function.

**Antioxidants**

An extensive, highly effective group of protective agents and defense mechanisms referred to collectively as the Antioxidant Defense System (ADS), acts to regulate oxidative reactions.

The ADS includes enzymes and antioxidants to prevent the start of oxidative damage and/or control its spread. There are also enzymes to repair oxidative damage, and mechanisms to target damaged molecules for destruction and replacement. Essential antioxidants are either endogenous (internally synthesized) or exogenous (consumed). They are typically categorized as scavenger antioxidants and prevention antioxidants.

Scavenger antioxidants remove ROS and include the following categories:

- Small molecule antioxidants include both water-soluble compounds such as Vitamin C or glutathione and lipid soluble compounds such as Vitamin E, carotenes, lipoic acid, and Coenzyme Q<sub>10</sub>.
- Large molecule "enzyme" antioxidants include superoxide dismutase (SOD) that detoxifies the superoxide ion, catalase, which deals with hydrogen peroxide (H<sub>2</sub>O<sub>2</sub>), and glutathione peroxidase (GPx) whose function is to detoxify cellular peroxides. These enzymes must be synthesized by cells and are subject to genetic and/or macromolecular regulatory mechanisms.

Preventative antioxidants hinder the formation of new ROS. These antioxidants are proteins that bind ROS to protect essential proteins. The group includes albumin, metallothionein, transferrin, ceruloplasmin, myoglobin, and ferritin.

**ROS and RNS today:**

Researchers are now making rapid progress in understanding the role of oxidative and nitrosative stress in cardiovascular diseases such as atherosclerosis, ischemia/reperfusion injury, restenosis and hypertension; cancer; inflammatory diseases such as acute respiratory distress syndrome (ARDS), asthma, inflammatory bowel disease (IBD), dermal and ocular inflammation and arthritis; metabolic disease such as diabetes; and diseases of the central nervous system (CNS) such as amyotrophic lateral sclerosis (ALS), Alzheimer's, Parkinson's, and stroke. The increased awareness of oxidative stress related to disease and the need to measure the delicate balance that exists between free radicals and the systems in place to regulate them has given rise to a demand for new research tools.

**II. Quantifying Oxidative and Nitrosative Stress**

Researchers wishing to assess the level and/or effect of oxidative or nitrosative stress on a given system typically have two basic methods at their disposal:

1. Measure levels of specific ROS and RNS present in the system
2. Measure the effect (including damage and ADS remodeling) of ROS and RNS on a given system.

Since ROS and RNS are generally highly reactive, exacting Spin Trap methodologies are typically required to adequately detect and quantify their existence in biological samples. For this reason the second method is the most often employed strategy in the study of oxidative and nitrosative stress. To accomplish this, researchers can look at modifications to a system's regulatory mechanisms (e.g. endogenously expressed antioxidants), or look at by-products of the applied stressor (e.g. damage/modification of lipids, proteins or DNA). Researchers may also wish to look at causative factors such as the events facilitating an inflammatory cascade such as neutrophil activation and subsequent release of myeloperoxidase and/or lactoferrin.

The OxisResearch® product line provides a “toolbox” for researchers to accomplish the basic goals as described earlier. Generally speaking, we offer researchers four basic types of tools...

- A. Assays
- B. Antibodies
- C. Enzymes
- D. Controls

...directed against four major areas of biomarkers:

- 1) Oxidative
- 2) Nitrosative
- 3) Antioxidant (regulatory)
- 4) Inflammatory (pro-oxidant)

**A. Assays**

Tools for assaying the presence and quantity of:


- 1. Oxidative Biomarkers:
  - a) Lipids (MDA, HAE, LOOH, 8-Isoprostane, 8-iso-metabolite)
  - b) Proteins (Aconitase,  $\alpha_1$ -Antiproteinase, Nitrotyrosine)
  - c) DNA (8-OHdG)
- 2. Nitrosative Biomarkers:
  - a) Nitric Oxide (Enzymatic and non-enzymatic Griess methods)
  - b) Nitric Oxide Synthase (Colorimetric and Radioactive methods)
  - c) Peroxynitrite (Nitrotyrosine)
- 3. Antioxidant Biomarkers:
  - a) Glutathione (GSH/GSSG, cGPx, pGPx, GR)
  - b) Superoxide Dismutase (SOD-525)
  - c) Catalase (Catalase-520)
- 4. Inflammatory Biomarkers:
  - a) Neutrophils (Myeloperoxidase, Lactoferrin)



**B. Antibodies**


Tools for semi-quantitative (e.g. blotting and/or IHC) detection of:

- 1. Oxidative Biomarkers:
  - a) Lipids (Anti-4-Hydroxynonenal)
  - b) Proteins (Anti-Nitrotyrosine)
  - c) DNA (Anti-8-Hydroxydeoxyguanosine)
- 2. Nitrosative Biomarkers:
  - a) Nitric Oxide Synthase (Anti-iNOS, Anti-cNOS, Anti-bNOS)
  - b) Peroxynitrite (Anti-Nitrotyrosine)
- 3. Antioxidant Biomarkers:
  - a) Glutathione (Anti-GST M1-1, Anti GST A1-1, Anti GST P1-1)
  - b) Superoxide Dismutase (Anti-Cu/ZnSOD, Anti-MnSOD)
  - c) Catalase (Anti-Catalase)
  - d) Cytochromes (Anti P450-2E1, Anti-P450 3A4)
- 4. Inflammatory Biomarkers:
  - a) Neutrophils (Anti-Myeloperoxidase, Anti-Lactoferrin)

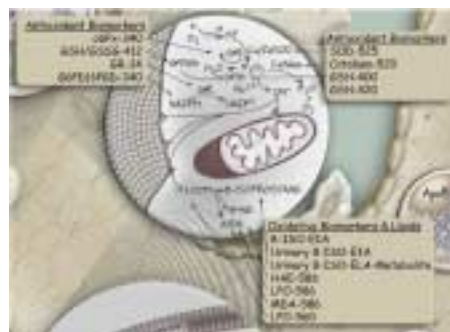


**C. Enzymes**

1. Oxidative Biomarkers:
  - a) Proteins (Anti-Nitrotyrosine)
  - b) DNA (Anti-8-hydroxydeoxyguanosine)
2. Nitrosative Biomarkers: (NO associated enzymes)
  - a) Nitric Oxide Synthase (iNOS, cNOS, bNOS)
3. Antioxidant Biomarkers:
  - a) Glutathione (GST M1-1, GST A1-1, GST P1-1)
  - b) Superoxide Dismutase (bovine Cu/ZnSOD, MnSOD)
  - c) Catalase (Catalase)
4. Inflammatory Biomarkers:
  - a) Neutrophils (Myeloperoxidase)

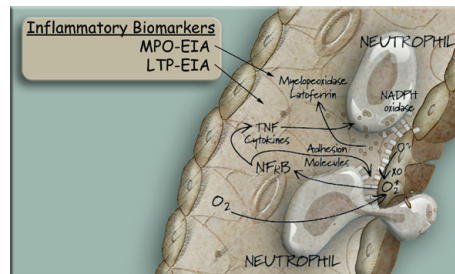

**D. Controls/Markers**

1. Oxidative Biomarkers:
  - a) Lipids (4-HNE, 4-HNE-diethylacetal, 4-HHE-diethylacetal)
  - b) Proteins (3-Nitro-L-Tyrosine)
  - c) DNA (8-OHG 8-OHdG)
2. Nitrosative Biomarkers:
  - a) Peroxynitrite (3-nitro-L-Tyrosine)
3. Antioxidant Biomarkers:
  - a) Glutathione (GR-control, cGPx-control)
  - b) Superoxide Dismutase (Cu/ZnSOD-control)
  - c) Catalase (Catalase-control)

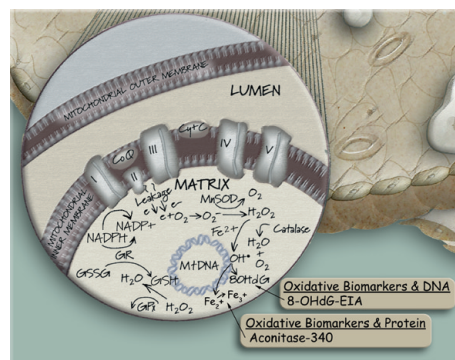

III. Selecting OxisResearch® Products

**Catalog Product Information**

OXIS International, Inc. through its OxisResearch® line of products has been addressing the need for measuring the balance that exists between free radicals, and the systems in place to regulate them. OXIS offers an array of innovative products for testing of Antioxidant, Oxidative, Nitrosative and Inflammatory Biomarkers. To help you more effectively evaluate our products for your particular system, this catalog thoroughly describes each OxisResearch® kit including useful applications and reference materials for all. Each assay listing shows the following information in detail:

- Product Description
- Method of Use
- Product References
- Summary of Specifications
- Selected Research Applications

If unable to locate information specific to your model or sample type, the OXIS application specialists are available for consultation and can recommend a product or technique best suited for your particular application.

OXIS products are playing a key role in what is now known about the role of reactive oxygen (ROS) and nitrogen (RNS) species in physiology and disease. Our products have been cited in over 150 peer reviewed international scientific journals. We hope the selected citations listed in this catalog will convey the multiplicity of current uses for our products and perhaps inspire you in finding new ones.

For more OxisResearch® product information and assistance, visit our web site at [www.oxisresearch.com](http://www.oxisresearch.com), email us at [information@oxis.com](mailto:information@oxis.com), or call OXIS International at 800.547.3686 (USA) to speak with one of our expert application specialists.



## Bioxytech® SOD-525 Assay

21010

*Quantitative Measurement of Superoxide Dismutase Enzyme Activity***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21010


The Superoxide dismutases (SOD) are a family of metalloproteins. These enzymes catalyze the decomposition of the superoxide radical into hydrogen peroxide and oxygen.



SOD constitutes the primary defense against the toxic effect of superoxide in aerobic organisms. Currently, four isoforms of SOD have been identified; three (Cu/Zn-SOD, Mn-SOD and EC-SOD) in eukaryotes and one (Fe-SOD) in Prokaryotes. In mammals, two of three SOD isoforms have Cu and Zn in their catalytic center and are localized intracellularly (Cu/Zn-SOD or SOD-1) or extracellularly (EC-SOD or SOD 3). SOD-1 (M.W. 32 kDa) has been found in the cytoplasm, nuclear compartments and lysosomes in mammals. SOD 3 (heterodimer, M.W. 135 kDa with high affinity to heparin) is the most recently discovered and least characterized member of the SOD family. It was first detected in plasma, lymph, ascites and cerebrospinal fluids. The expression of SOD-3 is highly restricted to specific cell types and tissues where its activity can exceed that of SOD-1. Mn-SOD or SOD-2, the third form of mammalian SOD, has manganese as a cofactor and is localized to the mitochondria of aerobic cells.

**SOD-525 Method**

The Bioxytech SOD-525 Assay takes advantage of a proprietary reagent whose alkaline autoxidation is accelerated by the presence of SOD. The autoxidation product absorbs light at 525 nm. Rates of auto-oxidation are determined spectrophotometrically in the presence and absence of an SOD sample.

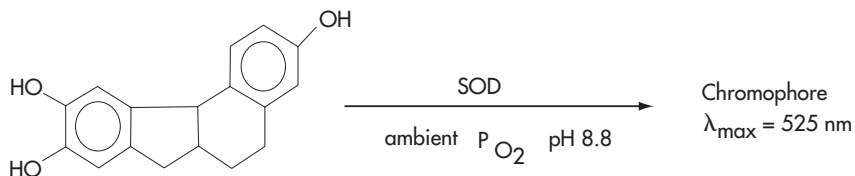

The ratio of the two rates measured can be converted directly to SOD concentration in the sample with the aid of a table (or equation) provided. No calibration is required however an SOD control (27617) is available for use with this kit.

U.S. Patent No. 5,543,298

**References**

1. Nebot C., Moutet M., Huet P., Xu J.Z., Yadan J. C., Chaudière J. Spectrophotometric assay of superoxide dismutase activity based on the activated autoxidation of a tetracyclic catechol. *Analytical Biochemistry* 1993; 214:442-451.
2. Zelko, I. Et al. Superoxide Dismutase Multigene Family: A Comparison of CuZn-SOD (SOD1), Mn-SOD (SOD2), And EC-SOD (SOD-3) Gene Structures, Evolution, and Expression. *2002 FRBM* 33:337-349

**SOD-525 Product Summary**

Intended Use:	Quantitative measurement of superoxide dismutase activity. For Research Use Only. Not for Use in Diagnostic Procedures.		
Format:	100 Test Colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• Chromogenic Reagent (R1)</li> <li>• Mercaptan Scavenger (R2)</li> <li>• Buffer</li> </ul>		
Storage and Stability:	18 months from date of manufacture when stored as specified.		
Specimen Requirements:	Tissue and/or cell homogenates. (e.g. cell culture, WBC, RBC, etc.)		
Assay Precision:			
Replicates = 30	Level 1	Level 2	Level 3
	<u>Day1/Day2</u>	<u>Day1/Day2</u>	<u>Day1/Day2</u>
Mean (U-525/mL)	0.82/0.81	2.85/2.86	13.60/13.22
Standard Dev	0.06/0.10	0.25/0.23	1.29/1.42
Total % CV	7.32/12.35	8.77/8.04	9.48/10.74
Assay Range:	Assay has best precision at Vs/Vc ratios less than 8.		
Sensitivity:	LLD (SOD-525 U/mL) = 0.1 in the assay		

**SOD-525 Selected Research Applications**

The Bioxytech SOD-525 assay has been used in research for measuring both SOD-1 and SOD-2 in a variety of models including melatonin pharmacology, cobalt and copper toxicity, preeclampsia related fetal hypoxia, diabetes, neurological diseases (amyotrophic lateral sclerosis disease or ALS), aging, gastrointestinal disorders, cystic fibrosis, pancreatic-fibrosis, asthma, organ transplantation and blood transfusion.

Perinatal encephalopathy induced by hypoxia or ischemia is a major cause of childhood neurological disability. Using OXIS' SOD-525 assay in a rat model, the Okatani group showed that the melatonin application to the mother increased antioxidant enzyme activities, including SOD, in the fetal brain and may thereby provide indirect protection against free radical injury (1). Ishi et al. showed that SOD activity is increased in the renal cortex in a streptozotocin induced diabetic rat model (2). In asthma patients, reduction of SOD activity has been associated with the severity of the disease (3). Improper antioxidant supplementation was shown to cause depletion of endogenous antioxidants (SOD, GSH and cGPx) and increased lipid peroxidation in cystic fibrosis patient blood (4). The Matsumura group studied the pancreatic injury in rats showing that SOD activity was significantly reduced in rats with pancreatic-fibrosis (5). In aging human populations, decreased SOD activity accompanied by increased lipid peroxidation has been reported (6). The Abe group showed that both "neo-red cells" (RBC substitutes) and RBC maintained the same SOD activity and other physical properties such as oxygen consumption and biocompatibility in an organ transplantation model. Thus, they suggested that "neo-red cells", which contain SOD, may be useful as an artificial oxygen carrier in blood transfusions and may reduce oxidative stress in organ transplantations (7).

1. Melatonin increases activities of glutathione peroxidase and superoxide dismutase in fetal rat brain Okatani, Y. et al. (2000) *J. Pineal Res.* 28: 89-96.
2. Nitric Oxide synthesis and oxidative stress in the renal cortex of rats with diabetes mellitus. Ishi, N et al. *J. Am. Soc. Nephrol.* 12:1630-39.
3. Lipid peroxidation as determined by plasma isoprostanes is related to disease severity in mild asthma. Wood LG et al. (2000). *Lipids* 35(9): 967-74.
4. Increased plasma fatty acid concentration after respiratory exacerbations are associated with elevated oxidative stress in cystic fibrosis. Wood LG et al. (2002). *Am. J. Clin. Nutr.* 75: 668-75.
5. Study of free radical and pancreatic fibrosis-pancreatic fibrosis induced by repeated injections superoxide dismutase inhibitor. Matsumura N et al. (2001). *Pancreas* 22 (1) 53-61.
6. Aging is associated with increased lipid peroxidation in human hearts, but not with mitochondrial respiratory chain enzyme defects. Mior, O. et.al (2000) *Cardiovasc Res.* 47: 624-31.
7. Superoxide generation from human polymorphonuclear leukocytes by liposome-encapsulated hemoglobin. Abe, H et al. (2002). *Artificial Cells, Blood Substitute and Immob. Biotechnology* 29 (4): 275-283

## Quantitative Measurement of Glutathione


For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21011


Glutathione ( $\gamma$ -glutamylcysteinylglycine or GSH) is a naturally occurring tripeptide whose nucleophilic and reducing properties play a central role in metabolic pathways, as well as in the antioxidant system of most aerobic cells. GSH plays a critical role as a coenzyme with a variety of enzymes including glutathione peroxidase, glutathione S-transferase, and thiol transferase. GSH also plays major roles in drug metabolism, calcium metabolism,  $\gamma$ -glutamyl cycle, blood platelet and membrane functions. In addition, GSH is crucial to a variety of life processes, including the detoxification of xenobiotics, maintenance of the -SH level of proteins, thiol-disulfide exchange, removal of hydroperoxides and free radicals as well as amino acid transport across membranes. Physiological values of intracellular GSH generally range from 1 to 10 mM. Although several methods have been described for the assay of GSH, reliable ones are often labor intensive and not easy to use. The GSH-400 from OxisResearch® provides a simple, specific test for GSH that does not require an enzyme as a reagent. Because of its simplicity, the GSH-400 Assay method is especially well adapted to the assay of glutathione in large series of biological samples.

## GSH-400 Method

The GSH-400 method is based on a chemical reaction that proceeds in two steps. The first step leads to the formation of substitution products (thioethers) between a patented reagent, R1 (4-chloro-1-methyl-7-trifluoromethyl-quinolinium methylsulfate), and all mercaptans (RSH) present in the sample:

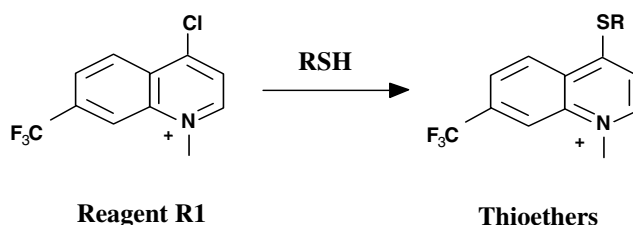

The second step is a  $\beta$ -elimination reaction which takes place under alkaline conditions. The reaction is mediated by an R2 reagent (30% NaOH) that transforms the specific substitution product (thioether) obtained with GSH into a chromophoric thione having a maximal absorbance wavelength at 400 nm:

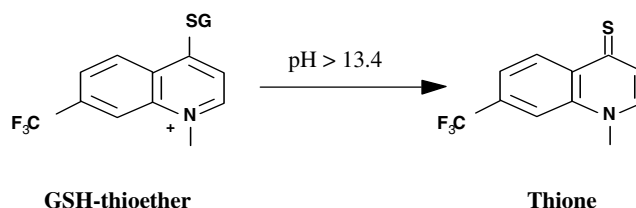

The GSH-400 method makes it possible to specifically assay glutathione with only one sampling and one colorimetric measurement. A modification of this method can be used to assay other mercaptans. This is based on the measurement of thioethers, which absorb light at 356 nm in the absence of reagent R2.

Researchers are cautioned that although only "reduced" GSH reacts with R1 to form specific GSH-thioethers, some GSSG will become reduced *in vitro* by normal means causing some detection of GSSG in the assay. For exacting measurements of "total" GSH, OXIS recommends the GSH-420 (21023) kit or for "ratio" measurement, the GSH/GSSG-412 (21040) kit.

U.S. Patent Number 5,817,520.

## References

1. Blume K. G., Paniker N. V., Beutler E. (1975) Glutathione (Flohé L., Benöhr H. Ch., Sies H, Waller H. D., Wendel A. eds.), pp. 157-165, Academic Press, New York.

**GSH-400 Product Summary**

Intended Use:	Quantitative measurement of glutathione in biological samples. For Research Use Only. Not For Use in Diagnostic Procedures.
Format:	100 Test Colorimetric
Kit Contents:	<ul style="list-style-type: none"> <li>• R1 Chromogenic Reagent</li> <li>• R2 (30% NaOH)</li> <li>• Buffer</li> </ul>
Storage and Stability:	18 months from date of manufacture when stored as specified.
Specimen Requirements:	Tissue homogenates and cell lysates
Accuracy	With a series of 30 measurements performed on the same day and under the same experimental conditions, using GSH (20-200 µM), the standard error of the mean value (SEM) was less than 2%.
Reproducibility	With the same experiments performed twice at three-day intervals, using the same samples, the new SEM calculated on the two measurement series was lower than 2%.
Sensitivity:	5 µmol/L in the final reaction mixture

**GSH-400 Selected Research Applications**

The OXIS GSH-400 has been on the market for approximately 10 years and the citations supporting its use are numerous. Here we present a few examples of the versatility of this assay as regards its use in various sample types and conditions.

Baumer et al. evaluated effect of ROS on various parameters of male fertility. They studied the influence of ROS generated by a xanthine (X)-xanthine oxidase system on equine sperm motility, viability, acrosomal integrity, membrane potential and membrane lipid peroxidation (1).

Elsewhere, Demols et al. studied cerulein-induced acute pancreatitis (AP) and the effect on GSH by NAC in prophylactic (NAC given before the induction) and therapeutic groups (NAC given after the induction). The authors showed that NAC administration abolished the decrease in pancreatic and hepatic tissue GSH that normally occurs during AP (2).

Lonigro, R et al. used L-buthionine-(S,R)-sulfoimine (BSO) to irreversibly inhibit  $\gamma$ -glutamylcysteine synthetase and tested FRTL-5 cells for the effects of diminished GSH. The authors showed a direct correlation between intracellular GSH concentration and the expression of thyroid specific genes: thyroglobulin (TO) thyroid peroxidases (TPO) which are mediated by the combination of TTF-1 (thyroid specific transcription factor), TTF-2 and Pax-8 (3).

Delcourt et al. evaluated GSH levels in erythrocytes of human subjects with cataracts showing that lower GSH levels correlate to increased incidence of cataracts (4).

Relating GSH content to immune function, Victor et al. performed an interesting timecourse study of GSH response to endotoxin mediated oxidative stress. They measured GSH levels in murine leukocytes isolated from peritoneum, thymus, spleen and axillary nodes showing that the endotoxin LPS causes decreased GSH in spleen and particularly axillary node immune cells.

1. The effect of reactive oxygen species on equine sperm motility, viability, acrosomal integrity, mitochondrial membrane potential, and membrane lipid peroxidation. Baumber J et al. (2000). *J. Androl.* 21: 895-902.

2. N-acetylcysteine (NAC) decreases severity of acute pancreatitis (AP) in mice. Demols A et al. (2000) *Pancreas* 20: 161-69.

3. Thyroid-specific gene expression is differentially influenced by intracellular glutathione level in FRTL-5 cells. Lonigro, R et al. (2000). *Endocrinology* 141: 901-909.

4. Risk factors for cortical, nuclear, and posterior sub-capsular cataracts. Delcourt C et al. (2000). *Am. J. Epidemiol.* 151: 497-504.

5. Changes in the antioxidant content of mononuclear leukocytes from mice with endotoxin-induced oxidative stress. Victor, VM et al. (2002). *Mol. Cell Biochem.* 229:107-111

## Bioxytech® Plasma GPx-EIA Assay


21014

*Immunoassay for Quantification of Human Plasma Glutathione Peroxidase Enzyme Expression*

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**


**Catalog Number** 21014

Glutathione peroxidases (GPx's) are selenoenzymes that catalyze the reduction of hydroperoxides ( $H_2O_2$  or  $ROOH$ ) in the presence of glutathione (GSH). Human blood contains, in addition to Se-GPx in erythrocytes, a plasma-specific glutathione peroxidase (pl-GPx). Glutathione peroxidase is a tetramer of approximately 94-100 kDA with each of the four identical subunits containing an active site with an essential selenocysteine residue. Plasma GPx differs from the other glutathione peroxidases by its primary sequence, its glycosylated N-terminal region, and its extra-cellular location. Initially purified from human plasma, pl-GPx has also been found in human milk. It has now been reported that pl-GPx is mostly synthesized and secreted by renal cells. pl-GPx may play a key role in the methionine sulfoxide reductase/thioredoxin system (see figure).

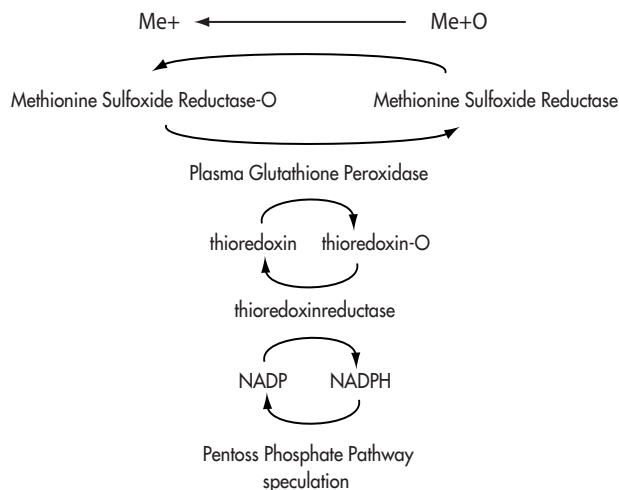

### pl-GPx-EIA Method

The Standard curve is obtained by fitting the standard absorbance at 405 nm to the concentration of pl-GPx by 4-parameter logistic curve fit.

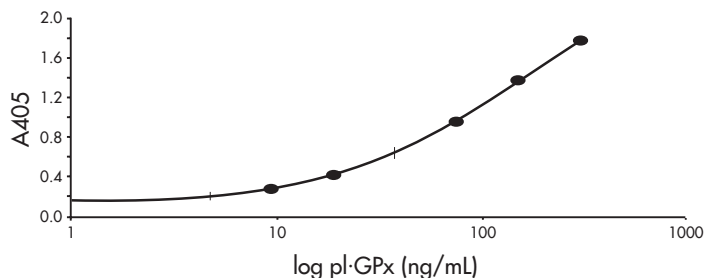

The Bioxytech pl-GPx-EIA assay system is a standard three (3) step "sandwich" ELISA. In step one, a plate bound (e.g. "solid phase") anti-human pl-GPx polyclonal antibody captures sample and/or standard pl-GPx. In step two, the bound pl-GPx antigen is detected using a biotin-labeled polyclonal anti-human pl-GPx. Finally, step three involves binding of a streptavidin/alkaline phosphatase conjugate to the biotinylated antibody. *p*-Nitrophenyl phosphate (pNPP) is then added causing the enzymatic hydrolysis of pNPP. The formation of yellow *p*-Nitrophenol is monitored at 405 nm and the amount of sample pl-GPx calculated against the color development of standards.

U.S. Patent Number US 5,861,262.

### References

1. Maddipati, K.R. and Marnett, L.J., (1987) *J. Biol. Chem.* 262:33, 17398-17403 Esworthy,
2. R.S., *Arch. Biochem. Biophys.* (1993) 307:1, 29-34.
3. OXIS unpublished results.

**pl•GPx-EIA Product Summary**

Intended Use:	Quantitative measurement of human plasma glutathione peroxidase. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	96 well ELISA
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 anti-pl•GPx Coated Microtiter Wells (6x16 strips) Plus Frame</li> <li>• Purified, Lyophilized pl•GPx Standard</li> <li>• Biotin Coupled Polyclonal Anti-pl•GPx</li> <li>• Streptavidin-Coupled Alkaline Phosphatase Solution</li> <li>• pNPP Tablets Plus Diluting Buffer</li> <li>• Stop Solution</li> <li>• Diluting and Wash Buffers</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Human serum or heparinized plasma.
Specificity:	No observed cross-reactivity with other GPx isoforms.
Sensitivity:	2.5 ng/mL

**pl•GPx-EIA Selected Research Applications**

Jorres et al. investigate the effect of repeated ozone exposure on inflammatory biomarkers. The OXIS plasma GPx-EIA is used to test for pl•GPx concentration in bronchioalveolar lavage fluid (BALF)

Delcourt et al studied the effect of sunlight exposure on human cataract pathogenesis.

Both physical and biological measurements (plasma glutathione peroxidase expression and superoxide dismutase activity, using the OXIS pl•GPx-EIA and SOD-525 assays respectively) were evaluated. Plasma was used for the pl•GPx-EIA and erythrocytes were tested for using the SOD-525 kit.

Wendland et al. studied plasma antioxidants, vitamins and lipid peroxidation among "healthy" patients and CD patients (smokers and malnourished subjects were eliminated from the study). Plasma GPx and plasma selenium concentrations were similar in both groups. They noted that previous reports, showing elevated plasma GPx in CD patients could be due to the involvement of smokers in those studies.

Atherosclerosis and cardiovascular complications are common features in end-stage renal failure patients (Glomerulonephritis, Renal hypoplasia, Polycystic kidney disease, Renal failure due to stasis and Bilateral nephrostomy). Donica et al. studied the relationships between lipid peroxidation products (MDA & 4-HNE) and plasma-GPx contents versus proinflammatory cytokines (TNF $\alpha$ , IL-6 and soluble receptor, sIL-6R) contents in plasma samples. Plasma-GPx levels was lower in patients, compared to the control, but not significant.

1. The effect of repeated ozone exposures on inflammatory markers in bronchioalveolar lavage (BAL) fluid and mucosal biopsies. Jorres, RA et al. (2000). *Am. J. Respir. Crit. Care. Med.* 161: 1855-61

2. Light Exposure and the risk of cortical, nuclear and posterior sub capsular cataracts. Delcourt, C et al. (2000). *Arch Ophthalmology* 118(3): 385-392.

3. Lipid peroxidation and plasma antioxidant micronutrients in Crohn's disease. Wendland, B.E., et. al., (2001) *Am. J. Nutr.* 74(2):259-264

4. Evaluation of lipids peroxidation products versus proinflammatory cytokines in hemodialized patients. Donica, H. (2001) *Renal Failure* 23(2): 231-38

Bioxytech® cGPx-340 Assay


21017

Quantitative Measurement of Cellular Glutathione Peroxidase Enzyme Activity

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21017


Cellular glutathione peroxidase (cGPx) is a member of a selenium containing family of GPx enzymes whose function is to detoxify peroxides in the cell. Since peroxides decompose to form highly reactive radicals, the GPx enzymes play a critical role in protecting the cell from free radical damage, particularly lipid peroxidation. The GPx enzymes catalyze the reduction of H<sub>2</sub>O<sub>2</sub> to water and organic peroxides (ROOH) to the corresponding stable alcohols (ROH) using glutathione (GSH) as a source of reducing equivalents.

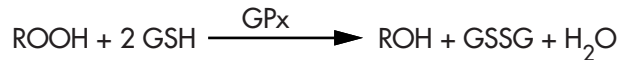


With few exceptions, GPx enzymes are comprised of 4 identical subunits. Each subunit contains a molecule of selenocysteine in the enzyme's active site. When associated with GSH, GPx selenocysteine is thought to participate directly in electron donation to the peroxide substrate and to become oxidized in the process. This oxidation of selenocysteine then strips electrons from 2 units of GSH generating, disulfide oxidized glutathione GSSG. Since cGPx activity depends on selenium the Bioxytech cGPx-340 assay is an ideal functional assay for detecting the effect of Se specific trace mineral deficiency.

**cGPx-340 Method**

The Bioxytech cGPx-340 method is an indirect measure of cGPx. Oxidized glutathione (GSSG), produced upon the reduction of organic peroxides by cGPx, is recycled to its reduced state by the enzyme glutathione reductase (GR).

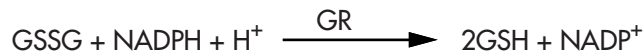

To assay for GPx activity, samples are added to a solution containing GSH, glutathione reductase (GR) and NADPH. An enzymatic reaction is initiated by addition of *tert*-butyl hydroperoxide substrate. By monitoring the NADPH consumption, as noted by a decrease in absorption at 340 nm, the Bioxytech cGPx-340 method allows for sensitive detection of cGPx activity. The graph below illustrates the usefulness of the Bioxytech assay as a functional enzyme biomarker. In this example, blood samples were taken from auction yard feeder calves and assayed for cGPx activity using the Bioxytech cGPx-340 test. After shipment to the feedyard, calves were monitored for signs and treatment of bovine respiratory disease complex (BRD). The figure below shows correlation between pre-shipment cGPx values and predisposition for BRD based on symptoms and incidence of treatment.

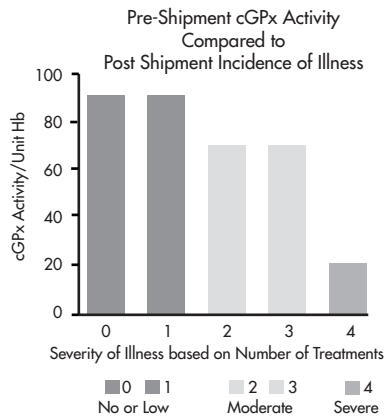

**References**

1. Paglia, D.E., Valentine, W.N., Studies on the quantitative and qualitative characterization of erythrocyte glutathione peroxidase. J. Lab. Clin. Med. July (1967) 70:158-169.

**cGPx-340 Product Summary**

Intended Use:	Quantitative measurement of cellular glutathione peroxidase activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• NADPH Reagent <i>Contains NADPH, GSH and GR</i></li> <li>• Assay Buffer</li> <li>• <i>tert</i>-Butyl Hydroperoxide</li> </ul>		
Storage and Stability:	21 months from date of manufacture when stored as specified.		
Specimen Requirements:	Cell and tissue lysates		
Assay Precision:	<u>Low</u>	<u>Med</u>	<u>High</u>
Inter- assay (% CV)	4.0	4.1	4.9
Sensitivity:	6 mU/mL ( <i>final assay concentration</i> )		

**cGPx-340 Selected Research Applications**

GPx-1 or cellular-GPx has been identified and studied in erythrocytes, lens, liver, brain, heart, platelet, lungs and other tissues. The cGPx-340 Assay has been used in research to study various conditions including ischemia/reperfusion injury, diabetes, neurological disorders, cystic fibrosis, cancer, preeclampsia, gastrointestinal, metal toxicity, and others.

Transplantation of any organ necessarily involves an unavoidable period of ischemia. In a research study using the OXIS cGPx-340 assay, Seth et al. showed that treatment with Picroliv (a natural antioxidant) protects rat renal tissues from ischemia-reperfusion injury by modulating antioxidant activities such as cGPx (1). Janciauskiene et al. examined cGPx levels in relation to LDL receptor negative and LDL receptor positive cell lines in a study targeting the detrimental cardiovascular effect of amyloid fibril deposits (IAPP) in Type 2 diabetes. They showed that elevated cGPx levels correlated with decreased lipid peroxidation and hence minimizing IAPP formation (2). Using a rat model, Parikh et al. demonstrated that some antipsychotic drugs (e.g., atypical antipsychotics: Clozapine, Risperidone, and Olanzapine) reduced oxidative stress and oxidative cell injuries by increasing the levels of antioxidants such as GPx, catalase and Cu/Zn-SOD and decreasing lipid peroxidation compared to the typical drug treatment, Haloperidol (3). Cystic Fibrosis is caused by the loss or dysfunction of the CFTR/Cl<sup>-</sup> channel. The resulting defective ion permeability leads to viscous epithelial secretion. In general, energy requirements of CF patients are 120-150% higher than the "normal" requirement, despite the administration of pancreatic enzyme supplement and others. However, it has been shown that a high fat diet without "proper" antioxidant supplementation may lead to higher degree of oxidation (e.g. lipid peroxidation) as well as endogenous antioxidant "depletion" such as GSH and cGPx (4). A research study in a murine model of somatic mutagenesis (cancer), using various OXIS kits (GPx-340, Catalase-520 and GSH/GSSG-412), the Felix group demonstrated that glucose-6-phosphate dehydrogenase (G6PD) is required for limiting oxidative mutagenesis in the mouse spleen (5). Preeclampsia is a leading cause of maternal and neonatal mortality and morbidity. Okatani et al. showed that melatonin could be useful in treating this condition and possible other clinical states involving excessive free radical production, such as intrauterine fetal growth retardation and fetal hypoxia (6). They showed GPx and SOD were significantly increased in chorionic homogenates from treated women, compared to the controls.


1. Prevention of renal ischemia-reperfusion-induced injury (IRI) in rats by Picroliv. Seth, PS et al. (2000) *Biochem. Pharmacol.* 59: 1315-22
2. Fibrillar islet amyloid polypeptide differentially affects oxidative mechanisms and lipoprotein uptake in correlation with cytotoxicity in two insulin-producing cell lines. Janciauskiene S and Bo Ahren (2000). *Biochem. BioPhy. Res. Com.* 267: 619-25.
3. Differential effects of antipsychotics on expression of antioxidant enzymes and membrane lipid peroxidation in rat brain. Parikh, V. et al.(2003) *J. Psychiatric Research* 37: 43-51
4. Increased plasma fatty acid concentration after respiratory exacerbations are associated with elevated oxidative stress in cystic fibrosis. Wood LG et al. (2002). *Am. J. Clin. Nutr.* 75: 668-75.
5. Redox imbalance and mutagenesis in spleens of mice harboring a hypomorphic allele of *Gpdx* encoding glucose 6-phosphate dehydrogenase. Felix, K et al.(2003) *Free Radical Biology & Medicine* 34 (2): 226-232.
6. Melatonin stimulates glutathione peroxide activity in human chorion. Okatani Y et al. (2001). *J. Pineal Res.* 30: 199-205




*Quantitative Measurement of Glutathione Reductase Enzyme Activity***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21018

Glutathione reductase (GR, EC 1.6.4.2) is a ubiquitous enzyme, which catalyzes the reduction of oxidized glutathione (GSSG) to glutathione (GSH). Glutathione reductase is essential for the glutathione redox cycle that maintains adequate levels of reduced cellular GSH. GSH serves as an antioxidant, reacting with free radicals and organic peroxides, in amino acid transport, and as a substrate for the glutathione peroxidases and glutathione S-transferases in the detoxification of organic peroxides and metabolism of xenobiotics, respectively. This homodimeric enzyme is a member of the family of flavoprotein disulfide oxidoreductases. Each subunit has four domains; beginning at the N-terminus: an FAD-binding domain, an NADPH-binding domain, a central domain, and an interface domain. The active site of GR is at the dimeric interface. Since the GSSG binding site is composed of residues from both subunits, only the dimeric form is active.

Oxidized glutathione is reduced by a multi-step reaction in which GR is initially reduced by NADPH forming a semiquinone of FAD, a sulfur radical and a thiol. The reduced GR ( $GR_{red}$ ) reacts with a molecule of GSSG, resulting in a disulfide interchange, which produces a molecule of GSH and the  $GR_{red}$ -SG complex. An electron rearrangement in  $GR_{red}$ -SG results in a second disulfide interchange, splitting off the second molecule of GSH and restoring the GR to the oxidized form.

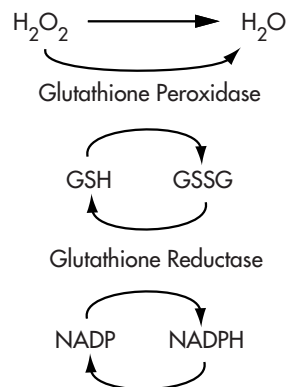**GR-340 Method**

The GR-340 assay is based on the oxidation of NADPH to  $NADP^+$  catalyzed by a limiting concentration of glutathione reductase.

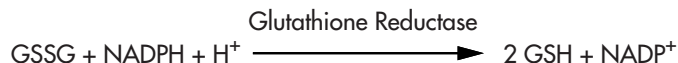

One GR activity unit is defined as the amount of enzyme catalyzing the reduction of one micromole of GSSG per minute at pH 7.6 and 25°C. As shown in the above reaction, one molecule of NADPH is consumed for each molecule of GSSG reduced. Therefore, the reduction of GSSG is determined indirectly by the measurement of the consumption of NADPH, as demonstrated by a decrease in absorbance at 340 nm (A340) as a function of time.

**References**

1. Bashir A, et al, *Biochem J* (1995) 312, 527–533.
2. Massey V, Williams CH, *J Biol Chem* (1965) 240(11), 4470–4480.
3. Worthington DJ, Rosemeyer MA, *Eur J Biochem* (1976) 67, 231–238.
4. García-Alfonso C, et al, *Int. J. Biochem* (1993) 25(1), 61–68.
5. Carlberg I, Mannervik B, *Meth Enzymol* (1985) 113, 485–490.
6. Peindo J, et al, *Molec Cell Biol* (1991) 101, 175–187.
7. García-Alfonso C, et al, *Int. J. Biochem* (1993) 25(4), 513–520.
8. Smith I, et al, *Anal Biochem* (1988) 75, 408–413.
9. Beutler E, *J Clin Invest* (1969) 48, 1957–1966.

**GR-340 Product Summary**

Intended Use:	Quantitative measurement of Glutathione Reductase activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• NADPH</li> <li>• Oxidized Glutathione (GSSG) in Potassium Phosphate Buffer.</li> <li>• Potassium Phosphate Buffer (K<sub>2</sub>PO<sub>4</sub>)</li> <li>• Diluent Buffer</li> </ul>		
Storage and Stability:	18 months from date of manufacture when stored as specified.		
Specimen Requirements:	Tissue homogenates and/or cell lysates.		
Assay Precision:			
n=20	<u>Low</u>	<u>Med</u>	<u>High</u>
Mean mU/mL	2.36	4.60	9.12
SD Total	0.22	0.40	0.72
Total (%CV)	9.20	8.60	7.80
Assay Range:	Minimum Net Rate = 0.0050 A <sub>340</sub> /Min (10x typical blank rate) Maximum Net Rate = 0.0625 A <sub>340</sub> /Min Corresponding to 0.8-10.0 mU SOD-525		
Sensitivity:	0.14 mU/mL (final concentration in the assay)		

**GR-340 Selected Research Applications**

The GR-340 assay has been used for many purposes including the study of emphysema related functional muscle impairment, Cystic Fibrosis related research and various liver toxicology studies.

Skeletal muscle functional impairment has been reported in emphysema (EMP) and COPD patients. Mattson et al. looked at markers of oxidative stress in skeletal muscles in a porcine elastase induced hamster model of emphysema. They showed that EMP may induce oxidative stress in peripheral skeletal muscle. The study reported elevated GR activities in gastrocnemius muscle of hamsters with emphysema compared to controls (1). Velsor et al studied the levels of endogenous antioxidants in Wild type vs CFTR knockout mice. They reported elevated levels of GR activity as measured with the OXIS GR-340 kit (2). Christova et al. looked at the effect of Cobalt induced oxidative stress on liver antioxidant enzymes. They showed that GR activities increase along with GSH and heme-oxygenase-1 under chronic cobalt exposure conditions (3). In a separate publication by the same author, acute effects of cobalt administration were compared in rat and guinea pig. Again they showed significant remodeling of the antioxidant biomarkers, including GR, as a result of acute cobalt induced oxidative stress.

1. Lipid peroxidation in the skeletal muscle of hamsters with emphysema. Mattson, J.P. et. al., (2002) Pathophysiology 8:215-221.
2. Antioxidant imbalance in the lungs of cystic fibrosis transmembrane conductance regulator protein mutant mice. Velsor et al. (2001) Am. J. Physiol. Lung Cell Mol. Physiol. 281:L31-38.
3. Heme oxygenase (HO) is the main protective enzyme in rat liver upon 6 day administration of cobalt chloride. Christova et al. (2001) Arch. Toxicology 75: 445-451.
4. Enhanced heme oxygenase (HO) activity increases the antioxidant defense capacity of guinea pig liver upon acute cobalt chloride loading: comparison with rat liver. Christova et al. (2002) Comparative Biochem, & Physiol. Part C 131:177-184.

## Bioxytech® GSH-420 Assay

21023


## Quantitative Measurement of Total Glutathione

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21023

Reduced glutathione is a ubiquitous tripeptide ( $\gamma$ -glutamyl-cysteinylglycine) which functions as a coenzyme in amino acid transport, detoxification of xenobiotics and carcinogens, synthesis of DNA precursors and as an antioxidant. The Bioxytech® GSH-420 is the next generation of the popular GSH-400 assay for the quantitative, colorimetric determination of glutathione in biological samples. This new assay allows for measurement of "total" GSH.

The GSH-420 compares favorably with other commercially available assay methods. The figure to the right shows the correlation between the GSH-420 and the DTNB-enzyme (Glutathione Reductase) recycling assay which is used in the GSH/GSSG-412 assay from OxisResearch®. The correlation coefficient obtained by linear regression is 0.9874.

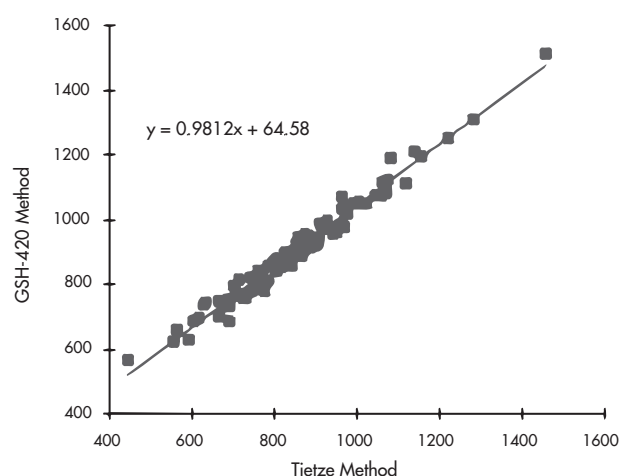

## GSH-420 Method

This assay is based on the formation of a chromophoric thione. There are three major steps to the reaction:

1. Sample is buffered and the reducing reagent tris (2-carboxyethyl) phosphine (TCEP) is added to reduce any oxidized glutathione (GSSG) present in the sample.
2. The chromagen, 4-chloro-1-methyl-7-trifluoromethyl-quinolinium methylsulfate, is added to the sample forming thioethers with all thiols.
3. The pH is raised by base addition resulting in a  $\beta$ -elimination reaction and formation of a chromophoric thione specific to GSH.


The measurement is made at 420 nm to negate the contribution by the thioether pool having peak absorbance at 356 nm. The figure below shows characteristic spectra generated after addition of base.

Improvements over existing GSH-400 assay:

- No ambiguity in "total" vs. "reduced" measurement.
- Greater specificity
- Precipitation reagent included

\* Note: Omission of reducing agent allows for approximation of reduced GSH as with the existing GSH-400 assay.

U.S. Patent Number 5,817,520

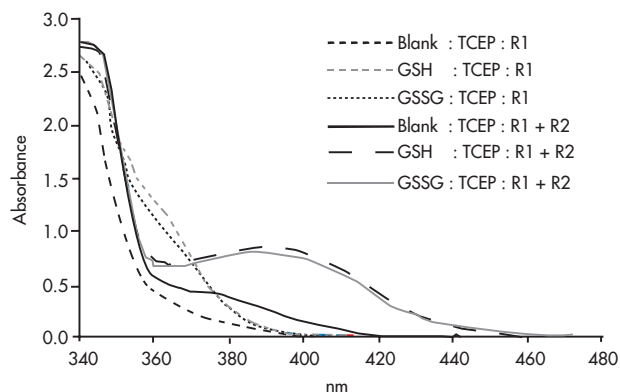

## References

1. Meister, A. and M.E. Anderson, (1983) *Glutathione*, Annual Review of Biochemistry, 52, 711-760.
2. Michelet, F. et al (1995), *Blood and Plasma Glutathione Measured in Healthy Subjects by HPLC: Relation to Sex, Aging, Biological Variables, and Life Habits*, Clinical Chemistry, 41(10), 1509-1517.
3. French Patent No. 9115868, United States Patent 5,817,520.
4. Burns, J.A. et al, (1991) *Selective Reduction of Disulfides by Tris(2-carboxyethyl)phosphine*, Journal of Organic Chemistry, 56, 2648-2650.
5. Richie Jr., J.P. (1996) et al, *Blood Glutathione Concentrations in a Large-Scale Human Study*, Clinical Chemistry, 42(1), 64-70.

**GSH-420 Product Summary**

Intended Use:	Quantitative measurement of "total" glutathione. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• Assay Buffer</li> <li>• Chromogen</li> <li>• Color Developer</li> <li>• Reducing Reagent</li> <li>• Precipitation Reagent</li> <li>• Calibrators (low and high)</li> </ul>		
Storage and Stability:	21 months from date of manufacture when stored in the dark at 2-4°C.		
Specimen Requirements:	Erythrocyte lysates, whole blood, tissue homogenates		
Precision:	<u>Low</u>	<u>High</u>	<u>Blood</u>
Mean	124.3	248.6	687.7
Intra-assay (% CV)	2.3	1.6	2.6
Inter-assay (% CV)	5.8	3.5	5.8
Total Precision (% CV)	6.0	3.7	6.1
Sensitivity:	9 µM in the sample		

**GSH-420 Selected Research Applications**

The GSH-420 has not been on the market as long as GSH-400, however it has already been used in a wide variety of studies.

In an animal feed supplement study, Williams et al. evaluated the safety and antioxidant potential of Lipoic acid (LA), a water-soluble and fat-soluble antioxidant, in horses undergoing light voluntary pasture exercise. They tested for endogenous antioxidant status (i.e. GSH and cGPx) using the GSH-420 and cGPx-340 kits on the automated OxyScan™ Analyzer. Lipid hydroperoxides were also evaluated using the OXIS LPO-560 Assay. They showed that LA supplementation of 10 mg/kg did not result in any adverse signs/effects over a 14-day period and moderately reduced the oxidative stress of horses that were allowed light activity (1).

In a performance animal related study Hargreaves et al. tested the antioxidant status in horses during prolonged and strenuous endurance exercise (2). They showed both erythrocyte GSH and Vitamin C were decreased during exercise. Similarly cellular GPx concentration tested with the OXIS GPx-340 assay was 4-fold higher before the exercise. They showed that changes of antioxidant status such as cellular enzymatic antioxidants (e.g. cGPx) are good indicators of muscle leakage associated with muscle damages and hypo-dehydration.

In one of the more innovative uses of our GSH-420 assay Downs et al. studied oxidative stress in corals and showed that coral (Zooxanthellae) bleaching may be a final strategy to defend coral from oxidative stress (3). They showed high levels of antioxidant enzymes and heat shock proteins were negatively correlated with levels of oxidative damage products.

1. Lipoic acid as an antioxidant in mature thoroughbred geldings: A preliminary study. Williams, C.A. et. al., (2002). J. Nutr. 132: 1628S-1631S
2. Antioxidant status of horses during two 80-km endurance races. Hargreaves, B.J. et al., (2002) J. Nutr. 132:1781S-1783S
3. Oxidative stress and seasonal coral bleaching. Downs, CA et al., (2002). Free Radic. Biol. Med. 33(4):533-543

Bioxytech® GSH/GSSG-412 Assay

21040


Quantitative Measurement of Total Oxidized Glutathione Ratio

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21040

OxisResearch® has developed an improved method to measure the ratio of reduced to oxidized glutathione (GSH/GSSG). This method features a more efficient mercaptan scavenger, 1-methyl-2-vinyl-pyridinium trifluoromethane-sulfonate (M2VP), to trap GSH. The figure below compares the scavenging abilities of the commonly used 2-vinylpyridine (2VP) with the OXIS M2VP reagent. One can easily see that M2VP is many times more effective in removing reduced GSH (99% within 1 minute) when compared to the 2VP reagent (11.7% within 1 minute).



The increased scavenging capability of the OXIS M2VP reagent allows for more accurate measurement of small amounts of GSSG without overestimation.

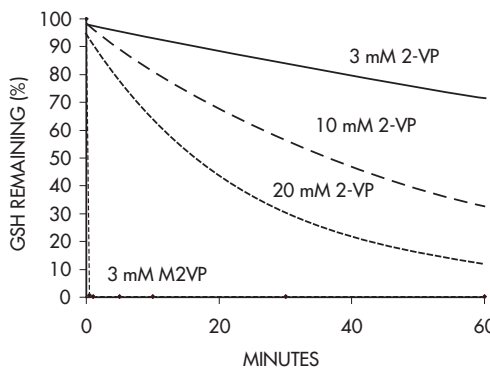

GSH/GSSG-412 Method

The test method is based on the DTNB-GSSG reductase recycling assay for GSH first described by Tietze in 1969. The method employs glutathione reductase to ensure all GSH is in the reduced form capable of reacting with DTNB (also called Ellman's reagent) to form a spectrophotometrically detectable product at 412 nm.

The data below represents a study that was done using the OXIS GSH/GSSG-412 assay. During a segmented charity run, whole blood was sampled from healthy OXIS volunteers. Blood was collected, with and without the OXIS M2VP GSH scavenging reagent, and tested with the GSH/GSSG-412 assay. Note the decreasing ratio as subject undergoes progressively more strenuous exercise. Also note the ratio recovery after a 1 hour rest period. This data represents the utility of the GSH/GSSG-412 assay in effectively assessing oxidative stress relative to the glutathione system.

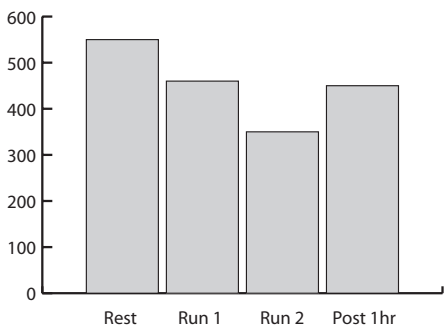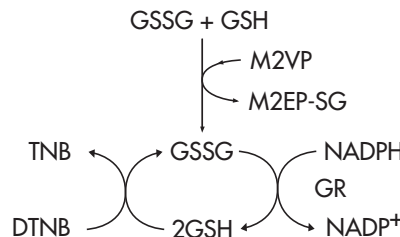

The GSH/GSSG-412 assay system contains reagents suitable for 100 tests representing the ratio of total glutathione to the oxidized form or 200 tests of either.

US Patent 5,543,298.

References

1. Tietze F (1969) Analytical Chemistry 27, 502-520.
2. Guntherberg H and Rost J (1966) Analytical Biochemistry 15, 205-210.
3. Griffith OW (1980) Analytical Biochemistry 106, 207-212.
4. Richie JP Jr., et al (1996) Clinical Chemistry 42, 64-70.
5. Anderson, M. (1996) Glutathione in Free Radicals, A Practical Approach, ed. N.A. Punchard and F.J. Kelly Oxford University Press, New York, p 213.

**GSH/GSSG-412 Product Summary**

Intended Use:	Quantitative measurement of total glutathione (GSH plus GSSG) and/or oxidized glutathione (GSSG) alone. This assay allows for determination of the GSH/GSSG ratio. For Research Use Only. Not Intended For Use in Diagnostic Procedures.			
Format:	100 test spectrophotometric cuvette (ratio) or 200 test spectrophotometric cuvette (GSH or GSSG)			
Kit Contents:	<ul style="list-style-type: none"> <li>• Assay Buffer</li> <li>• GSSG Buffer</li> <li>• Enzyme</li> <li>• NADPH</li> <li>• Scavenger</li> <li>• Chromogen</li> <li>• Standards</li> </ul>			
Storage and Stability:	12 months from date of manufacture when stored as specified			
Specimen Requirements:	Whole blood or tissue samples			
Precision:	GSSG in Buffer		Whole blood	
n=26	High	Low	GSH	GSSG
MeanRate at A <sub>412</sub>	0.4194	0.0427	0.2938	0.00490
Intra-assay (% CV)	0.59	3.50	0.96	6.45
Inter-assay (% CV)	2.84	2.96	3.11	7.61
Total Precision (% CV)	2.87	3.86	3.18	8.86
Sensitivity:	LLD in $\mu$ M reaction mixture = 0.009			
	LLD in $\mu$ M original sample = 0.54			

**GSH/GSSH-412 Selected Research Applications**

Our GSH assays have been and continue to be used in various research model systems including pulmonary diseases, alcoholism, fertility and cancer models. Using the GSH/GSSG-412 kit, Cavarra E et al. studied cigarette smoking associated oxidative stress on human protease enzymes (specifically anti-trypsin activity) in a mouse model of chronic obstructive pulmonary disease (COPD). They demonstrated that acute smoke exposure *in-vivo* can result in significant changes to the antioxidant capacity of bronchoalveolar lavage fluid (BAL). These changes were ascribed to depletion of many factors including the perturbation in the glutathione redox system as revealed by the increase in oxidized GSH (GSSG).

In an alcoholism and fertility study, the Emaluette group showed that when male rats were exposed to EtOH (8 wks of ethanol containing diet), there was increased paternal testicular oxidative injury demonstrated by decreased ratio of GSH/GSSG as well as enhanced lipid peroxidation (2).

In a diabetic model, Liang reported that metallothioneine-transfected diabetic mice exhibit significantly lower levels of GSSG compared to the non-transfected diabetic mouse and a similar to "normal"/control mouse (3). Many diabetic patients suffer from cardiomyopathy, even in the absence of vascular disease. Therefore, these results demonstrate that cardiomyocyte-specific expression of an antioxidant protein may reduce the damage to diabetic heart.

Most recently, researchers at the National Cancer Institute (NCI), National Institutes of Health (NIH) and a German group studied the effects of G6PD activity in terms of general metabolism, regulation of oxidative stress and cancer (somatic mutagenesis) in mice. They showed that the level of "total GSH" did not change significantly in the brain tissues of G6PD deficient (G6PD<sup>-</sup>) mice, compared to the "wild type". However, the GSH ratio (GSH/GSSG) was significantly increased (2-3 fold) in the "deficient" mice, as was their mutation rates (G6PD<sup>-</sup> =  $24 - 30 \times 10^{-5}$  vs. G6PD<sup>+</sup> =  $6.95 \pm 1.85 \times 10^{-5}$ ) (4). The mutagenesis was measured by examining the mutant frequency in Lac-Z gene, the target reporter gene of mutagenesis in the shuttle vector, pUR288.

1. Human SLPI (secretory leukoprotease inhibitor) inactivation after cigarette smoke exposure in new in vivo model of pulmonary oxidative stress. Cavarra E. et al. (2001). *Am. J. Physiol Lung Cell Mol. Physiol* 282: L412-L417.
2. Peripubertal paternal EtOH exposure: testicular oxidative injury, fertility and offspring. Emaluette NV et al., (2001) *Endocrine* 14: 213-19.
3. Over expression of metallothionein reduces diabetic cardiomyopathy. Liang Q et al. (2002) *Diabetes* 51 (1): 174-81.
4. Moderate G6PD deficiency increase mutation rates in the brain of mice. Felix, K et al. (2002) *Free Radical Biology & Medicine* 32 (7): 663-73

## Bioxytech® Catalase-520 Assay

21042

## Quantitative Measurement of Catalase Enzyme Activity

For Research Use Only, Not Intended For Use in Diagnostic Procedures


Catalog Number 21042

Catalase is a ubiquitous antioxidant enzyme that is present in most aerobic cells. It serves as one of the body's defense systems against  $H_2O_2$ , a strong oxidant that can cause intracellular damage. It is found in high concentration in erythrocytes and liver while lower concentrations are found in skeletal muscle, brain and heart. Measurement of catalase activity can be useful as a research tool for certain diseases such as acute pancreatitis and some liver diseases where values are elevated. Each unit of catalase decomposes 1  $\mu M$  of  $H_2O_2$  per minute at 25°C and pH 7.0. The common method for the determination of catalase involves measuring the change in absorbance at 240 nm (hydrogen peroxide decomposition due to catalase activity) in one minute. Criticisms of the method include interference from non-specific UV absorbing materials and the requirement of quartz cuvettes. Because of this, OxisResearch® has developed a colorimetric endpoint assay. The release of the new Catalase-520 assay adds to the popular Bioxytech® family of Antioxidant Biomarkers and provides another useful tool for oxidative stress investigations.


## Catalase-520 Method

The OxisResearch® Catalase-520 assay employs a two-step reaction scheme.


Reaction 1: Sample catalase reacts with a known quantity of  $H_2O_2$ . The reaction is stopped after 1 minute with sodium azide ( $NaN_3$ ).

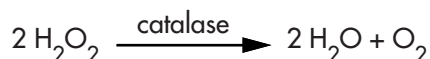

Reaction 2: In the presence of horseradish peroxidase (HRP), remaining  $H_2O_2$  reacts with 3,5-Dichloro-2-hydroxybenzenesulfonic (DHBS) acid and 4-aminophenazone (AAP) to form a chromophore with a color intensity inversely proportional to the amount of catalase in the original sample.

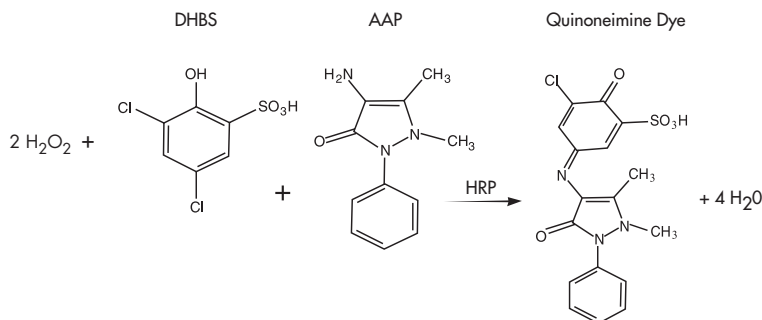

As a test of the assay's performance, Oxis compared red blood cell (RBC) catalase activity in a Thalassemia patient with that of a control population. In agreement with published values, the graph below shows increased activity in the patient relative to controls.

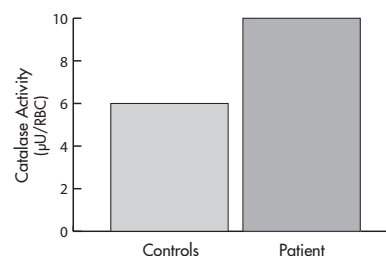

## References

1. Deisseroth, A. Dounce, A.L., (1970) Catalase: Physical and Chemical Properties, Mechanism of Catalysis, and Physiological Role, *Physiol. Rev.* 50, 319-375.
2. Zamocky, M. & Koller, F. (1999) Understanding the Structure and Function of Catalases: Clues from Molecular Evolution and In Vitro Mutagenesis, *Prog. Biophys. Mol. Biol.* 72, 19-66.
3. Kirkman, H.N., et al., (1987) The Function of Catalase-bound NADPH, *J. Biol. Chem.* 262, 660-666.
4. Aebi, H. & Suter, H. (1971) Acatlasemia, *Adv. Human Genet.* 2, 143-199.
5. Ogata, M. (1991) Acatlasemia, *Human Genet.* 86, 331-340.
6. Wakimoto, M., et al., (1998) Determination of Glutathione Peroxidase Activity and Its Contribution to Hydrogen Peroxide Removal in Erythrocytes, *Acta Med. Okayama* 52, 233-237.
7. Aebi, H. (1984) *Methods Enzymol* 105, 121-126
8. Fossati, P., et al. (1980) *Clin. Chem.* 26, 227-231

**Catalase-520 Product Summary**

Intended Use:	Quantitative measurement of catalase enzyme activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• Assay buffer</li> <li>• Substrate</li> <li>• Standards</li> <li>• Sample Diluent</li> <li>• Chromagen</li> <li>• HRP</li> <li>• Stop Reagent</li> </ul>		
Storage and Stability:	12 months from date of manufacture when stored as specified.		
Specimen Requirements:	Whole blood, serum, plasma, cell lysate and tissue		
Assay Precision:			
n=20	<u>Low</u>	<u>High</u>	<u>Cell Lysate</u>
Mean (U/mL)	14.1	56.7	38.7
Intra-assay (%CV)	2.39	0.78	1.29
Inter-assay (%CV)	4.03	2.13	5.55
Total Precision (%CV)	4.53	2.23	5.66
Sensitivity:	LLD in reaction mixture = 0.097 U/mL LLD in sample = 1.71 U/mL		

**Catalase-520 Selected Research Applications**

The OXIS Catalase-520 assay is relatively new but we are already beginning to collect information regarding its use in current research models. Here we list two such citations wherein the kit was used to monitor catalase activity in relation to NADPH availability and G6PD activity then in a second paper as related to anti-psychotic drug treatment.

Felix et al showed that catalase activity in the murine spleen is diminished in G6PD knockout mice as compared to controls (1). This correlates with catalase's dependency on NADPH to remain in the active state. Parikh et al studied endogenous antioxidant activities in the brain of Rats treated with various antipsychotic drugs and determined that chronic Haloperidol treatment had the effect of reducing Catalase, SOD and GPx activities as compared to animals treated with atypical antipsychotic drugs such as Clozapine, Risperidone and Olanzapine.

1. Redox imbalance and mutagenesis in spleens of mice harboring a hypomorphic allele of *Gpdx<sup>a</sup>* encoding glucose 6-phosphate dehydrogenase. Felix, K. et al. (2003) Free Radical Biology and Medicine 34 (2): 226-232.

2. Differential effects of antipsychotics on expression of antioxidant enzymes and membrane lipid peroxidation in rat brain. Parikh, V. et al. (2003) J. Psychiatric Research 37:43-51



Bioxytech® G6PD ; 6PGD-340 Assay


21045

Quantitative Measurement of Glucose-6-phosphate Dehydrogenase (G6PD) and 6-Phosphogluconate Dehydrogenase (6-PGD) Enzyme Activity

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21045

Since the amount of glutathione (GSH) present in the cell is finite, oxidized glutathione must be recycled to the reduced state to maintain protection against hydrogen peroxide. In cells experiencing oxidative stress, a continual supply of reducing equivalents, in the form of NADPH, is required in order to convert GSSG into GSH. The two reactions in the pentose phosphate pathway that produce NADPH are catalyzed by the enzymes glucose-6-phosphate dehydrogenase (G6PD) and 6-phosphogluconate dehydrogenase (6-PGD).


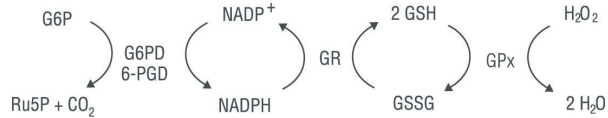


Of these two enzymes, G6PD is the predominant player in NADPH regulation and can play an important role in an organism's overall redox status and subsequent ability to handle oxidative stress. The easiest method for measuring G6PD activity is to monitor the production of NADPH in the presence of the appropriate substrate. A caveat of this method is that a product of the G6PD reaction produces substrate for the 6PGD reaction causing overestimation of G6PD when both G6PD and 6-PGD are present. In their failure to account for this contribution, other commercially available assays lack the specificity and subsequent sensitivity for studies relating G6PD activity and NADPH availability to oxidative stress. Because the G6PD ; 6PGD-340 assay from OxisResearch® does discriminate between the two enzymes, even subtle changes in G6PD activity can be detected.

G6PD ; 6PGD-340 Method

To avoid overestimation of G6PD and to obtain accurate measurement of both G6PD and 6-PGD activity, the OXIS G6PD ; 6PGD assay employs the method of Glock and McLean wherein two measurements of NADPH production are made: one with saturating concentrations of both glucose-6-phosphate (G6P) and 6-phosphogluconic acid (6-PGA) and one with only 6-PGA.

G6PD activity is then calculated as the difference in rate of NADPH production between the assays with both substrates and only 6-PGA. The figure below illustrates the over-estimation of red cell G6PD when both substrates are not used.

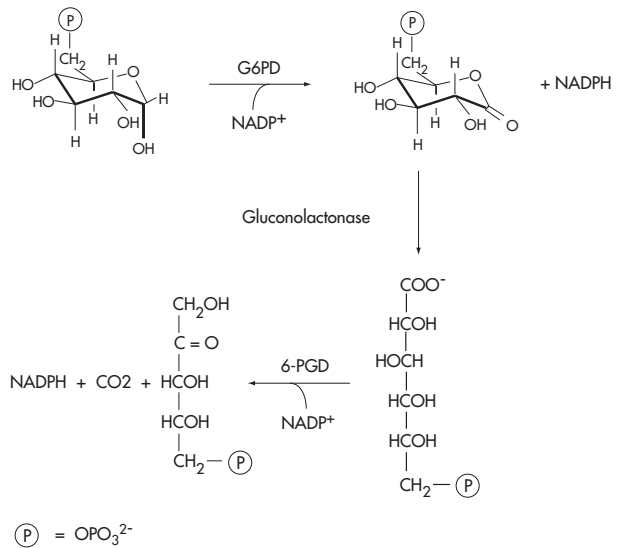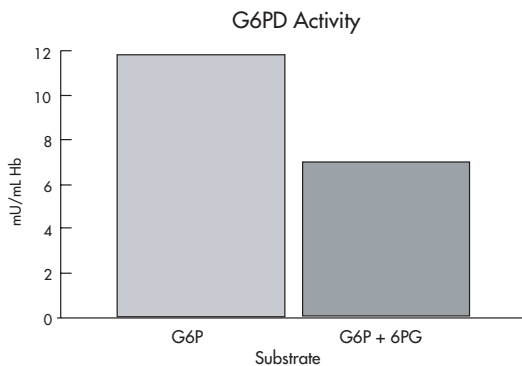

References

- Martini, G. & Ursini, M.V. (1996) A New Lease of Life for an Old Enzyme, *BioEssays* 18, 631-637.
- Beutler, E. (1993) Study of Glucose-6-Phosphate Dehydrogenase: History and Molecular Biology, *Am. J. Hematol.* 42, 53-58.
- Beutler, E. (1994) G6PD Deficiency, *Blood* 84, 3613-3636.
- Glock, G.E. & McLean, P. (1953) Further Studies on the Properties and Assay of Glucose-6-Phosphate Dehydrogenase and 6-Phosphogluconate Dehydrogenase of Rat Liver, *Biochem. J.* 55, 400-408.
- Beutler, E. (1984) *Red Cell Metabolism: A Manual of Biochemical Methods* pp. 68-71.

**G6PD ; 6PGD-340 Product Summary**

**Intended Use:** Quantitative measurement of G6PD and/or 6PGD enzyme activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.

**Format:** 100 test colorimetric (single enzyme)  
50 test colorimetric (both enzymes)

- Kit Contents:**
- Assay Buffer
  - NADP Reagent
  - G6P Substrate
  - 6-PGA Substrate
  - Sample Diluent

**Storage and Stability:** 12 months from date of manufacture when stored as specified

**Specimen Requirements:** Tissue, cell or RBC lysate

	G6PD			6-PGD		
	<u>Low</u>	<u>Med</u>	<u>High</u>	<u>Low</u>	<u>Med</u>	<u>High</u>
mU/mL enzyme	2.46	4.54	8.50	3.08	5.40	9.00
Mean ( $\Delta A_{340}/\text{min}$ )	0.0136	0.0403	0.0787	0.0192	0.0336	0.0560
Intra-assay (%CV)	2.07	1.14	2.45	1.65	1.38	1.30
Inter-assay (%CV)	1.62	1.05	2.42	1.84	1.25	1.42

<b>Sensitivity:</b>	<u>G6PD</u>	<u>6-PGD</u>
LLD in reaction mixture (mU/mL)	0.024	0.027
LLD in sample (mU/mL)	0.48	0.54

*Quantitative Measurement of Glutathione S-transferase Enzyme Activity***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21046

Glutathione S-transferase (GST) utilizes glutathione (GSH) to scavenge electrophilic xenobiotics as part of an organism's defense mechanism against the mutagenic, carcinogenic and toxic effects of such compounds. GST activity is present throughout the phylogenetic scale, generally as a family of enzymes whose major classes differ in their xenobiotic specificity. Each major class contains multiple isozymes. All GST isozymes catalyze the same reaction, shown schematically here, where X is a xenobiotic.



GST activity is present in most human tissues but the expression of the different isozymes is organ specific. It is especially prevalent in the liver, which plays a major role in detoxification. Due to its activity in detoxifying xenobiotics, GST plays an important role in the well-being of all living systems. Because of the interaction of GST with glutathione, a major component of the antioxidant system, and because of the implications relative to removal of oxidatively generated harmful substances, OxisResearch® has introduced the GST-340 assay for measurement of GST activity.

Because the various GST isozymes have widely varying  $K_m$  and  $V_{max}$  values towards different xenobiotics and GSH, the contribution of each to the overall GST activity will depend on the xenobiotic used as well as the reaction conditions. For this reason the substrate with the broadest range of isozyme detectability coupled with ease of monitoring was chosen for the GST-340 activity assay.

**GST-340 Method**

The Bioxytech® GST-340 assay is based on the GST-catalyzed reaction between 1-chloro-2,4-dinitrobenzene (CDNB) and GSH. Following addition of GST to the reaction vessel CDNB and GSH combine to form a dinitrophenyl thioether chromophore and a chloride ion (Figure 1).

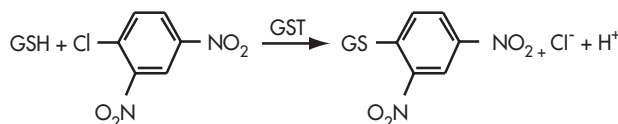
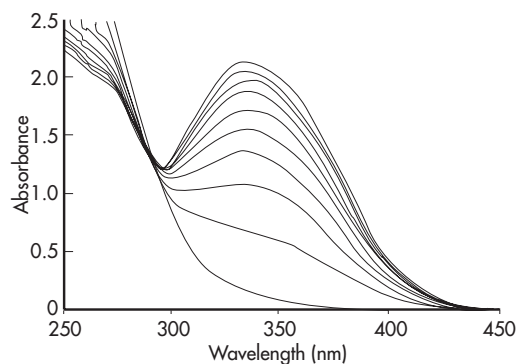


Figure 1 - The Reaction of CDNB with GSH

One unit of GST activity is defined as the amount of enzyme producing 1  $\mu\text{mol}$  of CDNB-GSH conjugate/min under the conditions of the assay. Figure 2 shows how addition of GST results in the gradual increase of the chromophore. The non-enzymatic reaction of CDNB and GSH under assay conditions is slow relative to the enzyme-catalyzed reaction.

Figure 2. (right) The reaction of CDNB and GSH in the presence of GST. The lowest scan at 340 nm is the CDNB/GSH mixture before addition of GST. Scans were taken at two minute intervals.

**References**

- Xie, C. et al. (2001) Expression of Glutathione S-Transferase Isozyme in the SY5Y Neuroblastoma Cell Line Increases Resistance to Oxidative Stress, *Free Rad. Biol. Med.* 31, 73-81.
- Hayes, J.D. & Strange, R.C. (2000) Glutathione S-Transferase Polymorphisms and Their Biological Consequences, *Pharmacology* 61, 154-166.
- Landi, S. (2000) Mammalian Class Theta GST and Differential Susceptibility to Carcinogens: A Review, *Mutation Res.* 463, 247-283.
- Galli, F. et al. (1999) Overexpression of Erythrocyte Glutathione S-Transferase in Uremia and Dialysis, *Clin. Chem.* 45, 1781-1788.
- Edwards, R. et al. (2000) Plant Glutathione S-Transferases: Enzymes with Multiple Functions in Sickness and in Health, *Trends Plant Sci.* 5, 193-198. (Additional references available upon request)

**GST-340 Product Summary**

Intended Use:	Quantitative measurement of glutathione S-transferase enzyme activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Specimen Requirements:	Tissue, cell or RBC lysate		
Storage and Stability:	9 months from date of manufacture when stored as specified		
Kit Contents:	<ul style="list-style-type: none"> <li>• Assay Buffer</li> <li>• CDNB Substrate</li> <li>• GSH</li> <li>• Sample Diluent</li> </ul>		
Precision:	<u>Low</u>	<u>Med</u>	<u>High</u>
mU/mL enzyme	1.42	4.20	8.20
Mean ( $\Delta A_{340}/\text{min}$ )	0.0136	0.0403	0.0787
Intra-assay (%CV)	0.89	0.80	1.95
Inter-assay (%CV)	1.05	0.89	2.13
Sensitivity:			
LLD in reaction mixture (mU/mL)	0.060		
LLD in sample (mU/mL)	1.2		

For Research Use Only, Not Intended For Use in Diagnostic Procedures


Catalog Number 21052

Four antioxidant strategies can be found in biological systems: (1) enzymatic (superoxide dismutase, glutathione peroxidase and catalase), (2) large molecules (albumin, ceruloplasmin and ferritin), (3) small molecules (ascorbic acid,  $\alpha$ -tocopherol,  $\beta$ -carotene, plasma ubiquinol, uric acid, and glutathione) and (4) some hormones (estrogen, melatonin, angiotensin and others). These antioxidants can be found either as water-soluble or fat-soluble molecules. OxisResearch® offers many assays for detection of specific, endogenously expressed enzymes and now offers the AOP-490 assay for detection of overall, antioxidant potential in aqueous biological samples.

This new assay compares well to the popular ORAC (Oxygen Radical Absorbance Capacity) assay without the need for fluorescence detection and requires only 3 minutes to perform compared to the 1 hour plus timeframe for ORAC. The AOP-490 assay also offers advantages over other existing assays such as FRAP (Ferric Reducing Ability) and TEAC (Trolox Equivalent Antioxidant Capacity) due to its ability to measure all antioxidant contributions (including thiol containing compounds) with little interference due to endogenous Reactive Oxygen Species (ROS).


### AOP-490 Method

The Bioxytech® AOP-490 Assay is based upon the reduction of  $\text{Cu}^{++}$  to  $\text{Cu}^+$  by the combined action of all antioxidants present in the sample. A chromogenic reagent, bathocuproine (2,9-dimethyl-4,7-diphenyl-1,10-phenanthroline), selectively forms a 2:1 complex with  $\text{Cu}^+$ . The resulting chromophore, having a maximal absorbance at 490 nm, provides the basis for quantification of antioxidant potential.

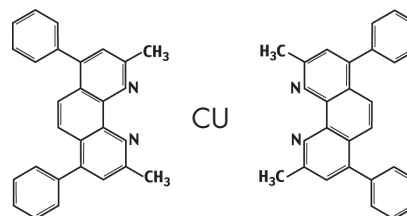

Using the standard provided, results of the assay may be expressed either as “ $\mu\text{M}$  Uric Acid Equivalents” directly from the plot of absorbance change versus mM uric acid concentration, or as “ $\mu\text{M}$  Copper Reducing Equivalents,” using a given conversion factor.

The Bioxytech® AOP-490 is capable of measuring of antioxidant potential in biological samples across a wide range of dilutions with excellent recovery.

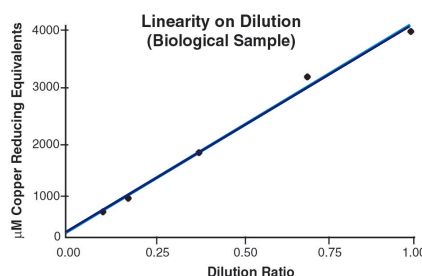

### References

1. Sies H, (1997) Oxidative stress: oxidants and antioxidants. Review. Exp. Physiol. 82: 291-5
2. Scott G (1997) Antioxidants in science, technology, medicine and nutrition. Coll House, UK Publishing.
3. Porter NA (1990) Auto-oxidation of polyunsaturated fatty acids: Initiation, propagation and product distribution (basic chemistry). Vigo-Pelfrey, C., ed Membrane lipid oxidation. Vol. 1 Boca Raton, FL, CRC Press.
4. Simic M, Karel M (1980) Auto oxidation in food and biological systems. New York, Plenum Press.
5. Winkler BS (1992) Unequivocal evidence in support of the nonenzymatic redox coupling between glutathione/glutathione disulfide and ascorbic acid/dehydroascorbic acid. Biochem. Biophys. Acta. 117: 287-90
6. Yamashita N, et al. (1998) Alpha-tocopherol induces oxidative damage to DNA in the presence of copper (III) ions. Chemical Res. Toxicol. 11: 855-858
7. Schilt AA (1966) Analytical application of 1-10-phenanthroline and related compounds. Pergamon Press, London, New York, Paris.
8. Prior RL, Cao P (1999) Antioxidant capacity and polyphenolic component of teas: implications for altering in vivo antioxidant status. Proc. Soc. Exp. Biol. Med. 220: 250-61
9. Ninfali P, Aluigi G (1998) Variability of oxygen radical absorbance capacity (ORAC) in different animal species. Free Radic. Res. 29: 299-408

**AOP-490 Product Summary**

Intended Use:	Quantitative measurement of antioxidant potential in biological samples. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	100 test colorimetric
Kit Contents:	<ul style="list-style-type: none"><li>• R1 (containing bathocuproine)</li><li>• R2 (containing Cu<sup>++</sup>)</li><li>• Uric Acid standard</li><li>• Stop solution</li></ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Serum, plasma, tissues and other biological samples
Assay Precision:	
Intra-assay (%CV)	2.22
Inter-assay (%CV)	4.20
Sensitivity:	30 µM uric acid equivalents

## Bioxytech® LPO-586 Assay

21012

## Quantitative Measurement of Malondialdehyde (MDA) Plus 4-hydroxyalkenals (4HAE)

For Research Use Only, Not Intended For Use in Diagnostic Procedures


Catalog Number 21012

Lipid peroxidation is a well-established mechanism of cellular injury in both plants and animals and is used as an indicator of oxidative stress in cells and tissues. Lipid peroxides are unstable and decompose to form a complex series of compounds including reactive carbonyl compounds. Polyunsaturated fatty acid peroxides generate malondialdehyde (MDA) and 4-hydroxyalkenals (HAE) upon decomposition. These substances in tandem provide a simple biomarker for measurement of general lipid peroxidation. The LPO-586 method is best suited for measuring MDA in combination with 4-hydroxyalkenals although the kit provides reagents for specific measurement of MDA alone as well. Researchers are advised that it is not possible to differentiate HAE from MDA using this kit due to slight differences in reaction stoichiometry when comparing the hydrochloric acid (HCl) reaction with the methanesulfonic acid (MSA) reaction.


We recommend this kit for detection of MDA plus HAE as a general indication of lipid peroxidation in tissue and cell samples. For MDA only we recommend the MDA-586 assay (21044) or for HAE only, we recommend the HAE-586 assay (21043). For measurement of lipid peroxidation in plasma we recommend researchers look at lipid hydroperoxides (LOOH) using the OXIS LPO-560 assay or MDA using the OXIS MDA-586 assay coupled with 3<sup>rd</sup> derivative spectroscopy.

## LPO-586 Method

The LPO-586 assay is based on the reaction of a chromogenic reagent, N-methyl-2-phenylindole (R1), with MDA and 4-hydroxyalkenals at 45°C. One molecule of either MDA or 4-hydroxyalkenal reacts with 2 molecules of reagent R1 to yield a stable chromophore with maximal absorbance at 586 nm.

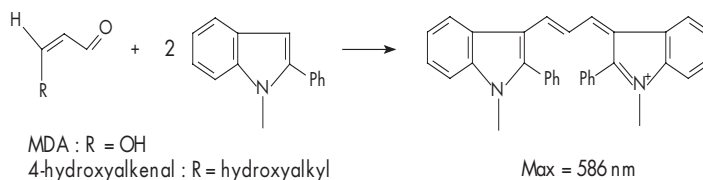

The graph below shows the relationship of LPO-586 lipid peroxidation measurement with other common methods including MDA and HNE by HPLC and lipid hydroperoxide (LOOH) using the OXIS LPO-560 method.

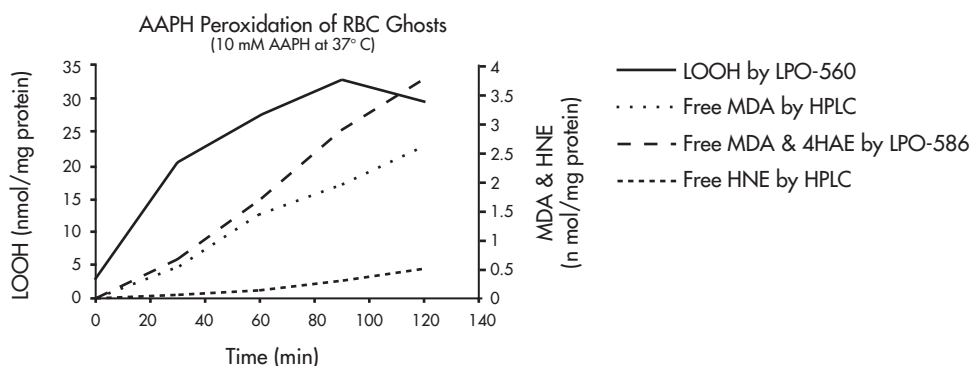

U.S. Patent Number 5,726,063.

## References

1. Esterbauer H., Schaur R.J., Zollner H. (1991) Chemistry and biochemistry of 4-hydroxynonenal, malonaldehyde and related aldehydes. *Free Radical Biology & Medicine* 11:81-128.

**LPO-586 Product Summary**

Intended Use:	Quantitative measurement of glutathione in biological samples. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	100 test colorimetric
Kit Contents:	<ul style="list-style-type: none"> <li>• R1 Chromogenic reagent</li> <li>• R2 Acid reagent</li> <li>• S1 Standard (4-HNE-diethylacetal)</li> <li>• S2 Standard (TMOP, Tetramethoxypropane)</li> </ul>
Storage and Stability:	24 months from date of manufacture when stored as specified.
Specimen Requirements:	Tissue homogenates and cell lysates.
Reproducibility:	Experiments in which standard samples (0-20 µM) were assayed using the same protocol over a period of 10 days established the standard error of the measurement (SEM) at less than 5%.
Sensitivity:	0.1 nmol/mL assay concentration 0.5 nmol/mL sample concentration

**LPO-586 Selected Research Applications:**

The ability of reactive oxygen species [e.g., hydroxyl, (OH<sup>•</sup>), peroxy (RO<sup>•</sup>)] and advanced lipid peroxidation end products (ALE) to react with amino acids/proteins, enzymes and DNA molecules can cause a broad range of pathophysiological conditions in plants, animals and humans. These conditions include fruit ripening disorders (plants), environmental toxicity, fertility, diabetes, cardiovascular disease (e.g. atherosclerosis), neurodegenerative diseases (e.g. Parkinson's and Alzheimer's), arthritis, cancer, inflammation, ischemia/reperfusion, stroke and organ transplantation.

Using the OXIS LPO-586 assay, the Jimenez group researched the oxidative process during tomato fruit ripening. Their data indicated that, under "off-the-vine" ripening conditions (tomatoes were harvested at mature green stage and left at room temperature to ripen), lipid peroxidation in the fruit was increased compared to the vine ripened fruit (1). Ethanol consumption has been reported to cause testicular oxidative injury as measured by lipid peroxidation in a rat model. The condition was positively correlated with decreased fertility and diminished performance of offspring (2). Patients with diabetes mellitus have high incidence of coronary heart disease (CHD) and over 50% of their mortality rate may be due to CHD. In a hyperlipidemic-diabetic hamster model, El-Swefy et al. reported that advanced lipid peroxidation end products (MDA & 4-HAE) were found in LDL and RBC of NIDDM and IDDM hamsters (3). Parkinson's disease is a progressive degenerative disorder of the central nervous system. Using cabergoline (a dopamine agonist), the Finotti group demonstrated in a rat model that lipid peroxidation can be reduced (4). In the brain, cholesterol metabolism can produce LDL and HDL is synthesized by glial cells and associated with amyloid plaques in Alzheimer's disease (AD). Thus, lipid peroxidation is a central feature of vascular illness and may have an important role in AD. Recently, the White group reported that the ability of amyloid precursor protein products from fruit fly, worm, frog, puffer fish, electric ray and humans to induce direct or oxidized lipoprotein mediated neuronal damage *in vivo* will depend on the availability of both Cu and lipoprotein (5). Phenobarbitone (PB) is used for preventing intraventricular hemorrhage, a major complication of pre-birth. The Hodges group used a rat model to show that PB could promote the formation of liver tumors indirectly due to the enhanced production of lipid peroxy radicals. They showed elevated lipid peroxidation levels positively correlated with transient induction of epidermal growth factor inducing S-phase DNA synthesis and subsequent cell proliferation (6). Reperfusion of ischemic tissues causes the generation of reactive oxygen species that contribute to tissue injury. The Ozaki group hypothesized that Rac1, (from the rho family of small GTP-ases) expression increases due to ischemia/reperfusion (I/R) induced necrosis and apoptosis during organ transplantation, myocardial infarction and stroke. Rac1 regulates the production of ROS by an NADPH oxidase in phagocytic and non-phagocytic cells. Using the Bioxytech® LPO-586 assay and others, they demonstrated that targeting inhibition of NADPH oxidase provides a new avenue for *in vivo* therapy aimed at protecting organs at risk from I/R (7).

1. Changes in oxidative processes and components of the antioxidant system during tomato fruit ripening. Jimenez, A. et al (2002). *Planta* 214: 751-58.
2. Peripubertal paternal EtOH exposure.: testicular oxidative injury, fertility and offspring. Emaluele NV et al., (2001) *Endocrine* 14: 213-19
3. The effect of Vitamin E, probucol and lovastatin on oxidative status and aortic fatty lesions in hyperlipidemic-diabetic hamster. El-Swefy S et al. (2000). *Atherosclerosis*. 149: 277-86.
4. Reduction of lipid peroxidation in different rat brain areas after cabergoline treatment. Finotti, N et al. (2000). *Pharmacol. Res.* 42 (4): 287-87
5. Contrasting species-dependent modulation of copper-mediated neurotoxicity by the Alzheimer's disease amyloid precursor protein. White, AR et al. (2002). *J. Neuroscience* 22 (2): 365-76.
6. Potentiation of Epidermal growth factor or EGF-induced DNA synthesis in rat hepatocytes by phenobarbitone (PB): possible involvement of oxidative stress and kinase activation. Hodges NJ et . (2000). *Carcinogenesis* 21: 2041-47.



## Bioxytech® 8-Isoprostane Assay


21019

Immunoassay For 8-epi-Prostaglandin  $F_{2\alpha}$ 

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21019

Prostaglandin  $F_{2\alpha}$  ( $PGF_{2\alpha}$ ) is formed naturally from arachidonic acid via the cyclooxygenase (COX I and II) mediated pathways. It functions as a potent vasoconstrictor *in vivo* under the control of COX I and II. Inhibition of COX activity and subsequent production of  $PGF_{2\alpha}$  provides the mechanism for the anti-inflammatory properties of common non-steroidal anti-inflammatory drugs (NSAIDs) such as aspirin, ibuprofen and others.

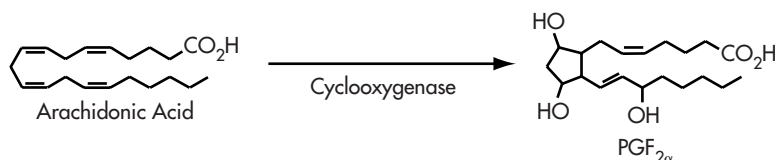


Isoprostanes are prostaglandin-like compounds that are produced upon peroxidation of arachidonic acid. They arise as a result of oxidative insult both *in vivo* and *ex vivo*. Because of this non-enzymatic, oxidative pathway, 8-epi-prostaglandin  $F_{2\alpha}$  (8-epi- $PGF_{2\alpha}$ ) has become a popular biomarker relating oxidative stress to various disease pathologies. Similar to the COX derived  $PGF_{2\alpha}$ , 8-epi- $PGF_{2\alpha}$  has been shown to have biological activity as a potent pulmonary and renal vasoconstrictor. The Bioxytech® 8-Isoprostane-EIA provides a simple method for determination of 8-isoprostane (8-epi- $PGF_{2\alpha}$ ) in biological samples such as tissue, plasma and urine.

## 8-Isoprostane-EIA Method


The 8-Isoprostane-EIA assay is a competitive enzyme-linked immunosorbent assay (ELISA) for determining levels of 8-epi- $PGF_{2\alpha}$  in biological samples. Briefly, samples or standards are added simultaneously with an unbound 8-epi-horseradish peroxidase (HRP) conjugate. Sample and standard 8-epi compete with 8-epi-HRP for binding to the stationary anti-8-epi bound to the plate wells. After washing, a Tetramethyl Benzidine (TMB) substrate is added, resulting in color development. Color intensity is proportional to the amount of 8-epi-HRP bound and inversely proportional to the amount of 8-epi in the samples or standards. The Bioxytech® 8-isoprostane-EIA has a high degree of correlation with the gold standard gas chromatograph mass spectrometry (GC/MS) methods.

The Bioxytech® 8-isoprostane EIA is recommended for use in tissue, serum or plasma samples and requires solid phase extraction using a  $C_{12}$  or silica column prior to sample assay. Although suitable for measurement of 8-isoprostane in urine, OXIS recommends our Urinary 8-isoprostane-EIA method (21048) for this purpose. A third alternative is our new Urinary 8-isoprostane metabolite assay (21049).

Comparison of isoprostane levels in human urine by GC/MS and ELISA

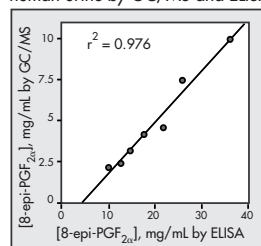

Typical Standard Curve

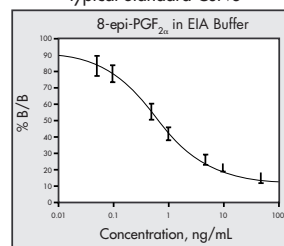

## References

1. Hazbun, M.E., et al. Cell Mol. Biol. (1993) 9:568-572.
2. Morrow, J.D., Harris, T.M., and Roberts, L.J., Analyt. Biochem., 184:1-10 (1990).
3. Morrow, J.D., Hill, K.E., Burke, R.F., Nammour, T.M., Badr, K.F., and Roberts, L.J., Proc. Natl. Acad. Sci. U.S.A. 87:9383-9387 (1990).
4. Morrow, J.D., and Roberts, L.J., Free Rad Biol. Med, 10:195-200 (1991).
5. Morrow, J., et al, Proc. Natl. Acad. Sci. USA 89:10721-10725 (1992).

**8-Isoprostane-EIA Product Summary**

Intended Use:	Quantitative measurement of 8-Isoprostane in biological samples. Also known as 8-epi-Prostaglandin F <sub>2α</sub> and 15-F <sub>2t</sub> -isoprostane. For Research Use Only. Not Intended For Use in Diagnostic Procedures.	
Format:	96 Test Colorimetric (ELISA)	
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 Well Anti-8-epi Coated Microtiter Plate</li> <li>• 8-epi Standard</li> <li>• Wash Buffer</li> <li>• Dilution Buffer</li> <li>• TMB Substrate</li> <li>• 8-epi-HRP Conjugate</li> <li>• Reagent Troughs</li> </ul>	
Storage and Stability:	12 months from date of manufacture when stored as specified.	
Specimen Requirements:	Tissue, plasma, and urine. (Recommended for plasma and tissue)	
Specificity:	Cross-reactivity at 50% B/B <sub>0</sub>	
	8-epi-PGF <sub>2α</sub>	100%
	9α,11b-Prostaglandin F <sub>2α</sub>	4.1%
	13,14-Dihydro-15-keto PGF <sub>2α</sub>	3.0%
	9b,11α-Prostaglandin F <sub>2α</sub>	<.01%
	Prostaglandin F <sub>2α</sub>	<.01%
	Prostaglandin E <sub>2</sub>	<.01%
	Prostaglandin D <sub>2</sub>	<.01%
	Arachidonic Acid	<.01%
Sensitivity:	0.05 ng/mL (precision best on steeper, linear portion of the curve)	

**8-Isoprostane-EIA Selected Research Applications**

The OXIS 8-isoprostane assay has been used to study a variety of diseases and parameters including circadian rhythms as related to uric acid and lipid peroxidation in diabetic and "healthy" men and the effects of cigarette smoking on oxidative stress parameters related to Chronic Obstructive Pulmonary Disease (COPD) related to emphysema. In other cardiovascular related studies, researchers have used the OXIS kit to look at the effects of certain lipid lowering therapies (e.g. HMG-CoA inhibitor/Cerivastatin) on plasma lipid peroxidation.

Kanabrocki, E.L., et. al., showed that oxygen-based free radicals, such as  $\cdot\text{OH}$ ,  $\text{O}_2^{\cdot-}$ , may be involved in the peroxidation of lipids, which was enhanced in the diabetic subjects, but that  $\text{ONOO}^-$  is unlikely to make a major contribution. Although higher levels of uric acid in the diabetics helped to scavenge out  $\text{ONOO}^-$ , it was ineffective in preventing lipid peroxidation (1).

SLPI, like  $\alpha_1$ -anti-proteinase loses its activity due to oxidative stress. This results in enhanced proteolysis of lung connective tissue (e.g., elastin), a key event in the pathogenesis of emphysema. Cavarra E. et al. showed that 8-isoprostane levels were correlated with decreased antitrypsin activity among smokers illustrating some of the manifestation of smoking induced oxidative stress (2). Lipid-lowering therapy, using HMG-CoA inhibitor/Cerivastatin decreases the risk of coronary events in both primary and secondary prevention. Tsunekawa T et al. showed that plasma 8-isoprostane was decreased with the treatment (3).

1. Circadian variation in oxidative stress markers in healthy and type II diabetic men. Kanabrocki, E.L., et. al., (2002) *Chronobiology International* 19(2):423-39

2. Human SLPI (secretory leukoprotease inhibitor) inactivation after cigarette smoke exposure in new in vivo model of pulmonary oxidative stress. Cavarra E.. et al. (2001). *Am. J. Physiol Lung Cell Mol. Physiol* 282: L412-L417

3. Cerivastatin, a hydroxymethylglutaryl coenzyme A reductase (HMG-CoA) inhibitor, improved endothelial function in elderly diabetic patients within 3 days. Tsunekawa T et al. (2001). *Circulation*. 104(4):376-379

Bioxytech® H<sub>2</sub>O<sub>2</sub>-560 Assay

21024


*Quantitative Measurement of Hydrogen Peroxide in Aqueous Solutions***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21024

Hydrogen peroxide is a powerful oxidant at the center of many redox pathways.

It is also a key player in the healing process as neutrophils gather at a wound site post trauma and release bactericidal reactive oxygen species (ROS) and H<sub>2</sub>O<sub>2</sub> to kill bacteria and prevent infection. The ongoing interest in understanding the role of hydrogen peroxide in biological systems including its role as a second messenger prompts the need for a rapid and quantitative measurement tool. Researchers also require a method of hydrogen peroxide measurement for laboratory reagents and detergents in order to prevent the unintended contribution of peroxidation by reagent use. The Bioxytech® H<sub>2</sub>O<sub>2</sub>-560 assay allows for simple colorimetric measurement of hydrogen peroxide.

**H<sub>2</sub>O<sub>2</sub>-560 Method**

The Bioxytech® H<sub>2</sub>O<sub>2</sub>-560 Assay is based on the oxidation of ferrous ions (Fe<sup>2+</sup>) to ferric ions (Fe<sup>3+</sup>) by hydrogen peroxide under acidic conditions, equation (1). The ferric ion binds with the indicator dye xylenol orange (3,3'-bis[N,N-di(carboxymethyl)-aminomethyl]-o-cresolsulfone-phthalein, sodium salt) to form a stable colored complex which can be measured at 560 nm, equation (2).

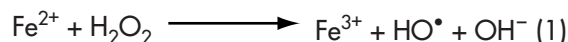

The Fe<sup>3+</sup>-xylenol orange complex has an extinction coefficient of 1.5 X 10<sup>4</sup> M<sup>-1</sup>·cm<sup>-1</sup> at 560nm when ferric ions are added in the absence of hydrogen peroxide. However, in the presence of sorbitol there is substantial chain oxidation of ferrous ion, increasing the color yield. This results in an apparent extinction coefficient of 2.67 X 10<sup>5</sup> M<sup>-1</sup>·cm<sup>-1</sup>, indicating that approximately 18 moles of Fe<sup>2+</sup> are oxidized to Fe<sup>3+</sup> for every mole of hydrogen peroxide present, increasing the sensitivity of the assay.

**H<sub>2</sub>O<sub>2</sub>-560 Product Summary**

Intended Use:	Quantitative measurement of hydrogen peroxide. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	100 Test Colorimetric
Kit Contents:	<ul style="list-style-type: none"> <li>• R1, Ammonium Iron Sulfate (Fe<sup>2+</sup>)</li> <li>• R2, Sorbitol/Xylenol Orange solution</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified. Working reagent stable 12 hours refrigerated.
Specimen Requirements:	Aqueous samples
Sensitivity:	1 µM in the assay

**H<sub>2</sub>O<sub>2</sub>-560 Selected Research Applications**

The OXIS H<sub>2</sub>O<sub>2</sub>-560 assay has been used to assess endogenous production of the pro-oxidant hydrogen peroxide as well as to monitor levels of exogenously applied H<sub>2</sub>O<sub>2</sub> in various research applications.

Joo et al. studied the relationship of gravitropism in plant roots to ROS production. They showed increased H<sub>2</sub>O<sub>2</sub> production in various regions of the root and suggested this could play a role in early gravitropic induced root response (1).

Kaul et al. found that the endogenous antioxidant activities (e.g. SOD, catalase and GSH) are not inhibited by rhinovirus infection, suggesting that oxidative stress is a result of increased production of pro-oxidants. The OXIS H<sub>2</sub>O<sub>2</sub>-560 kit was used to measure the H<sub>2</sub>O<sub>2</sub> production in cell culture supernatant as a means of monitoring pro-oxidant production (2).

LeVine et al. evaluated hydrogen peroxide production by macrophages in rat bronchoalveolar lavage (BAL) using the OXIS H<sub>2</sub>O<sub>2</sub>-560 assay (3). They utilized murine surfactant knockout models to show that both surfactant protein A and D play an apparent role in reducing inflammation as measured by decreased H<sub>2</sub>O<sub>2</sub> in SP-A and SP-D deficient mice.

Avshalumov et al. tested Long-Evans Rat Hippocampal slices for evoked potential in the presence of exogenously provided H<sub>2</sub>O<sub>2</sub>. Hydrogen peroxide was evaluated in the applied media using the OXIS H<sub>2</sub>O<sub>2</sub>-560 assay (4).

1. Joo JH et al. Role of auxin-induced reactive oxygen species in root gravitropism. *Plant Physiology*, (2002) 126:1055-1060.
2. Rhinovirus-induced oxidative stress and interleukin-8 elaboration involves p47-phox but is independent of attachment to intercellular adhesion molecule-1 and viral replication. Kaul, P et al. (2000). *J. Infec. Dis.* 181:1885-90.
3. Distinct effects of surfactant protein A or D deficiency during bacterial infection on the lung. Le Vine, AM et al. (2000). *J. Immunol.* 165:3934-40.
4. Mechanism underlying H<sub>2</sub>O<sub>2</sub>-mediated inhibition of synaptic transmission in rat hippocampal slices. Avshalumov MV et al., (2000). *Brain Research* (2000) 882: 86-94.


*Quantitative Measurement of Lipid Hydroperoxides (LOOH)***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21025

Lipid peroxidation (LPO), which is associated with oxidative stress, contributes to pathological processes in aging and many diseases such as atherosclerosis, diabetes, arthritis and Alzheimer's disease. Oxidation of polyunsaturated lipids involves an allylic hydrogen abstraction, insertion of molecular oxygen, and subsequent abstraction of hydrogen by the resulting hydroperoxyl radicals to form lipid hydroperoxides (LOOHs). The measurement of LOOHs are extensively used as an indicator of the LPO reaction. Some techniques, such as HPLC-chemiluminescence, thiobarbituric acid conjugation of MDA, iodide oxidation, and conjugated diene formation etc., have been used for the determination of LOOH. However, these methods either require sophisticated instrumentation or lack specificity and sensitivity. Therefore, the LPO-560 assay is designed to detect total LOOHs in plasma by a simple and sensitive method. An advantage of the LPO-560 assay is that total available plasma or serum LOOHs can be determined without involvement of a lipid extraction step.

**LPO-560 Method**

The design concept is based on the theory that hydroperoxides can be determined by the Ferrous Oxidation-Xylenol Orange (FOX) assay in conjunction with the reductant tris(2-carboxyethyl) phosphine (TCEP). Catalase is first used to decompose the existing hydrogen peroxide in biological samples.

The graph below shows the relationship of LPO-560 lipid peroxidation measurement with other common methods including MDA and 4HNE by HPLC and MDA and HAE using the OXIS LPO-586 method.

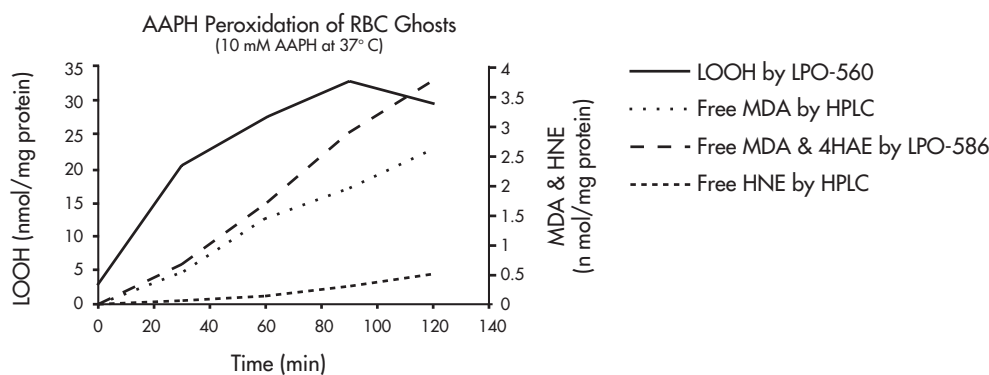**References**

1. Nourooz-Zadeh, J. et. al. *Anal Biochem* 220:403-409, 1994.
2. Nourooz-Zadeh, J. et. al. *Diabetologia* 40:647-653, 1997.
3. Wolff, S.P. *Methods in Enzymology* 223:182-189, 1994.
4. Tateishi, T. et. al. *Exp Gerontol* 22(2):103-111, 1987.
5. Anderson, D.J. *Clin Chem* 35:2152-2153, 1989.

**LPO-560 Product Summary**

Intended Use:	Quantitative measurement of total LOOH. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• Enzyme</li> <li>• Reducing Agent</li> <li>• Chromagen</li> <li>• Color Developer</li> <li>• Buffer</li> </ul>		
Storage and Stability:	24 months from date of manufacture when stored in the dark at 2° – 8° C.		
Specimen Requirements:	Heparinized plasma, serum or tissue homogenates		
Assay Precision:	<i>tert</i> -Butyl Hydroperoxide		
n=26	<u>Low</u>	<u>Med</u>	<u>High</u>
Average (µM)	1.94	3.80	10.0
Intra-assay (% CV)	3.64	2.82	2.41
Inter-assay (% CV)	4.67	6.08	3.10
Total Precision (% CV)	5.31	6.39	3.53
Sensitivity:	1.2 µM in original sample		

**LPO-560 Selected Research Applications**

This popularity of this assay has increased among researchers seeking an alternative marker other than the more commonly looked at malondialdehyde (MDA). While originally marketed as an acute LPO plasma assay, researchers have found the kit useful for samples as diverse as cell culture supernatant to coral polyps.

In a cardiovascular related study investigating the effect of remnant lipoproteins (RLPs) on plasma redox status Doi et al. evaluated levels of lipid hydroperoxides (LOOH) in cell culture supernatants. They showed that LOOH levels were dramatically increased in the cell culture supernatant of HUVECs incubated with RLPs compared to controls.

Williams et al. evaluated the safety and antioxidant potential of Lipoic acid (LA), a water-soluble and fat-soluble antioxidant, in horses undergoing light voluntary pasture exercise. They tested for endogenous antioxidant status (i.e. cGPx and GSH) using the OXIS cGPx and GSH-420 kits with the automated OxyScan™ Analyzer and lipid hydroperoxides using the OXIS LPO-560 Assay (3). They showed that LA supplementation of 10 mg/kg did not result in any adverse signs/effects over a 14-day period and moderately reduced the oxidative stress of horses that were allowed light activity.

Downs, et al. studied the oxidative stress in corals and showed that coral (Zooxanthellee) bleaching may be a final strategy to defend coral from oxidative stress. High levels of antioxidant enzymes and heat shock proteins were negatively correlated with levels of oxidative damage products including LOOH measured using the OXIS LPO-560 assay.

Wang et al. investigated the shortened lifespan of insect cell lines due to viral infection, a necessary part of recombinant protein expression. Using the LPO-586 assay they showed that oxidative stress is increased upon infection leading a postulation that lipid peroxidation can cause "leaky membranes" leading to cell death (4).

1. Remnant lipoprotein induce proatherothrombogenic molecules in endothelial cells through a redox-sensitive mechanism. Doi, H et al. (2000) *Circulation* 102: 670-6

2. Lipoic acid as an antioxidant in mature thoroughbred geldings: A preliminary study. Williams, C.A. et. al., (2002). *J. Nutr.* 132: 1628S-1631S

3. Oxidative stress and seasonal coral bleaching. Downs, CA et al.,. (2002). *Free Radic. Biol. Med.* 33(4):533-543

4. Evidence of oxidative stress following the viral infection of two lepidopteran insect cell lines. Wang et al. (2001) *Free Radical Biology and Medicine* 31:1448-1455.

## Bioxytech® 8-OHdG Assay

21026

## Immunoassay for 8-Hydroxy-2'-deoxyguanosine (8-OHdG)

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21026


In recent years a number of studies have correlated elevated levels of oxidative mediated DNA damage to various disease conditions. This has created a demand for easy to measure biomarkers that can be used to monitor oxidation of DNA in biological samples. 8-Hydroxy-2'-deoxyguanosine (8-OHdG) is a modified base that occurs in DNA after attack by reactive oxygen species (ROS), most notably hydroxyl (OH) radical. 8-OHdG modified bases are normally repaired by excision of the modified base resulting in formation of 8-hydroxy-2'-deoxyguanosine (8-OHdG).

8-OHdG can be identified through Southern blotting and tissue IHC. It can be found in plasma and in higher concentrations in the urine using HPLC or other means. For this reason it has become a very popular biomarker used to correlate oxidative stress with damage to DNA. The OXIS Kit allows for quantitative measurement of 8-hydroxy-2'-deoxyguanosine (8-OHdG) in tissue, serum, plasma, urine and cell culture supernatant.

## 8-OHdG-EIA Method

The Bioxytech® 8-OHdG-EIA kit is a competitive enzyme-linked immunosorbent assay (ELISA). Briefly, 8-OHdG antibody (anti-8-OHdG) is added to wells along with sample or standard. "Free" 8-OHdG in the sample or standards competes with "stationary" (plate bound) 8-OHdG for the anti-8-OHdG binding sites. Therefore, *increased* sample 8-OHdG concentrations leads to *decreased* plate bound anti-8-OHdG.

After "free" anti-8-OHdG/8-OHdG complexes are washed from the plate, the stationary 8-OHdG/anti-8-OHdG complexes are then treated with an enzyme labeled secondary antibody. Subsequent addition of a chromagenic substrate allows quantification of plate bound anti-8-OHdG. The resulting color is inversely proportional to the original sample 8-OHdG.

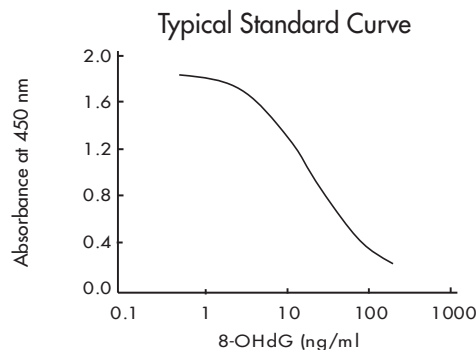

## References

1. R.G. Cutler, Am. J. Clin. Nutr. 50:373S-379S, 1991.
2. H. Kasai, P.F. Crain, Y. Kuchino, S. Nishimura, A. Ootsuyama and H. Tanooka. Carcinogenesis 7:1849-51, 1986.
3. C.G. Fraga, M.K. Shigenaga, J.W. Park, P. Degan and B.N. Ames. Proc.Natl.Acad.Sci.USA 87:4533-4537, 1990.
4. M.K. Shigenaga, and B.N. Ames. Free Radical Biology and Medicine 10:211-6,1991.
5. S. Okamoto and H. Ochi. Chemical Abst. 129859a, 1992.
6. S.S. Kantha, S. Wada, H. Tanaka, M. Takeuchi, S. Watabe and H. Ochi. Biochem.Biophys.Res.Comm. 233:278-82, 1996.
7. S. Toyokuni, T. Tanaka, Y. Hattori, Y. Nishiyama, A. Yoshida, K.Uchida, H.Hiai, H. Ochi and T. Osawa. Lab.Invest 76:365-74, 1997.
8. M. Erhola, S. Toyokuni, K.Okada, T. Tanaka, H. Hiai, H. Ochi, K. Uchida, T. Osawa, M.M. Nieminen, H. Alho and P.K. Lehtinen. FEBS Letters 40:287-91, 1997.
9. H. Ochi. Rouka seigyo syokuhin no kaihatsu (Development of foods for control of aging) Korin Press, p. 294, 1995.

**8-OHdG-EIA Product Summary**

Intended Use:	Quantitative measurement of 8-hydroxy-2'-deoxyguanosine. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	96 Test colorimetric (ELISA)
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 well 8-OHdG coated microtiter wells</li> <li>• Primary 8-OHdG antibody</li> <li>• Primary antibody dilution buffer</li> <li>• Secondary antibody</li> <li>• Secondary antibody dilution buffer</li> <li>• Chromagen</li> <li>• Chromagen dilution buffer</li> <li>• Wash buffer</li> <li>• Standard</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Serum, plasma, urine and DNA that has been extracted and digested.
Specificity:	Specific for DNA adduct 8-hydroxydeoxyguanosine. No cross-reactivity with other known 8-OHdG analogues was observed.
Sensitivity:	Threshold detection limit of 0.6 ng/mL

**8-OHdG-EIA Selected Research Applications**

The OXIS 8-OHdG ELISA assay kit has been used for a number of research applications in both human and veterinary related topics including studying the effect of antioxidant supplementation in high performance animals (e.g. sled dogs). In human studies the kit has been used to look at circadian variations in markers of oxidative stress in diabetic men as well as the effect of prolonged aerobic exercise on human immune function.

Baskin et al. showed that dietary antioxidant supplementation for exercising sled dogs can decrease levels of oxidative stress and subsequent levels of DNA damage as measured by decreased 8-OHdG in supplemented animals vs increased 8-OHdG in control animals (1). Kanabrocki et al used urinary 8-OHdG as an index of peroxynitrite mediated damage to DNA. They showed that oxygen-based free radicals, such as  $\cdot\text{OH}$ ,  $\text{O}_2^{\cdot-}$ , may be involved in the peroxidation of lipids which was enhanced in the diabetic subjects, but that ONOO<sup>-</sup> (causes oxidative damage in DNA or 8-OHdG pools) is unlikely to make a major contribution. Although higher levels of uric acid in the diabetics helped to scavenge out ONOO<sup>-</sup> as measured by decreased 8-OHdG, it was ineffective in preventing lipid peroxidation. Overall there appeared to be a statistically significant circadian variation of fluctuating baseline 8-OHdG levels and uric acid levels (i.e elevated uric acid levels correspond to decreased basal 8-OHdG levels) (2). Tsai et al. showed that 8-OHdG is increased during prolonged human aerobic exercise (e.g. as measured in marathon runners) and remains elevated for up to a week post exercise. The authors note that this may be responsible in part for exhaustive related immune dysfunction (3).

1. Effects of dietary antioxidant supplementation on oxidative damage and resistance to oxidative damage during prolonged exercise in sled dogs. Baskin CR et al. (2000) Am. J. Vet Res. 61(8): 886-91
2. Circadian variation in oxidative stress markers in healthy and type II diabetic men. Kanabrocki, E.L., et. al., (2002) Chronobiology International 19(2):423-39
3. Oxidative DNA damage in human peripheral leukocytes induced by massive aerobic exercise. Tsai, K. et al. (2001). Free Radic. Biol. Med. 31(11):1465-72



## Aconitase-340 Assay

21041

## Quantitative Measurement of Aconitase Enzyme Activity

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21041

There are three biological components most commonly looked at in assessing oxidative insult: lipid peroxidation, protein oxidation and DNA oxidation. While excellent means for assessing lipid peroxidation and DNA oxidation exist, assessment of protein oxidation has been problematic. Carbonylated proteins, once thought to be a very useful marker, have not proven adequate due to sample instability and reproducibility problems.

According to recent literature, decreased aconitase enzyme activity is a sensitive and specific indicator of oxidative damage during aging, Parkinson's and other disease progression. The function of aconitase is to isomerize citrate to isocitrate a key intermediate of the citric acid cycle. Because of its role in cellular energy production, aconitase enzyme function is well positioned as an important marker relative to biological decline.



Two forms of aconitase are known (mitochondrial and cytosolic). The cytosolic form is a bifunctional protein. The holo (4Fe/4S) form has aconitase activity while the fully apo (iron-free) form, known as Iron Regulatory Protein-1 (IRP-1) is devoid of enzymatic activity but plays an important role in regulating the intracellular iron level. In a complete cellular homogenate, both mitochondrial and cytosolic forms will contribute to the total aconitase activity, with the mitochondrial aconitase being predominant in most tissue types.

## Aconitase-340 Method

The OxisResearch® Aconitase-340 assay utilizes the coupled reaction of citrate to isocitrate and isocitrate to  $\alpha$ -ketoglutarate as the basis for quantitating aconitase enzyme activity.

Under the assay conditions, the rate of NADH production is a measure of aconitase activity. To test the performance of the OXIS kit on biological samples, aconitase activity was measured in liver samples from a parenterally fed rat model. Using the Aconitase-340 assay, activity levels of parenterally fed animals were significantly ( $P = 0.0002$ ) reduced compared to controls.

Aconitase Activity in Total Parenteral Rat Model

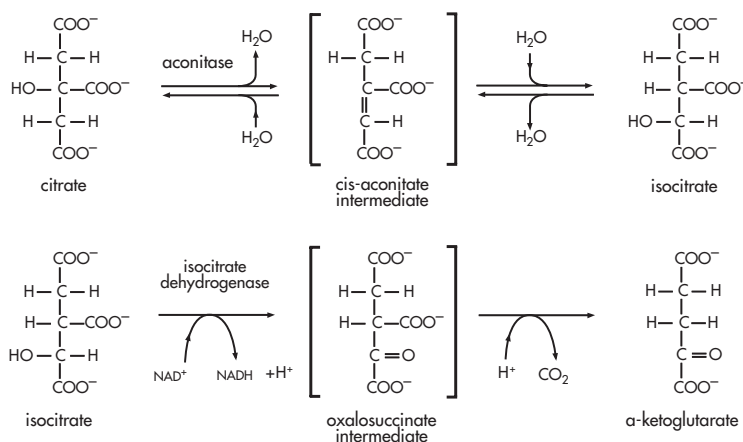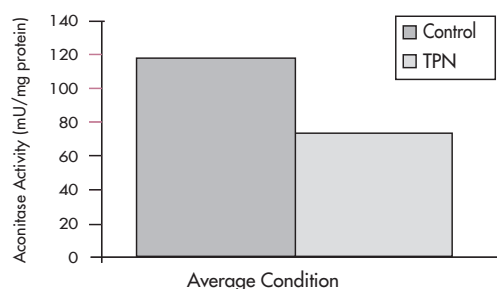

## References

- Gardner, PR, (1997) Superoxide-Driven Aconitase Fe-S Center Cycling, *Bioscience Reports*, 17: 33-42.
- Gardner, PR and Fridovich, I, (1992) Inactivation-Reactivation of Aconitase In *E. Coli*: A Sensitive Measure of Superoxide Radical, *J. Biol. Chem.* 267: 8757-8763.
- Kennedy, MC, et. al., (1997) An EPR Investigation of the Products of the Reaction of Cytosolic and Mitochondrial Aconitases with Nitric Oxide, *J. Biol. Chem.* 272: 20340-20347.
- Yan, LJ, et. al., (1997) Oxidative Damage During Aging Targets Mitochondrial Aconitase, *PNAS*, 94: 11168-11172.
- Rose, IA and O'Connell, EL, (1967) Mechanism Of Aconitase Action, *J. Biol. Chem.* 242: 1870-1879.
- Hentz, MW and Kühn, LC, (1996) Molecular Control of Vertebrate Iron Metabolism: mRNA based Regulatory Circuits Operated by Iron, Nitric Oxide and Oxidative Stress, *PNAS*, 93:8175-8182.

**Aconitase-340 Product Summary**

Intended Use:	Quantitative measurement of aconitase enzyme activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric.		
Kit Contents:	<ul style="list-style-type: none"> <li>• Assay Buffer</li> <li>• Substrate</li> <li>• Enzyme</li> <li>• NAD<sup>+</sup></li> <li>• Standards</li> </ul>		
Storage and Stability:	12 months from date of manufacture when stored as specified.		
Specimen Requirements:	Tissue or cultured cells.		
Assay Precision:	<u>Aconitase in buffer</u>		<u>Biological</u>
n=18	<u>Low</u>	<u>High</u>	<u>Rat Lung</u>
Mean (A 340 /min)	0.0052	0.0254	0.0049
Intra-assay (% CV)	5.11	1.63	3.64
Inter-assay (% CV)	4.82	3.53	8.63
Total Precision (% CV)	6.00	3.71	9.01
Sensitivity:	0.625 mU/ mL in the reaction mixture. 2.5mU/ mL in original sample.		

**Aconitase-340 Selected Research Applications**

The OXIS assay for aconitase enzyme activity is gaining favor among researchers looking for novel ways to quantify the oxidation of proteins. The assay has not been on the market long enough to show many citations but here we highlight one that has come to our attention.

Doxorubicin is an anthracyclin antibiotic sometimes often used in treatment of human neoplasmas such as lymphomas, acute leukemias, stomach, breast, ovarian and cancers of the bone. Its clinical use is limited in that it may also induce cardiotoxicity. Childs, Leeuwenburgh et al. investigated the mechanisms of this cardiotoxicity employing the OXIS Aconitase-340 assay in the process. Although not statistically significant, they showed aconitase activity was decreased in the heart of Doxorubicin treated rats compared to controls (1). They also showed elevated serum levels of free iron. It is interesting to note that aconitase is also an iron storage protein but it is unknown whether its deactivation contributed to the elevated free iron in the serum.

1. Doxorubicin treatment *in-vivo* causes cytochrome c release and cardiomyocyte apoptosis, as well as increased mitochondrial efficiency, superoxide dismutase activity and Bcl-2:Bax ratio. Childs, et al. (2002) Cancer Research 62:4592-4598.

## Bioxytech® HAE-586 Assay

21043

## Quantitative Measurement of 4-Hydroxyalkenals

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21043



Lipid peroxidation, a well established mechanism of cellular injury in both plants and animals, is used as an indicator of oxidative stress in various biological samples. Lipid peroxides, derived from polyunsaturated fatty acids, are unstable and decompose to form a complex series of compounds including malondialdehyde (MDA) and 4-hydroxyalkenals (4-HAE). The most abundant 4-hydroxyalkenal formed by lipid peroxidation is 4-hydroxynonenal (4-HNE). HNE adducts, which are cytotoxic and mutagenic, can alter signal transduction and gene expression. 4-HNE as a biomarker has implications in the areas of neurodegenerative disease, atherosclerosis and nutrition research relative to food freshness and product stability.

The Bioxytech® LPO-586 from OxisResearch® has traditionally been used to measure either the sum of MDA and 4-hydroxyalkenal or MDA only. In response to an increased interest in 4-HNE, OxisResearch® has developed a new assay specific for the HAE component of lipid peroxidation by-products.


## HAE-586 Method

The OxisResearch® HAE-586 assay employs an extraction step that allows quantitative separation and detection of HAE in the presence of excess MDA. After sample extraction, reaction of N-methyl-2-phenylindole (NMPI) with hydroxyalkenals yields a stable carbocyanine dye visible at 586 nm.

The by-products of the reaction also react to form a stable chromophore visible at 505nm. The graph below shows the characteristic spectra obtained from the HAE-586 assay.

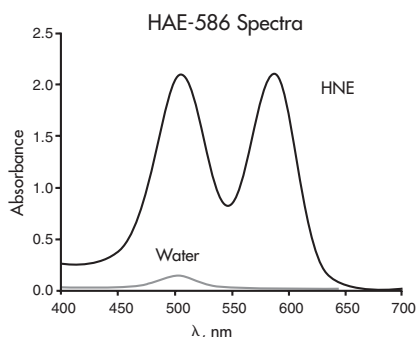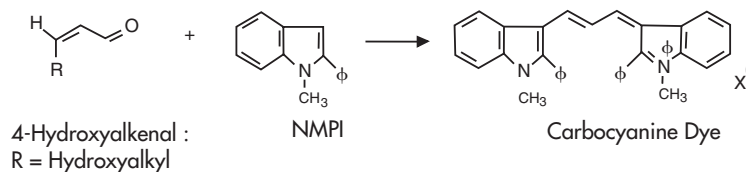

As an example of the assay's usefulness, a study was designed to test the HAE content of various cooking oils and their susceptibility to lipid peroxidation. The data shows the ability of the HAE-586 assay to detect sample HAE content.

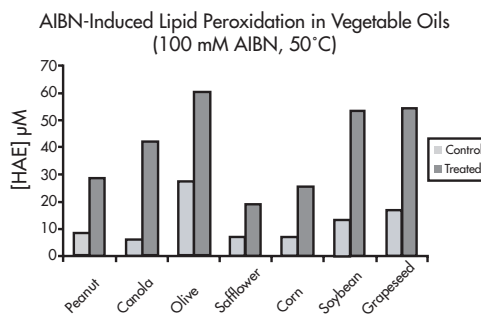

## References

1. Esterbauer, H. et al., (1991) Chemistry and Biochemistry of 4-Hydroxynonenal, Malondialdehyde and Related Aldehydes; Free Rad. Biol. Med. 11, 81-128.
2. Gérard-Monnier et al., (1998) Method of Colorimetric Analysis of Malonic Dialdehyde and 4-Hydroxy-2-enaldehydes as Indexes of Lipid Peroxidation, Kits for Carrying Out Said Method, Substituted Indoles for Use in Said Method and Their Preparation, US Patent No. US5726063.
3. Gérard-Monnier, D., Erdelmeier, I., Régnard, K., Moze-Henry, N., Yadan, J-C., and Chaudierè, J. (1997) Reactions of N-Methyl-2-phenylindole with Malondialdehyde and 4-Hydroxyalkenals. Analytical Applications to a Colorimetric Assay of Lipid Peroxidation, Chemical Research in Toxicology 11, 1176-1183.
4. OXIS Health Products, unpublished data. 5. Liu, J. et al, (1997) Assay of Aldehydes from Lipid Peroxidation: Gas Chromatography-Mass Spectrometry Compared to Thiobarbituric Acid; Analyt. Biochem. 245, 161-166.
6. Erdelmeier, I. et al., (1997) Reactions of N-Methyl-2-phenylindole with Malondialdehyde and 4-Hydroxyalkenals. Mechanistic Aspects of the Colorimetric Assay of Lipid Peroxidation, Chemical Research in Toxicology 11, 1184-1194.

**HAE-586 Product Summary**

Intended Use:	Quantitative measurement of 4-hydroxyalkenals. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric.		
Kit Contents:	<ul style="list-style-type: none"> <li>• N-Methyl-2-phenylindole</li> <li>• Methanesulfonic acid</li> <li>• HNE Standard</li> <li>• Butylated Hydroxytoluene (BHT)</li> <li>• Methanol</li> <li>• Dichloromethane</li> </ul>		
Storage and Stability:	12 months from date of manufacture when stored as specified.		
Specimen Requirements:	Cell Lysate & Tissue		
Assay Precision:			
n=10	<u>Low</u>	<u>Med.</u>	<u>High</u>
Mean (µM)	0.99	1.79	4.30
Intra-assay (%CV)	5.1	2.7	2.9
Total Precision (%CV)	5.9	4.2	3.3
Sensitivity:			
LLD in reaction mixture	0.0743 µM		
LLD A586 in sample	0.0078		

**HAE-586 Selected Research Applications**

Parikh et al studied endogenous antioxidant activities and lipid peroxidation (noting that EPUFAs in the brain are unstable and decompose, preferentially forming HAE's) in the brain of rats treated with various antipsychotic drugs. They determined that chronic Haloperidol treatment had the effect of reducing catalase, SOD and GPx activities (1). The OXIS HAE-586 kit was used to show that hydroxyalkenal (HAE) formation was increased in parallel with the decreased endogenous antioxidant activities in Haloperidol treated animals. These trends were not observed under treatment with atypical antipsychotic drugs such as Clozapine, Risperidone and Olanzapine.

1. Differential effects of antipsychotics on expression of antioxidant enzymes and membrane lipid peroxidation in rat brain. Parikh, V. et al. (2003) J. Psychiatric Research 37:43-51

## Bioxytech® MDA-586 Assay

21044

## Quantitative Measurement of Malondialdehyde (MDA)

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21044

Lipid peroxidation, a well-established mechanism of cellular injury in both plants and animals, is used as an indicator of oxidative stress. Lipid peroxides, derived from polyunsaturated fatty acids, are unstable and decompose to form a complex series of compounds including reactive carbonyl compounds. The most abundant of these is malondialdehyde (MDA), which is commonly used as a biomarker of lipid peroxidation. The MDA-586 from OxisResearch® is designed to measure free or total MDA (i.e. free and protein bound Schiff base conjugates) under assay conditions that minimize interference from other lipid peroxidation products such as 4-hydroxyalkenals. The assay improves upon our popular Bioxytech® LPO-586 assay system.


Improvements over existing method:

- Acid hydrolysis defined for "total MDA" measurement.
- Antioxidant reagent included.
- Enhanced sensitivity via 3<sup>rd</sup> derivative spectroscopy.
- Acid reagent included.

The MDA-586 allows for simple, rapid measurement of MDA with superior specificity and reproducibility compared to other commonly used methods such as the thiobarbituric acid reactive substance assay (TBARS).

## MDA-586 Method

The OxisResearch® MDA-586 method is based on the reaction of a chromogenic reagent, N-methyl-2-phenylindole (NMPI), with MDA at 45°C to form an intensely colored carbocyanine dye with a maximum absorption at 586 nm.

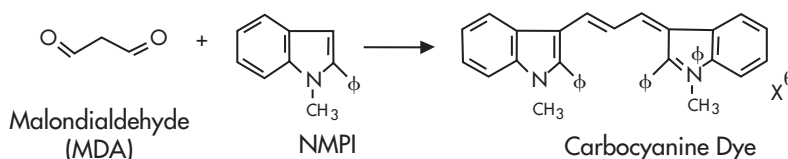


The reaction is carried out in hydrochloric acid with the addition of an antioxidant to further minimize the reaction of 4-hydroxyalkenals. Under these conditions, there is little absorbance at 586 nm from HNE, the most common 4-hydroxyalkenal produced in cells subjected to lipid peroxidation.

The table below shows improved standard deviation (SD) and lower limit of detection (LLD) values for MDA assayed in plasma utilizing the MDA-586 3<sup>rd</sup> derivative spectroscopy procedure.

	Zero Derivative	Third Derivative
SD	0.65 $\mu$ m	0.06 $\mu$ m
LLD	3.0 $\mu$ m	0.2 $\mu$ m *

\*LLD is 0.1  $\mu$ m when four replicates used.

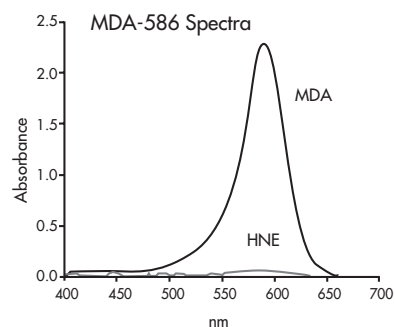

## References

1. Gérard-Monnier, et. al., (1998) Method of Colorimetric Analysis of Malonic Dialdehyde and 4-Hydroxy-2-enaldehydes as Indexes of Lipid Peroxidation, Kits for Carrying Out Said Method, Substitued Indoles for Use in Said Method and Their Preparation, US Patent No. US5726063.
2. Gérard-Monnier, et.al., (1997) Reactions of N-Methyl-2-phenindole with Malondialdehyde and 4-Hydroxyalkenals. Analytical Applications to a Colorimetric Assay of Lipid Peroxidation, Chemical Research in Toxicology 11:10, 1176-1183.
3. Erdelmeier, I., et. al., (1997) Reactions of N-Methyl-2-phenindole with Malondialdehyde and 4-Hydroxyalkenals. Mechanistic Aspects of the Colorimetric Assay of Lipid Peroxidation, Chemical Research in Toxicology 11:10, 1184-1194.
4. OxisResearch®, unpublished data.
5. Carbonneau, M.A. et. al., (1991) Free and Bound Malondialdehyde Measured as Thiobarbituric Acid Adduct by HPLC in Serum and Plasma, Clin. Chem. 37, 1423-1429.
6. Liu, J. et. al., (1997) Assay of Aldehydes from Lipid Peroxidation: Gas Chromatography-Mass Spectrometry Compared to Thiobarbituric Acid, Analyt. Biochem. 245, 161-166. (Additional references available upon request)

**MDA-586 Product Summary**

Intended Use:	Quantitative measurement of malondialdehyde. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• N-methyl-2-phenylindole</li> <li>• Concentrated hydrochloric acid</li> <li>• MDA Standard (tetramethoxypropane, TMOP)</li> <li>• Butylated hydroxytoluene (BHT)</li> <li>• Probucol</li> <li>• Methanol</li> </ul>		
Storage and Stability:	12 months from date of manufacture when stored as specified.		
Specimen Requirements:	Serum, plasma, cell lysate & tissue		
Assay Precision:			
n=10	<u>Low</u>	<u>Med.</u>	<u>High</u>
Mean ( $\mu\text{M}$ )	1.04	1.99	4.79
Intra-assay (%CV)	3.4	1.2	1.6
Total Precision (%CV)	3.7	1.6	1.7
Sensitivity:			
LLD in reaction mixture	0.0801 $\mu\text{M}$		
LLD $A_{586}$ in sample	0.0088		

**MDA-586 Selected Research Applications**

The OXIS MDA-586 assay is still rather new but fast gaining popularity among researchers needing an easy, highly specific method of malondialdehyde (MDA) determination. Here we present a very nice piece of literature by two prominent groups in which the assay was used to evaluate plasma MDA levels.

Dr. Jason D. Morrow and Dr. Lester Packer groups lead by Dr. Block carried out a nationwide study on oxidative stress in human populations. Along with other biomarkers, they looked at plasma MDA levels among different population sets: smokers/non-smokers, age, sex, ethnicity, alcoholic/non-alcoholic etc.

The authors showed that differences in plasma MDA levels were statistically significant in comparisons of smokers vs. nonsmokers, race and the sexes. No significant differences were reported comparing age, weight and alcohol consumption. The authors stated that the OXIS MDA-586 method for measuring MDA in plasma was equally competitive to HPLC method (1).

1. Factors associated with oxidative stress in human populations. Block G et al. (2002). Am. J. Epidemiology 156(3):274-285

$\alpha_1$ AP-410 Assay

21047

Quantitative Measurement of  $\alpha_1$ Antiproteinase Enzyme Activity


For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21047


The most common indicators of oxidative damage are lipid peroxidation, protein oxidation and DNA oxidation. Until recently, the main test for protein oxidation has been carbonylated protein assessment. While sometimes useful, tedious procedures and reproducibility issues have lead researchers to seek alternative methods. For this reason OxisResearch® has sought to provide sensitive assays allowing assessment of specific protein oxidation through measurement of enzyme activities shown to be affected by reactive oxygen species (ROS) and Reactive Nitrogen Species (RNS). The first kit focused on the enzyme aconitase. That assay has the most utility in mitochondria rich tissues. The need and demand for a systemic marker indicative of protein oxidation has prompted the development of the  $\alpha_1$ anti-proteinase ( $\alpha_1$ AP) activity assay. Excess ROS and RNS can inactivate  $\alpha_1$ AP through oxidation of essential methionines.  $\alpha_1$ AP has been shown to be very sensitive to peroxyntirite, hydroxyl radicals and hypochlorous acid. It is well positioned as a key biomarker in the assessment of oxidative stress due to inflammation, disease and other at risk factors in biological systems.

 $\alpha_1$ AP-410 Method

The Bioxytech®  $\alpha_1$ AP-410 Assay is based on the inactivation of elastase by  $\alpha_1$ AP. Since  $\alpha_1$ AP is an irreversible, equimolar inhibitor of elastase, the concentration of active  $\alpha_1$ AP is defined as the molar decrease in elastase activity. The difference between control (elastase alone) and test (elastase plus sample  $\alpha_1$ AP), in the presence of a substrate is used to calculate  $\alpha_1$ AP activity in the samples. Elastase inactivation is measured by monitoring the rate of cleavage of the substrate N-succinyl-(Ala)<sub>3</sub>-*p*-nitroanilide (NSAN). Cleavage of NSAN results in production of a chromagen, *p*-nitroaniline, having a maximal absorbance wavelength at 410nm.

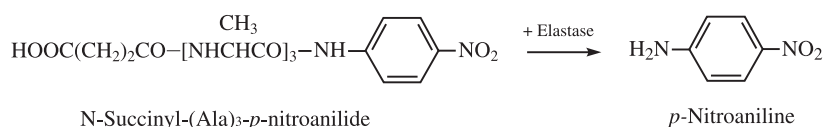

Serum, plasma and other bodily fluids including sputum, synovial fluid and fluid from sites of inflammation are sources of  $\alpha_1$ AP. Some tissues such as lung, liver and pancreas also contain  $\alpha_1$ AP making the assay useful in these tissues as well. The concentration of  $\alpha_1$ AP activity in normal human plasma is 20- 50  $\mu\text{M}$ . The graph to the right shows the assay's utility in monitoring free radical activity in a time-course exercise related study using plasma samples.

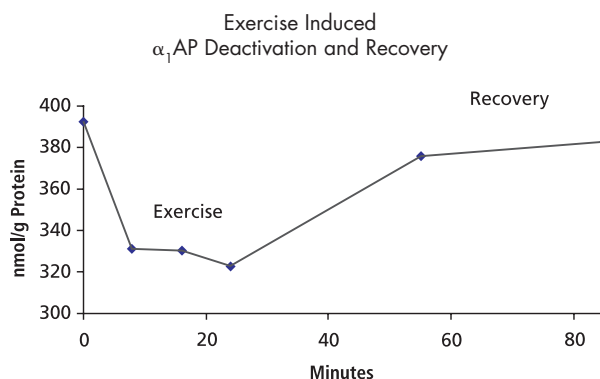

## References

1. Klumpp T and Bieth J (1979) Automated measurement of the elastase-inhibitory capacity of plasma with a centrifugal analyzer *Clin. Chem.* 25, 969-972.
2. Fujita J *et al.* (1990) Evaluation of elastase and antielastase balance in patients with chronic bronchitis and pulmonary emphysema *Am Rev Respir Dis* 142, 57-62.
3. Boudier C and Bieth J (1994) Oxidized mucus proteinase inhibitor: a fairly potent neutrophil elastase inhibitor *Biochem J* 303, 61-68.
4. Whiteman M *et al.* (1999) Modulation of peroxyntirite- and hypochlorous acid-induced inactivation of  $\alpha_1$ -antiproteinase by mercaptoethylguanidinel *British J of Pharmacology* 126, 1646-1652.
5. Browne R *et al.* (1996) Alpha1-antitrypsin deficiency deaths in the United States from 1979-1991: an analysis using multi-cause mortality data *Chest* 110, 78-83.
6. Skosey J and Chow D (1986) *Handbook of Methods for Oxygen Radical Research*, Ed. Robert A. Greenwald, CRC Press Inc. 413-416.
7. Bieth J and Wermuth C (1973) The action of elastase on *p*-nitroanilide substrates *Biochemical and Biophysical Research Communication* 53, 383-390.

**α<sub>1</sub>AP-410 Product Summary**

Intended Use:	Quantitative measurement of α <sub>1</sub> AP enzyme activity. For Research Use Only. Not Intended For Use in Diagnostic Procedures.		
Format:	100 test colorimetric		
Kit Contents:	<ul style="list-style-type: none"> <li>• Assay Buffer</li> <li>• Substrate</li> <li>• Enzyme</li> </ul>		
Storage and Stability:	9 months from date of manufacture when stored as specified.		
Specimen Requirements:	Plasma, serum, tissue or cultured cells.		
Assay Precision:			
n=23	<u>Low</u>	<u>Med</u>	<u>High</u>
Mean (μM)	0.50	0.79	1.29
Intra-assay (% CV)	5.88	5.30	1.08
Inter-assay (% CV)	4.46	4.00	0.23
Total precision (% CV)	6.67	6.01	1.09
Sensitivity:			
LLD in Reaction Mixture	0.025 μM		
LLD in Sample	0.5 μM		




Bioxytech® Urinary 8-EPI-PGF<sub>2α</sub> Assay

21048

*Immunoassay for 8-epi-prostaglandin F<sub>2α</sub> in Urine Samples***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21048

Prostaglandin F<sub>2α</sub> (PGF<sub>2α</sub>) is formed naturally from arachidonic acid via the cyclooxygenase (COX I and II) mediated pathways. It functions as a potent vasoconstrictor *in vivo* under the control of COX I and II. Inhibition of COX activity and subsequent production of PGF<sub>2α</sub> provides the mechanism for the anti-inflammatory properties of common non-steroidal anti-inflammatory drugs (NSAIDs) such as aspirin, ibuprofen and others.

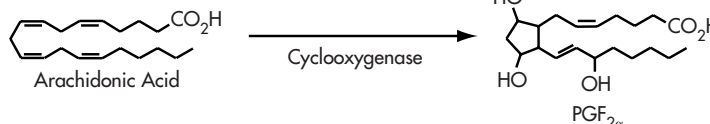


Isoprostanes are prostaglandin-like compounds that are produced upon peroxidation of arachidonic acid. They arise as a result of oxidative insult both *in vivo* and *ex vivo*. Because of this non-enzymatic, oxidative pathway, 8-isoprostane (8-epi-PGF<sub>2α</sub>) has become a popular biomarker relating oxidative stress to various disease pathologies. Similar to the COX derived PGF<sub>2α</sub>, 8-epi-PGF<sub>2α</sub> has been shown to have biological activity as a potent pulmonary and renal vasoconstrictor. In particular urinary isoprostane is well positioned as a non-invasive test for lipid peroxidation associated with oxidative stress. The Bioxytech® Urinary 8-Isoprostane-EIA provides a simple method for determination of 8-isoprostane (8-epi-PGF<sub>2α</sub>) in urine without the need for sample extraction.

**Urinary 8-Isoprostane-EIA Method**

The Urinary 8-Isoprostane-EIA assay is a competitive enzyme-linked immuno-sorbent assay (ELISA) for determining levels of 8-epi-Prostaglandin F<sub>2α</sub> (8-epi) in biological samples. The kit utilizes a pretreatment reagent that does away with the need for sample extraction prior to assay. Briefly, pre-treated samples or standards are added simultaneously with a "free" 8-epi-horseradish peroxidase (HRP) conjugate. Sample or standard 8-epi competes with 8-epi-HRP for binding to the stationary anti-8-epi bound to the plate wells. After washing, Tetramethyl Benzidine (TMB) substrate is added resulting in color development. Color intensity is proportional to the amount of 8-epi-HRP bound and inversely proportional to the amount of 8-epi in the samples or standards. The Bioxytech® 8-isoprostane-EIA has a high degree of correlation with the gold standard Gas Chromatograph/Mass Spectrometry (GC/MS) methods.

The Bioxytech® Urinary 8-isoprostane EIA is strongly recommended for use in urine and does not require solid phase extraction prior to sample assay. For measurement of 8-isoprostane parent compound please refer to our product 21019, 8-isoprostane EIA.

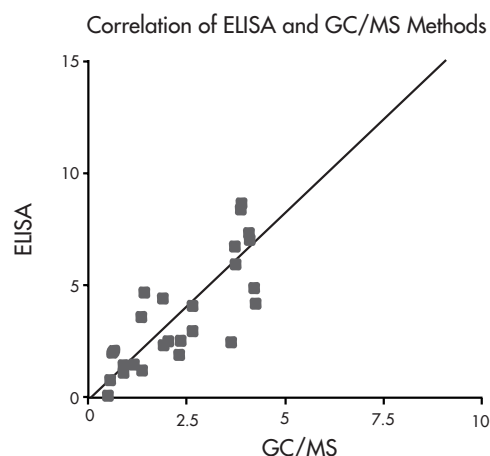**References**

1. Roberts LJ, Morrow JD (1997). The generation and action of isoprostanes. *Biochim Biophys Acta* 1345: 121-135.
2. Morrow JD, Roberts LJ (1997). The isoprostanes: unique bioactive products of lipid peroxidation. *Prog. Lipid Res.* 36(1): 1-21.
3. Kanji VK, Wang C, Salahudeen AK (1999). Vitamin E suppresses cyclosporin A-induced increase in the urinary excretion of arachidonic acid metabolites inducing F<sub>2</sub>-isoprostanes in the rat model. *Transplant Proc.* 31(3): 1724-28.
4. Banerjee M, Kang KH, Morrow JD, Roberts LJ, Newman JH (1992). Effects of a novel prostaglandin 8-epi PGF<sub>2</sub> alpha in rabbit lung in situ. *Am. J. Physiol.* 263: H660-H663.
5. Gniwotta C, Morrow JD, Roberts LJ II, Kuhn H (1997). Prostaglandin F<sub>2</sub>-like compounds, F<sub>2</sub>-isoprostanes, are present in increased amounts in human atherosclerotic lesion. *Arterioscler. Thromb. Vasc. Biol.* 17: 3236-3241.
6. Tangirala RK et al. (2001). Reduction of isoprostanes and regression of advanced atherosclerosis by apolipoprotein E. *J. Biol. Chem.* 276(1): 261-6.
7. Roberts LJ and Morrow JD (2000). Measurement of F<sub>2</sub>-isoprostanes as an index of oxidative stress *in vivo*. *Free Rad. Biol. Medicine.* 28(4): 505-513.

### Urinary 8-Isoprostane-EIA Product Summary

Intended Use:	Quantitative measurement of 8-Isoprostane in urine samples. Analyte also known as 8-epi-Prostaglandin F <sub>2α</sub> and 15-F <sub>2t</sub> -isoprostane. For Research Use Only. Not Intended For Use in Diagnostic Procedures.	
Format:	96 Test Colorimetric (ELISA)	
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 Eell Snti-8-epi Coated Microtiter Plate</li> <li>• 8-epi Standard</li> <li>• Pretreatment Reagent</li> <li>• Wash Buffer</li> <li>• Dilution Buffer</li> <li>• TMB Substrate</li> <li>• 8-epi-HRP Conjugate</li> <li>• Reagent Troughs</li> <li>• Stop Solution</li> </ul>	
Storage and Stability:	12 months from date of manufacture when stored as specified.	
Specimen Requirements:	Urine	
Specificity:	Cross-reactivity at 50% B/B <sub>0</sub>	
	8-epi-PGF <sub>2α</sub>	100%
	9α,11b-Prostaglandin F <sub>2α</sub>	4.1%
	13,14-Dihydro-15-keto PGF <sub>2α</sub>	3.0%
	9b,11α-Prostaglandin F <sub>2α</sub>	<.01%
	Prostaglandin F <sub>2α</sub>	<.01%
	Prostaglandin E <sub>2</sub>	<.01%
	Prostaglandin D <sub>2</sub>	<.01%
	Arachidonic Acid	<.01%
Sensitivity:	0.05ng/mL (precision best on steeper, linear portion of the curve)	

Bioxytech® Urinary F<sub>2</sub>-Isoprostane Metabolite Assay

21049



*Immunoassay for F<sub>2</sub>-Isoprostane Metabolite in Urine***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21049

F<sub>2</sub>-isoprostanes formed as a result of free radical-mediated non-enzymatic peroxidation of membrane-bound arachidonic acid are largely metabolized before excretion. They can be found in esterified tissues, plasma lipids and other body fluids (1). Thus, these isoprostanes can be used to evaluate local or systemic lipid peroxidation *in vivo*. Four classes have been described, each potentially present in 16 isomers. 8-epi-prostaglandin-F<sub>2α</sub> (8-epi-PGF<sub>2α</sub> or sometimes called iP F<sub>2α</sub>-III), a major component of the F<sub>2</sub>-isoprostane family with mitogenic and vasoconstrictor capability, can be noninvasively measured in urine to assess *in vivo* lipid peroxidation (2). In humans, increased F<sub>2</sub>-isoprostane levels have been found in different physiopathological conditions such as atherothrombotic disease, diabetes, hypercholesterolemia, Alzheimer's disease and cigarette smoking.

However, not all F<sub>2</sub>-isoprostanes detected in urine may be of systemic lipid peroxidation origin. Three major urinary and plasma metabolites of 8-epi-PGF<sub>2α</sub> have been reported in humans and rats: 2,3-dinor-5,6-dihydro-8-epi-PGF<sub>2α</sub>, 2,3-dinor-8-epi-PGF<sub>2α</sub> and 2,3,4,5-tetranor-15-keto-13,14-dihydro-8-epi-PGF<sub>2α</sub>. In addition, autoxidation of α-linolenic acid found in plants will also produce 2,3-dinor-5,6-dihydro-8-epi-PGF<sub>2α</sub>, a major urinary metabolite in humans. Thus, the measurement of major metabolites of endogenous 8-epi-PGF<sub>2α</sub> in addition to the parent compound may be useful for both (a) allowing the researcher to obtain a more accurate evaluation of the overall production of the biomarker *in vivo* while adding significance to individual measurements; and (b) provide a compound that can be measured without the risk of artifactual production *ex vivo*. This assay may be used for the quantification of metabolites of 8-epi-PGF<sub>2α</sub> in samples without the need for prior purification or extraction.

**F<sub>2</sub>-Isoprostane Metabolite Method**

The Bioxytech® F<sub>2</sub>-Isoprostane Metabolite Assay is a competitive enzyme-linked immunoassay (ELISA) for determining levels of F<sub>2</sub>-Isoprostane Metabolite. Briefly, the samples are mixed with a pre-treatment reagent. The F<sub>2</sub>-Isoprostane Metabolite in the sample or standard then competes with F<sub>2</sub>-Isoprostane Metabolite conjugated to horseradish peroxidase (HRP Conjugate) for binding to a polyclonal antibody specific for F<sub>2</sub>-Isoprostane Metabolite coated on the microplate. Following substrate addition, the intensity of the color is proportional to the amount of F<sub>2</sub>-Isoprostane Metabolite HRP Conjugate bound and inversely proportional to the amount of unconjugated F<sub>2</sub>-Isoprostane Metabolite in the sample or standard.

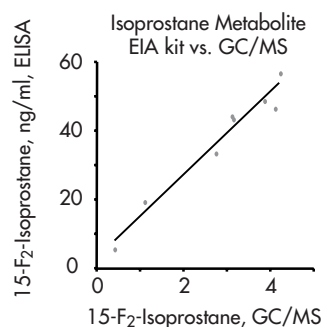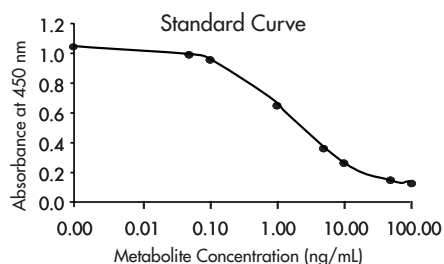

The Bioxytech® Urinary F<sub>2</sub>-Isoprostane Metabolite EIA is strongly recommended for use in urine and does not require solid phase extraction prior to sample assay. For measurement of the 8-isoprostane parent compound, please refer to our products 21019 and 21048.

**References**

- Morrow JD et al. (1990) "A series of prostaglandin F<sub>2</sub>-like compounds are produced *in vivo* in human by non-cyclooxygenase, free radical-catalyzed mechanism." *Proc. Natl. Acad. Sci. USA*. 87: 9383-9387.
- Reilly M et al. (1997) "Modulation of oxidant stress *in vivo* in chronic cigarette smokers." *Circulation* 94: 19-25.
- Reilly M et al. (1997) "Increased formation of distinct F<sub>2</sub>-isoprostanes in hypercholesterolemia." *Circulation* 98: 2822-2828.
- Paratico D et al. (1998) "Increased F<sub>2</sub>-isoprostanes in Alzheimer's disease: evidence for enhanced lipid peroxidation *in vivo*." *FASEB J* 12: 1777-1783.
- Burke A et al. (2000) "Specific analysis in plasma and urine of 2,3-dinor-5,6-dihydro-isoprostane F<sub>2α</sub>-III, a metabolite of isoprostane F<sub>2α</sub>-III and an oxidation product of α-linolenic acid." *J. Biol. Chem.* 275(4): 2499-2504.
- Chiabrando C et al. (1999) "Identification and measurement of endogenous α-oxidation metabolites of 8-epi-prostaglandin-F<sub>2α</sub>." *J. Biol. Chem.* 274(3): 1313-1319.

**F<sub>2</sub>-Isoprostane Metabolite Product Summary**

Intended Use:	Quantitative measurement of F <sub>2</sub> -Isoprostane Metabolite. For Research Use Only. Not Intended For Use in Diagnostic Procedures.	
Format:	96 test colorimetric (ELISA)	
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 well anti-F<sub>2</sub>-Isoprostane Metabolite Coated Microtiter Plate</li> <li>• F<sub>2</sub>-Isoprostane Metabolite Standard</li> <li>• Pretreatment Reagent</li> <li>• Wash Buffer</li> <li>• Dilution Buffer</li> <li>• TMB Substrate</li> <li>• HRP Conjugate</li> </ul>	
Storage and Stability:	12 months from date of manufacture when stored as specified.	
Specimen Requirements:	Urine	
Specificity:	Cross-reactivity at 50% B/B <sub>0</sub>	
	F <sub>2</sub> -Isoprostane Metabolite	100%
	F <sub>2</sub> -Isoprostane (8-epi-Prostaglandin F <sub>2α</sub> )	0.98%
	Prostaglandin F <sub>1α</sub>	0.75%
	11β-Prostaglandin F <sub>2α</sub>	0.29%
	6-Keto Prostaglandin F <sub>1α</sub>	< 0.25%
	Thromboxane B <sub>2</sub>	< 0.25%
	Arachidonic Acid	< 0.25%
Sensitivity:	0.05ng/mL with best precision on steeper, linear portion of the curve above 0.1 ng/mL.	

## Bioxytech® MPO-EIA Assay

21013


## Immunoassay for Human Myeloperoxidase

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21013


Human myeloperoxidase (MPO) is a hemoprotein with a molecular weight of 140 kDa. It is composed of two heavy subunits of 53 kDa and of light subunits of 15 kDa. Each MPO molecule contains two prosthetic porphyrins, which play an essential role in the catalytic cycle.

MPO is stored in primary granules (azurophilic) of neutrophils. It is a major component of the bactericidal armamentarium of neutrophils, due to its capacity to catalyze the production of hypochlorous acid (HOCl) a powerful oxidant. HOCl is derived from the chloride ion ( $\text{Cl}^-$ ) and hydrogen peroxide ( $\text{H}_2\text{O}_2$ ). MPO is a specific marker for polymorphonuclear cells (PMNs). It is released extracellularly via degranulation after PMN activation *in vitro* or *in vivo*. Extracellular MPO can be monitored as an index of PMN activation in a variety of clinical conditions where neutrophils contribute to inflammatory processes and subsequent disease pathology. These include rheumatoid arthritis, adult respiratory distress syndrome (ARDS), septic shock, psoriasis, gout, inflammatory bowel disease (IBD), emphysema and cardiovascular disease.

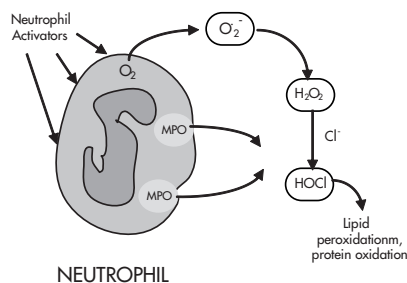

## MPO-EIA Method

The Bioxytech® MPO-EIA assay system is a standard three (3) step "sandwich" ELISA. In step one, a plate bound (e.g. "solid phase") anti-human MPO monoclonal antibody captures sample and/or standard MPO. In step two, the bound MPO antigen is detected using a biotin-labeled polyclonal anti-human MPO. Finally, step three involves binding of an avidin/alkaline phosphatase conjugate to the biotinylated antibody. p-nitrophenyl phosphate (pNPP) is then added causing the enzymatic hydrolysis of pNPP. The formation of yellow p-nitrophenol is monitored at 405 nm and the amount of sample MPO is calculated against the color development of on-board standards.

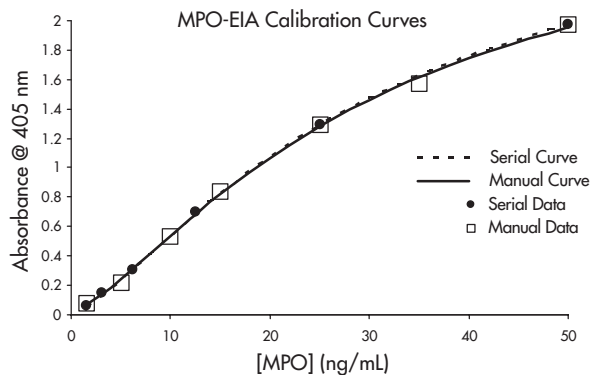

## References

1. Agner, K. (1958) Acta Chem. Scand. 58, 89-94.
2. Olsen, R. L. & Little, c. (1983) Biochem. J. 209, 781-787.
3. Klebanoff, S. J. (1991) in Peroxidases in Chemistry and Biology (J. Everse, K. E. Everse and M. B. Grisham eds.), pp. 1-35, CRC Press, Boston.

**MPO-EIA Product Summary**

Intended Use:	Quantitative measurement of human myeloperoxidase. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	96 Test Colorimetric (ELISA)
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 Anti-Hu-MPO Coated Microtiter Wells (6x16 strips) plus frame</li> <li>• Purified, Lyophilized MPO Standard</li> <li>• Biotin-Coupled Polyclonal Anti-MPO</li> <li>• Avidin-Coupled Alkaline Phosphatase Solution</li> <li>• pNPP Tablets Plus Diluting Buffer</li> <li>• Stop Solution</li> <li>• Diluting and Wash Buffers</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Human samples: serum, plasma, urine, bronchoalveolar lavage (BAL), cerebrospinal fluid (CSF), and cell culture supernatant.
Specificity:	Less than 2% cross reactivity with eosinophil peroxidase.
Sensitivity:	Threshold detection limit of 1.5 ng/mL

**MPO-EIA Selected Research Applications**

The Bioxytech® MPO-EIA has been used in researching various conditions including cardiovascular, fertility, pulmonary, RBC abnormalities, toxicity, inflammatory bowel disease, aging, "critically ill patient status", and transplantation.

Under certain circumstances, an excess of the reactive species produced by MPO (e.g., halides) can overwhelm local defenses and lead to oxidative stress and tissue injury. A process implicated in this pathogenesis is atherosclerosis where MPO is one pathway leading to protein and lipoprotein oxidation. Using OXIS' MPO-EIA kit, Bekesi's (1) showed that MPO could be used as an indicator of age and gender dependent cardiovascular risk. Menopause is positively correlated with decreased particle size following increased oxidation of LDL leading to cardiovascular disease. Hermenegildo (2) studied the effect of MPO level in the plasma of estradiol (E2) treated women. They found that MPO levels might undergo "bimodal function" where low dose E2 treatment promotes decreased levels of MPO (beneficial) and high dose E2 increases MPO levels. Altered functions of neutrophils have been demonstrated in blood from asthmatic patients both before and after challenge with antigens. For example, plant pollen (allergen) sensitized asthmatics were found to exhibit significantly higher levels of MPO released by neutrophils than non-sensitized controls (3). Red blood cell and platelet abnormalities cause essential thrombocythemia (ET) and polycythemia vera (PV). The Falanga group showed that MPO levels in white blood cells of ET and PV patients were significantly increased compared to controls (4). In a toxicity study, Alexis et al. (5) demonstrated that constitutive airway CD-14 (primary receptor for inhaled endotoxin such as lipopolysaccharide (LPS)) expression can predict the magnitude of post LPS inhalation neutrophil response. The authors showed increased CD-14 expression results in increased levels of MPO (5). Increased MPO production has also been reported in patients with inflammatory bowel syndrome. Using a rat model, Pacheco et al. (6) showed that corticoid treatments (budesonide and prednisone) significantly reduced the level of MPO, compared to the control subjects.

1. Plasma concentration of myeloperoxidase enzyme in pre- and post-climacterical people: related superoxide anion generation. Bekesi G et al. (2001). *Experimental Gerontology* 37: 137-148

2. Transdermal estradiol reduces plasma myeloperoxidase levels without affecting the LDL resistance to oxidation or the LDL particle size. Hermenegildo C et al. (2002). *Menopause*. 9 (2): 107-111.

3. IgE-dependent release of MPO by neutrophils from allergic patients Monteseirin, J et al., (2001). *Clin. Exp. Allergy* 31: 889-92.

4. Polymorphonuclear leukocyte activation and hemostasis in patients with essential thrombocythemia and polycythemia vera. Falanga A et al. (2000). *Blood* 96(13): 4261-6.

5. CD-14-dependent airway neutrophil response to inhaled LPS: role of atopy. Alexis N et al. (2001). *J. Allergy Clin Immunol.* 107(1): 31-5.

6. Corticosteroid pretreatment prevents small intestinal mucosal lesion induced by acetic acid-perfusion model in rats. Pacheco I et al. (2000). *Diseases and Sciences* 45: 2337-46

## Bioxytech® Lactof-EIA Assay


21015

*Immunoassay for Human Lactoferrin***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 21015

Human lactoferrin (LTF) is an 80 kDa glycoprotein which was first isolated from human milk. LTF is found in most of the body fluids and secretions, in the nose, genital tract and tear (Autiero, 1991). LTF in blood is secreted from the neutrophils and its plasma concentration is positively related to the total pool of neutrophils and to the rate of neutrophil turnover (Birgens, 1985). Because of its ability to strongly bind iron, LTF is considered to be bactericidal (Baynes, 1986). In a number of cases of inflammation, lactoferrin is released from secondary granules of neutrophilic leukocytes into the extracellular medium (Lash, 1983; Kähler, 1988). Its extracellular concentration can therefore be used as an index of neutrophil activation, especially in blood samples containing anti-myeloperoxidase antibodies.

**Lactof-EIA Method**

The Lactof-EIA method is an enzyme-linked immunosorbent assay (ELISA). Samples are incubated in the wells of a sectionable microplate which have been coated with a monoclonal antibody to LTF. The LTF-MAb complex is detected by a biotinylated monoclonal antibody to LTF. The final step of the assay is an amplification based on a biotin-avidin coupling in which avidin has been covalently linked to horseradish peroxidase. The amount of LTF is measured enzymatically upon addition of orthophenylenediamine (OPD) at 420 nm.

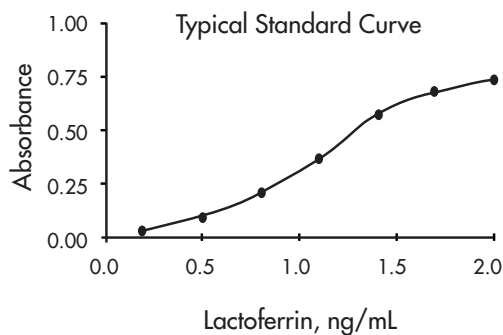**References**

1. Autiero M, Sansone G, Abrescia P. (1991) "Relative ratios of lactoferrin, albumin, and acid phosphatase seminal levels as sperm quality markers in fertile and infertile men" *J. Androl.* 12: 191-200.
2. Birgens HS (1985) "Lactoferrin in plasma measured by an ELISA technique: Evidence that plasma lactoferrin is an indicator of neutrophil turnover and bone marrow activity in acute leukemia" *Scand. J. Haematol.* 34: 326-331.
3. Baynes RD, Bezwoda WR, Khan Q, Mansoor N (1986) "Relationship of plasma lactoferrin content to neutrophil regeneration and bone marrow infusion" *Scand. J. Haematol.* 36: 79-84.
4. Lash JA, Coates TD, Lafuze J, Baehner RL, Boxer L (1983) "Plasma lactoferrin reflects granulocyte activation *in vivo*" *Blood.* 61: 885-888.
5. Kähler S, Christophers E., Schröder JM (1988) "Plasma lactoferrin reflects neutrophil activation in psoriasis" *Brit. J. Derm.* 119: 289-293.

**Lacto<sup>f</sup>EIA Product Summary**

Intended Use:	Quantitative measurement of human lactoferrin. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	96 test colorimetric (ELISA)
Kit Contents:	<ul style="list-style-type: none"> <li>• 96 Anti-LTF Coated Microtiter Wells (6x16 strips) Plus Frame</li> <li>• Purified, Lyophilized LTF Standard</li> <li>• Biotin-Coupled Monoclonal Anti-LTF</li> <li>• Avidin-Coupled Alkaline Horseradish Peroxidase Solution</li> <li>• OPD Tablets Plus Diluting Buffer</li> <li>• Stop Solution</li> <li>• Diluting and Washing Buffers</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Human samples: serum, EDTA plasma, urine, bronchoalveolar lavage (BAL), cerebrospinal fluid (CSF), and cell culture supernatant.
Sensitivity:	Threshold detection limit of 1.0 ng/mL

**Lacto<sup>f</sup>EIA Selected Research Applications**

The OXIS LTF-EIA assay has been used to investigate subjects as diverse as pregnancy, exercise and immunology including AIDS among others.

Pancora et al. studied the concentrations of lactoferrin with gestational age, infection, labor and rupture. Human amniotic fluid and plasma were evaluated for lactoferrin content using the OXIS LTF-EIA kit (1).

Santos-Silva et al. studied oxidative stress in highly competitive swimmers (20 hrs/week) versus moderately exercising (1-2 hr/week) adolescents (12 to 16 yrs) in terms of leukocyte activation and lipid profile. The authors reported elevated lactoferrin, elastase and GM-CSF in competitive swimmers, though only elastase presented a statistically significant difference (2). Note: OXIS now has an  $\alpha_1$ -antiproteinase kit (colorimetric assay).  $\alpha_1$ AP-410 can be used to indirectly evaluate the function of elastase activity. When  $\alpha_1$ -antiproteinase is inactivated (as a result of oxidative stress-protein oxidation) elastase will be higher.

Da Silva et al. studied the effect of the antibody CD11a mAb (ODULIMONAB) in preventing renal ischemia-reperfusion injury during kidney transplantation. A primate model was used wherein histological studies in kidney and creatinine clearance did not provide good indicators for glomerular function/neutrophil infiltration. Lactoferrin, however, did provide a good marker for neutrophil infiltration/retention. It was significantly higher in untreated animals than the ODULIMONAB treated ones (3).

In other studies, Dapsanse et al. showed that plasma levels of LTF in HIV 1-infected ("AIDS") patients was decreased in correlation with the progression of the disease (4). This profile was not found in HIV-2 infected patients (HIV-2 progression is far slower than HIV-1). In parallel with the drop in LTF, a marked increase in circulating nitrogen derivatives (total nitrites) was observed in HIV-1 infected patients whereas a low level was found in normal donors and in HIV-2 infected patients. Thus, they suggest that impairment of circulating LTF could be an important element in HIV- induced oxidative imbalance.

1. Lactoferrin in intrauterine infection, human parturition, and rupture of fetal membrane. Pacora, P et al. (2000). *Am. J. Obstetrics and Gynecology* 283: 904-10

2. Leukocyte activation, erythrocyte damage, lipid profile and oxidative stress imposed by high competition physical exercise in adolescents. Santos-Silva A et al. (2001). *Clinica Chimica Acta* 306:119-126

3. A primate model of renal ischemia-reperfusion injury for pre-clinical evaluation of the anti-leukocyte function associated antigen 1 monoclonal antibody ODULIMONAB. Da Silva, M et al. (2001). *J. Urology*. 168:1915-1919

4. Differential pattern in circulating nitrogen derivatives, lactoferrin, and anti-lactoferrin antibodies in HIV Type 1 and HIV Type 2 infection. Dapsanse V et al. (2001) *AIDS Research and Retroviruses* 17(11): 1041-45



## Bioxytech® Nitrotyrosine-EIA Assay

21055

## Quantitative Measurement of Nitrotyrosine


For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 21055

Nitrotyrosine is a relatively stable product formed from various reaction pathways. Perhaps most notable is the reaction of peroxynitrite (formed from superoxide and nitric oxide radicals) with protein bound tyrosine.

As a strong oxidant and nitrating agent, peroxynitrite mediates tyrosine nitration reactions on proteins resulting in inactivation of certain housekeeping enzymes (e.g.  $\alpha_1$ -antiproteinase) as well as endogenous antioxidant enzymes such as catalase and SOD.


This form of protein nitration has been shown to be involved in the pathology of several human conditions such as chronic myocardial dysfunction, respiratory distress syndrome, inflammatory bowel disease, lung injury, asthma, atherosclerosis, rheumatoid arthritis, chronic renal failure, ALS, septic shock and others.

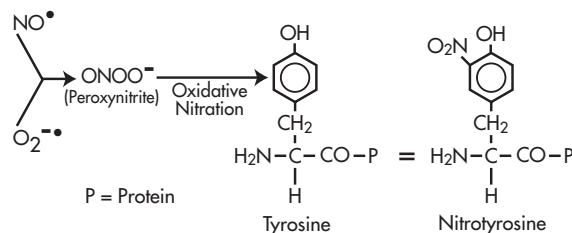

## Nitrotyrosine EIA Method

The Nitrotyrosine-EIA is a "sandwich" ELISA. Antigen captured by a solid phase monoclonal antibody is detected with a biotin-labeled goat polyclonal anti-nitrotyrosine. A streptavidin peroxidase conjugate then binds to the biotinylated antibody. The tetramethylbenzidine (TMB) substrate is added and the yellow product is monitored at 450 nm.

Demonstration of nitrotyrosine in biological samples is useful in determining oxidative stress level and specifically infers the presence of peroxynitrite or related nitrogen-centered oxidants. The new Nitrotyrosine ELISA kit from OxisResearch® provides a simple and reproducible method for NT detection in tissues and plasma.

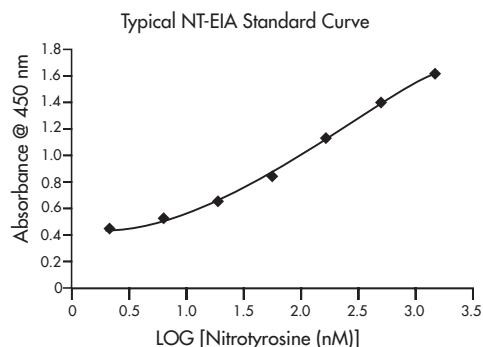

## References

1. van der Vliet A, et al. (2000) "Nitric oxide: a proinflammatory mediator in lung disease?" *Rep. Res.* 1:67-72.
  2. Wu W, et al. (1999) "Eosinophil peroxidase nitrates protein tyrosyl residues" *J. Biol. Chem.* 274(36):25933-44.
  3. van der Vliet A, et al. (1999) "Reactive nitrogen species and tyrosine nitration in the respiratory tract: Epiphenomena or a pathobiological mechanism of disease" *Am. J. Respir. Crit. Care Med.* 160: 1-9.
  4. Whiteman M., et al. (1996) "Protection against peroxynitrite-dependent tyrosine nitration and  $\alpha_1$ -antiproteinase inactivation by oxidized and reduced lipoic acid" *FEBS Lett.* 379: 74-6.
  5. Keng T., et al. (2000) "Peroxynitrite formation and decreased catalase activity in autoimmune MRL-lpr/lpr mice" *Mol. Med.* 6(9): 779-92.
  6. MacMillan-Crow LA, et al. (1996) "Nitration and inactivation of manganese superoxide dismutase in chronic rejection of human renal allografts" *Proc. Natl. Acad. Sci. USA* 93: 11853-8.
  7. Monezen SL, Augustus O (2001) "EPR detection of glutathionyl and protein-tyrosyl radicals during the interaction of peroxynitrite with macrophages (J774)" *J. Biol. Chem.* 276(43): 37679-84.
  8. Kooy NW, et al. (1997) "Extensive tyrosine nitration in human myocardial inflammation: evidence for the presence of peroxynitrite" *Crit. Care Med.* 25(5): 812-9.
  9. Beal MF, et al. (1997) "Increased 3-nitrotyrosine in both sporadic and familial amyotrophic sclerosis" *Ann. Neurol.* 42: 646-54.
  10. Sittipunt C, et al. (2001) "Nitric oxide and nitrotyrosine in the lungs of patients with acute respiratory distress syndrome" *Am. J. Res. Crit. Care Med.* 163: 503-510.
  11. Singer II, et al. (1996) "Expression of inducible nitric oxide and nitrotyrosine in colonic epithelium in inflammatory bowel disease" *Gastroenterology* 111: 8771-85.
  12. Chang L-Y, Crapo JM, (2002) "Inhibition of airway inflammation and hyper-reactivity by an antioxidant mimetic" *Free Radic. Biol. Med.* 33(3): 379-85.
- (Additional references available upon request)

**Nitrotyrosine EIA Product Summary**

Intended Use:	Quantitative measurement of nitrotyrosine in biological samples. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	2 X 96 well plates
Kit Contents:	<ul style="list-style-type: none"> <li>• Sample Diluting Buffer</li> <li>• Nitrotyrosine Standard</li> <li>• Nitrotyrosine Antibody</li> <li>• Streptavidin Peroxidase</li> <li>• TMB Substrate</li> <li>• Substrate Buffer</li> <li>• Substrate Diluting Buffer</li> <li>• Wash Buffer</li> <li>• Stop Solution</li> <li>• Pre-Coated Plates w/ Frame</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Tissue, plasma
Assay Precision:	
Intra-assay (%CV)	2.32
Inter-assay (%CV)	11.17
Sensitivity:	LLD = 2 nM

## Bioxytech® Nitric Oxide Assay (enzymatic)

22110

*Quantitative Measurement of Total Nitrite in Aqueous Samples***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 22110

Nitric Oxide (NO; also known as endothelial derived relaxation factor or EDRF) is produced in trace quantities in neurons, endothelial cells, platelets and neutrophils in response to homeostatic stimuli. Under these conditions, NO acts in a paracrine fashion to transduce cellular signals. NO is also produced by other cell types (macrophages, fibroblasts, hepatocytes) in micromolar concentrations as a response to inflammatory or mitogenic stimuli. In this case, the biological role is defense against pathogens through oxidative toxicity. NO reacts readily with other radicals, metal ions and protein thiols resulting in a short biological half-life. High NO levels can lead to the formation of peroxynitrite (ONOO), destruction of iron-sulfur clusters, thiol nitrosation and nitration of protein tyrosine residues.

The Bioxytech® NO Enzymatic assay provides reagents for use in the determination of nitrite as an indicator of nitric oxide production in biological samples. Nitric oxide is rapidly converted to nitrite and nitrate in typical oxygenated aqueous solutions. Because an excellent colorimetric reagent (the Griess reagent) exists for the determination of nitrite, it is common practice to use enzymatic or chemical reduction to convert all nitrate to nitrite in a sample and measure total nitrite as an indicator of nitric oxide production. This assay provides for enzymatic reduction of nitrate by nitrate reductase, followed by spectrophotometric analysis of total nitrite using Griess reagent. In addition to providing all necessary components in a microtiter format, this kit employs affinity purified nitrate reductase and NADH, thereby circumventing some of the potential problems reported for nitrite measurement using NADPH dependent nitrate reductases.

**NO-Enzymatic Method**

In aqueous solution, nitric oxide rapidly degrades to nitrate and nitrite. Spectrophotometric quantitation of nitrite using Griess Reagent is straightforward and sensitive, but does not measure nitrate, causing a possible underestimation of nitric oxide. This kit employs the NADH-dependent enzyme nitrate reductase for enzymatic reduction of nitrate to nitrite (Figure 1) prior to quantitation of nitrite using Griess reagent. In acid solution, nitrite is converted to nitrous acid ( $\text{HNO}_2$ ) which diazotizes sulfanilamide. This sulfanilamide-diazonium salt is then reacted with N-(1-Naphthyl)-ethylenediamine (NED) to produce a chromophore which is measured at 540 nm.

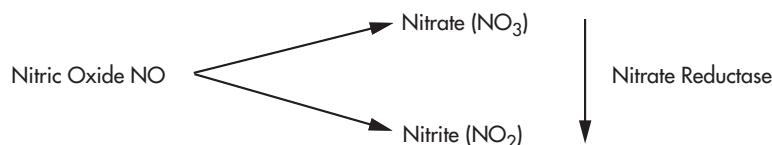


Figure 1 - Enzymatic Reduction of Nitrate to Nitrite

This kit can be used to accurately measure as little as 1  $\mu\text{M}$  of nitrite (final concentration in the assay). Very little sample is required (5  $\mu\text{L}$  to 85  $\mu\text{L}$  for most samples) so it is the assay we recommend for testing of samples such as rodent tail bleeds etc. For larger samples we recommend our NO-540 (22116) cadmium reductant assay.

**References**

- Schmit, H.H., et. al. *Biochemica* 2:22-23 (1995).

**NO-Enzymatic Product Summary**

Intended Use:	Quantitative measurement of total nitrite in aqueous samples. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	96 Test Microplate
Kit Contents:	<ul style="list-style-type: none"> <li>• Nitrate Reductase</li> <li>• NADH</li> <li>• Color Reagent #1 (Sulfanilamide)</li> <li>• Color Reagent #2 (N-(1-Naphthyl) ethylenediamine dihydrochloride)</li> <li>• Nitrate Standard</li> <li>• Microtitre Plate</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Tissue homogenates, plasma, cell lysate and culture supernatant .
Sensitivity:	1.0 µM

**NO-Enzymatic Selected Research Applications**

The authors studied circadian rhythms in urinary MDA, urinary 8-Isoprostane, urinary 8-OHdG and serum's total nitrite and uric acid in diabetic (4) and "healthy" men (7), using NO-22110, 8-OHdG-EIA, and 8-Isoprostane-ELISA.

1. Circadian variation in oxidative stress markers in healthy and type II diabetic men. Kanabrocki, E.L., et. al., (2002) Chronobiology International 19(2):423-39

## Bioxytech® Nitric Oxide Assay (non-enzymatic)

22111

*Quantitative Measurement of Total Nitrite in Aqueous Samples***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 22111

Nitric Oxide (NO; also known as endothelial derived relaxation factor or EDRF) is produced in trace quantities in neurons, endothelial cells, platelets and neutrophils in response to homeostatic stimuli. Under these conditions, NO acts in a paracrine fashion to transduce cellular signals. NO is also produced by other cell types (macrophages, fibroblasts, hepatocytes) in micromolar concentrations as a response to inflammatory or mitogenic stimuli. In this case, the biological role is as a defense against pathogens through oxidative toxicity. NO reacts readily with other radicals, metal ions and protein thiols resulting in a short biological half-life. High NO levels can lead to the formation of peroxynitrite (ONOO), destruction of iron-sulfur clusters, thiol nitrosation and nitration of protein tyrosine residues.

The Bioxytech® NO-Non-Enzymatic assay provides reagents for use in the determination of nitrite as an indicator of nitric oxide production in biological samples. Nitric oxide is rapidly converted to nitrite and nitrate in typical oxygenated aqueous solutions. Because an excellent colorimetric reagent (the Griess reagent) exists for the determination of nitrite, it is common practice to use enzymatic or chemical reduction to convert all nitrate to nitrite in a sample and measure total nitrite as an indicator of nitric oxide production. This assay provides for chemical reduction of nitrate by granulated cadmium, followed by spectrophotometric analysis of total nitrite using Griess reagent. In addition to providing all necessary components in a microtiter format, this kit provides for elimination of possible interference by proteins in concentrated biological samples by ZnSO<sub>4</sub> precipitation. The granulated cadmium beads provide for a quantitative and inexpensive assay.

**NO-Non-Enzymatic Method**

In aqueous solution, nitric oxide rapidly degrades to nitrate and nitrite. Spectrophotometric determination of nitrite using Griess reagent is straight forward and sensitive, but does not measure nitrate, causing a possible underestimation of nitric oxide. This kit employs granular cadmium metal for chemical reduction of nitrate to nitrite (Figure 1) prior to quantitation of nitrite using Griess reagent. In acid solution, nitrite is converted to nitrous acid (HNO<sub>2</sub>) which diazotizes sulfanilamide. This sulfanilamide-diazonium salt is then reacted with N-(1-Naphthyl)-ethylenediamine (NED) to produce a chromophore which is measured at 540 nm.

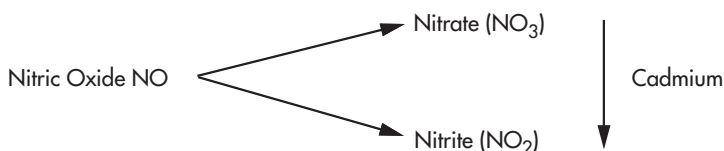


Figure 1 - Chemical Reduction of Nitrate to Nitrite

This kit can be used to accurately measure as little as 0.25 µM of nitrite (final concentration in the assay). Very little sample is required (10 µL to 100 µL for most samples).

**References**

- Schmidt, H.H., et. al. *Biochemica* 2:22-23 (1995).

**NO-Non-Enzymatic Product Summary**

Intended Use:	Quantitative measurement of total nitrite in aqueous samples. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	96 test microplate
Kit Contents:	<ul style="list-style-type: none"> <li>• Granulated Cadmium (25g Cd beads)</li> <li>• ZnSO<sub>4</sub> Solution (30% wt/vol)</li> <li>• Microcentrifuge Tubes (50 X 1 mL)</li> <li>• Color Reagent #1 (Sulfanilamide)</li> <li>• Color Reagent #2 (N-(1-Naphthyl) ethylenediamine dihydrochloride)</li> <li>• Nitrite standard</li> <li>• Microtitre Plate</li> <li>• Cadmium Bead Washing Solution</li> </ul>
Storage and Stability:	12 months from date of manufacture when stored as specified.
Specimen Requirements:	Tissue homogenates, plasma, cell lysate and culture supernatant.
Sensitivity:	0.25 µM

**NO-Non-Enzymatic Selected Research Applications**

Hyperglycemia is linked to vascular and glomerular dysfunction by many mechanisms, including a glucose-dependent abnormality in nitric oxide production and action. The authors investigated NO biosynthesis and action associated with glomerular infiltration and persistent micro albuminuria in type 1 diabetes patients.

1. Increased circulating nitric oxide in young patients with type 1 diabetes and persistent micro albuminuria: relation to glomerular hyperfiltration. Chiarelli F et al., (2000). *Diabetes* 49: 1258-63.

## Bioxytech® Nitric Oxide Synthase Assay (radioactive arginine) 22112

## Quantitative Assay for the Determination of NOS Activity

Catalog Number

22112

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Three distinct isoforms of NOS have been identified: inducible (iNOS), endothelial (eNOS) and neuronal (nNOS or bNOS). Nitric Oxide Synthase is a remarkably complex enzyme that acts on molecular oxygen and arginine in neurons, endothelial cells, platelets, neutrophils and other cells to produce nitric oxide (NO). Nitric oxide is a unique second messenger molecule that readily diffuses through cell membranes to exert a variety of biological actions in mammalian cells. Excess generation of NO can lead to the formation of peroxynitrite, destruction of iron-sulfur clusters, thiol nitrosylation and nitration of protein tyrosine residues.

Measuring NOS activity by monitoring the conversion of arginine to citrulline is a widely accepted assay for NOS activity in both crude and purified enzyme preparations. Advantages of the NOS assay kit include the use of radioactive substrates ( $^3\text{H}$ -arginine or  $^{14}\text{C}$ -arginine), and an easy method for separating neutrally charged citrulline from positively charged arginine.


The citrulline assay has been used to quantify levels of NOS activity in tissue homogenates from numerous sources including blood vessels, immune cells, visceral organs, nervous tissue and cultured cells. Nitric oxide synthase is relatively unstable; therefore, tissues should be harvested quickly after animal euthanasia. If enzyme assays are to be performed at a later time, it is best to freeze intact tissues or harvested cultured cells prior to homogenization. Wrap the tissues in aluminum foil, flash freeze in liquid nitrogen and then store at  $-70^\circ\text{C}$ .

The subcellular distribution of NOS is tightly regulated in tissues. Endothelial NOS is largely membrane associated as a result of N-terminal myristoylation. bNOS is found primarily in the cytoplasmic fractions in adult rat brain, yet in skeletal muscle, it is predominately associated with membrane fractions. Nitric oxide synthase in soluble and membrane-associated fractions can be separated by centrifuging the homogenized tissues at  $100,000 \times g$  for 60 minutes. The supernatant contains soluble NOS, while the pellet, which can be re-suspended in homogenization buffer, will contain membrane-associated NOS.

## NOS Arginine Method

This NOS assay kit is based on the biochemical conversion of L-arginine to L-citrulline by NOS. This reaction involves a five-electron oxidation of a guanidinonitrogen of L-arginine to nitric oxide (NO), together with the stoichiometric production of L-citrulline. The reaction consumes 1.5 equivalents of reduced nicotinamide adenine dinucleotide phosphate (NADPH) and also requires molecular oxygen, calcium, calmodulin and tetrahydrobiopterin.

For routine assays, radioactive arginine is added to tissue or cellular extracts. After incubation, the reactions are stopped with a buffer containing EDTA, which chelates the calcium required by NOS and, consequently, inactivates the NOS. Equilibrated resin, which binds to arginine, is added to the samples and they are centrifuged in spin cups. The citrulline, being ionically neutral at pH 5.5, passes through the spin cup column. The NOS activity is then reported by quantitating the radioactivity in the eluate. Nitric oxide synthase activity in the citrulline assay is defined as counts per minute (cpm) in an incubated test sample as compared to an appropriate blank.

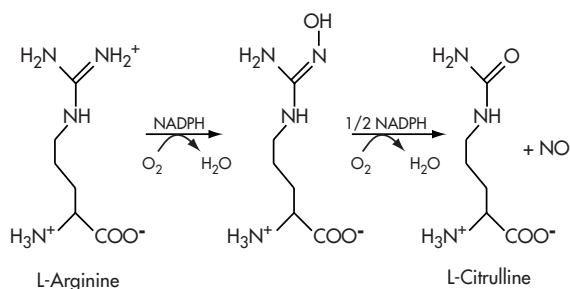

NOS catalyzes a 5-electron oxidation of an amidine nitrogen of L-arginine to generate NO and L-citrulline. L-Hydroxyarginine is formed as an intermediate that is tightly bound to the enzyme. Both steps in the reaction are calcium and calmodulin dependent.

## References

1. Bredt, D.S., and Snyder, S.H. 1994. *Annu. Rev. Biochem.* 63, 175.
2. Marletta, M.A. 1993. *J. Biol. Chem.* 268, 12231.
3. Moncada, S., and Higgs, A. 1994. *N. Engl. J. Med.* 329, 2002.
4. Nathan, C., and Xie, Q.W. 1994. *J. Biol. Chem.* 269, 13725.
5. Schmidt, H.H., et al. 1993. *Biochim. Biophys. Acta* 1178, 153.
6. Stuehr, D.J., et al. 1991. *Proc. Natl. Acad. Sci. U.S.A* 88, 7773.
7. Schmidt, H.H., et al. 1991. *Proc. Natl. Acad. Sci. U.S.A* 88, 365.
8. Busconi, L., and Michel, T. 1993. *J. Biol. Chem.* 268, 8410.
9. Kobzik, L., et al. 1994. *Nature* 372, 546.

**NOS Arginine Product Summary**

Intended Use:	Determination of NOS activity by monitoring the conversion of tritiated L-arginine to L-Citrulline. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	50 tests
Kit Contents:	Part 1, room temperature storage - <ul style="list-style-type: none"> <li>• Homogenization Buffer</li> <li>• Stop Buffer (stops cNOS reactions)</li> <li>• Equilibrated Arginine-Binding Resin</li> <li>• Calcium Chloride (CaCl<sub>2</sub>)</li> <li>• Elution Buffer</li> <li>• Spin Cups and Cup Holders</li> </ul> Part 2, -70°C storage - <ul style="list-style-type: none"> <li>• Cerebellum Extract</li> <li>• Calmodulin</li> <li>• Reaction Buffer</li> <li>• NADPH Tetrasodium Salt</li> <li>• N<sup>G</sup>-Nitro-L-arginine Methyl Ester, HCl, (L-NAME, HCl),</li> </ul>
Storage and Stability:	2 years from date of manufacture when stored as specified.
Specimen Requirements:	Tissue homogenates and cell lysates.

**NOS Arginine Selected Research Applications**

Role of nitric oxide relative to gastric mucosal injury is examined. Nitric oxide appears to play a protective role in the gastric mucosa as an endogenous vasodilator and its hyperemic response to damaging agents. The authors used NO modulators (L-arginine and NG-nitro-L-arginine methyl ester or L-NAME) in rats and measured the NOS activity of gastric tissue using the OXIS NOS 22112 kit (1).

Nishima et al evaluated the effect of a new NOS inhibitor in a hamster model of diaphragmatic dysfunction. Hamster diaphragmatic muscle was tested for nitric oxide using the OXIS NO assay (22112).

Propofol, a peroxynitrite scavenger and antioxidant, is often used in critically ill patients as a sedative or anesthetic. A potential negative effect of the antioxidant activity of this product includes increased severity of sepsis. Mikawa et al. investigate whether propofol may suppress the lipid peroxidation and NO species formation. They evaluated the effect of a new iNOS inhibitor on a rat model of pancreatitis wherein diaphragmatic muscle homogenate was tested for nitric oxide using the OXIS NO assay (22112).

1. Role of nitric oxide and mucus in ischemia/reperfusion-induced gastric mucosal injury in rats. Kim, H et al. (2001). *Pharmacology* 62: 200-207.

2. ONO1714, a new inducible nitric oxide synthase inhibitor, attenuates sepsis-induced diaphragmatic dysfunction in hamsters. Nishima K, et al. (2001). *Anesthesia and analgesia* 92: 959-66.

3. ONO-1714, a new inducible nitric oxide synthase inhibitor, attenuated diaphragmatic dysfunction associated with cerulein-induced pancreatitis in rats. Mikawa K et al. (2001). *Crit. Care Med.* 29(6) 1215-21.



## Bioxytech® Nitric Oxide Synthase Assay (colorimetric)

22113

## Colorimetric Assay for the Measurement of NOS Activity

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 22113

Three distinct isoforms of NOS have been identified: inducible (iNOS), endothelial (eNOS) and neuronal (nNOS or bNOS). Nitric Oxide Synthase is a remarkably complex enzyme that acts on molecular oxygen and arginine in neurons, endothelial cells, platelets, neutrophils and other cells to produce nitric oxide (NO). Nitric oxide is a unique second messenger molecule that readily diffuses through cell membranes to exert a variety of biological actions in mammalian cells. Excess generation of NO can lead to the formation of peroxynitrite, destruction of iron-sulfur clusters, thiol nitrosylation and nitration of protein tyrosine residues. The final products of NO *in vivo* are nitrate ( $\text{NO}_3^-$ ) and nitrite ( $\text{NO}_2^-$ ).


The Bioxytech® NOS Colorimetric Assay Kit is intended for the quantitative determination of total nitrite as an indicator of nitric oxide synthase (NOS) activity in biological samples. It may be used for the assay of total nitrite in urine, plasma, serum and tissue culture medium.

The subcellular distribution of NOS is tightly regulated in tissues. Endothelial NOS is largely membrane associated as a result of N-terminal myristoylation. nNOS is found primarily in the cytoplasmic fractions in adult rat brain, yet in skeletal muscle, it is predominately associated with membrane fractions. Nitric oxide synthase in soluble and membrane-associated fractions can be separated by centrifuging the homogenized tissues at 100,000 x g for 60 minutes. The supernatant contains soluble NOS, while the pellet, which is resuspended in homogenization buffer, contains membrane-associated NOS.

## NOS-Colorimetric Method

NO exists primarily as  $\text{NO}_2$  and  $\text{NO}_3$  in biological systems. Its measurement can be a useful tool in estimating NOS activity. The principle of this assay is based on the measurement of  $\text{NO}_2^-$  produced in the sample during a timed reaction compared with a heat inactivated control sample. NOS mediated NO production is halted by heat inactivation. Subsequent comparison of heat inactivated with non heat inactivated samples, over time, yields an indication of NOS activity.

Nitrate reductase is utilized for the enzymatic reduction of nitrate to nitrite.

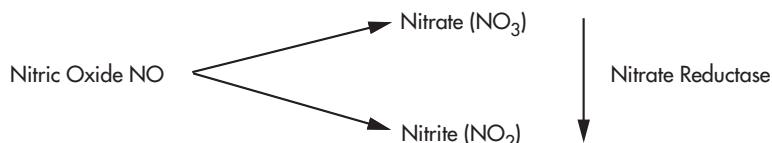

NADPH is an essential cofactor for the function of the NOS enzyme. Unfortunately, NADPH interferes with the Griess Reagents, the most commonly used reagents for nitrite detection. There are two common methods to reduce the NADPH interference. Very small concentrations of NADPH can be used, in conjunction with a catalytic system for recycling of the spent  $\text{NADP}^+$  back to NADPH; or the excess NADPH is removed or destroyed. This assay uses lactate dehydrogenase (LDH) to destroy the excess NADPH. Spectrophotometric quantitation of nitrite using Griess Reagents is straightforward and sensitive. In acidic solution, nitrite is converted to nitrous acid ( $\text{HNO}_2$ ), which diazotizes sulfanilamide. This sulfanilamide-diazonium salt is then reacted with N-(1-Naphthyl)-ethylenediamine to produce a chromophore which is measured at 540 nm.

## References

1. Busconi, L and Michel, T. 1993. *J. Biol. Chem.* 268, 8410.
2. Green, L.C., et al. 1982. *Anal. Biochem.* 126, 131.
3. Kobzik, L., et al. 1994. *Nature* 372, 546.
4. Moncada, S. 1992. *Acta Physiol. Scand.* 145, 201.
5. Nathan, C. 1992. *FASEB J.* 6, 3051.
6. Nims, R.W., et al. 1995. *Methods* 7, 48.
7. Pollock, J.S., et al. 1991. *Proc. Natl. Acad. Sci. U.S.A.* 88, 10480.

**NOS-Colorimetric Product Summary**

Intended Use:	Indirect determination of NOS activity by comparison of NO produced in active samples vs. heat inactivated samples. For Research Use Only. Not Intended For Use in Diagnostic Procedures.
Format:	80 microtitre tests
Kit Contents:	<ul style="list-style-type: none"><li>• Assay Buffer</li><li>• Nitrate Reductase</li><li>• Co-factor Preparation</li><li>• Nitrate Standard</li><li>• Lactate Dehydrogensae</li><li>• Color Reagent #1 (Sulfanilamide)</li><li>• Color Reagent #2 (N-(1-Naphthyl) Ethylenediamine Dihydrochloride)</li><li>• NADPH Tetrasodium Salt</li><li>• Microtitre Plate</li></ul>
Storage and Stability:	Seven months from date of manufacture when stored as specified.
Specimen Requirements:	Tissue homogenates and cell lysates.
Sensitivity:	5 nmol/mL

## L-Ergothioneine


26511  
26524

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number    26511 - 100 mg  
                           26524 - 250 mg

L-Ergothioneine is a natural thiol antioxidant which has been reported to have a number of different functions. It is present in tissues which are subject to high oxidative stress including the lens of the eye, the liver and red blood cells. It has been reported to block lipid peroxidation and protect against  $\gamma$ , UV radiation and singlet oxygen *in vitro*.

Form:                    White crystalline  
 Packaging Size:      100 mg and 250 mg  
 Source:                Synthetic  
 CAS Number:        497-30-3  
 Formula Weight:    229.3  
 Molecular Formula:  $C_9H_{15}N_3O_2S$   
 Purity:                >98% (HPLC)  
 Storage and Stability: Refrigerate. This product is stable for 1 year as supplied; storage and stability studies are ongoing.  
 References:         Hartman, P.E. 1990. *Meth. Enzy.* 186, 310-318.

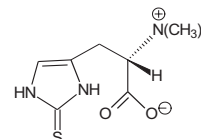

U.S. Patent No. 5,438,151



## Bioxytech® Cellular Glutathione Peroxidase Control

27617

*Control for Glutathione Peroxidase Enzyme Assay*

**Catalog Number** 27617 **For Research Use Only, Not Intended For Use in Diagnostic Procedures**

This companion product to the Bioxytech® GPx-340 Assay System enables investigators to assess the performance of the assay, assess recovery of enzyme activity in biological samples and normalize measured glutathione peroxidase (GPx) activity.

**Contents:**

- GPx control (2 vials, minimum 2500 mU/mL)
- Diluent

**Storage and Stability:** Freeze. This product is stable for six months from date of manufacture when stored frozen.

## Bioxytech® Glutathione Reductase Control

27618

*Control for Glutathione Reductase Enzyme Assay*

**Catalog Number** 27618 **For Research Use Only, Not Intended For Use in Diagnostic Procedures**

This companion product to the Bioxytech® GR-340 Assay System enables investigators to assess the performance of the assay, assess recovery of enzyme activity in biological samples and normalize measured glutathione reductase (GR) activity.

**Contents:**

- GR control (2 vials, minimum 200 mU/vial)
- Diluent

**Storage and Stability:** Freeze. This product is stable for six months from date of manufacture when stored frozen.

## Bioxytech® Cu/Zn SOD Control

27619

*Control for Cu/Zn SOD*

**Catalog Number** 27619 **For Research Use Only, Not Intended For Use in Diagnostic Procedures**

This companion product to the Bioxytech® SOD-525 Assay System enables investigators to assess the performance of the assay, assess recovery of enzyme activity in biological samples and normalize measured superoxide dismutase (SOD) activity. This control can also be used with other methodologies.


**Kit Contents:** Cu/Zn SOD Control; 5 mg, minimum 3000 U/mg protein (McCord-Fridovitch).

**Storage and Stability:** Freeze. This product is stable for at least one year from date of manufacture when stored frozen.

Catalog Number 27622

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalase (EC 1.11.1.6, 2 H<sub>2</sub>O<sub>2</sub> oxidoreductase) is a widely distributed enzyme that destroys hydrogen peroxide by dismutation (catalatic activity, Rxn 1) and also demonstrates peroxidatic activity (Rxn 2) (1,2). Hydrogen peroxide, a reactive oxygen species (ROS), is a toxic product of both normal aerobic metabolism and pathogenic ROS production. In the peroxidatic reaction, low molecular weight alcohols can serve as electron donors. The physiological significance of the peroxidatic activity of catalase is controversial (2).

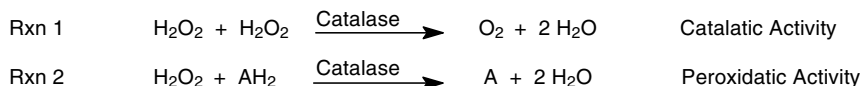

Eukaryotic catalases are heme enzymes but a manganese catalase has been described from prokaryotes (2). In humans, the highest levels of catalase are found in liver, kidney, and erythrocytes, where it is believed to account for the majority of hydrogen peroxide decomposition. Eukaryotic catalases bind NADPH, which stabilizes the enzyme and prevents the formation of Compound II, a form of catalase outside the normal catalytic cycle (3). Acatalasemia is a rare inherited condition in which there is little or no catalase produced (4,5). It is found most commonly in Asian populations, although European and other cases have been reported. Acatalasemic individuals occasionally present with oral ulcers (Takahara's Disease), but the condition is largely asymptomatic. Studies of a mouse model of acatalasemia have demonstrated the importance of catalase in preventing methemoglobin formation in erythrocytes (6).

### Catalase Control Method

- Contents:
- Control Diluent Surfactant in Phosphate Buffer, 30 mL.
  - Control Catalase, ≥ 200 U/vial, lyophilized, 2 vials.
- Items Required But NOT Provided:
- Deionized water.
  - Adjustable pipettes with disposable pipette tips.
- Control Preparation:
- Reconstitute the lyophilized contents of a Control vial with deionized water according to the vial label. See the Control vial label for the concentration of the resulting solution. Further dilute the Control with the provided Control Diluent to concentrations relevant to the study. Reconstituted and diluted catalase controls are stable for one week when stored at 4°C.

### References

1. Deisseroth A, Dounce AL (1970) Catalase: Physical and Chemical Properties, Mechanism of Catalysis, and Physiological Role, *Physiol. Rev.* 50, 319-375.
2. Zamocky M, Koller F (1999) Understanding the Structure and Function of Catalases: Clues from Molecular Evolution and In Vitro Mutagenesis, *Prog. Biophys. Mol. Biol.* 72, 19-66.
3. Kirkman HN, et al., (1987) The Function of Catalase-bound NADPH, *J. Biol. Chem.* 262, 660-666.
4. Aebi H, Suter H (1971) Acatalasemia, *Adv. Human Genet.* 2, 143-199.
5. Ogata M (1991) Acatalasemia, *Human Genet.* 86, 331-340.
6. Wakimoto M, et al., (1998) Determination of Glutathione Peroxidase Activity and Its Contribution to Hydrogen Peroxide Removal in Erythrocytes, *Acta Med. Okayama* 52, 233-237.

## Anti-Myeloperoxidase

24310

*Mouse Anti-human MPO Monoclonal Antibody*

<b>Catalog Number</b>	24310	<b>For Research Use Only, Not Intended For Use in Diagnostic Procedures</b>
Description:	Physiology and pharmacology of human PMN cells (neutrophils/secondary granules).	
Source:	Mouse	
Specificity:	<3% cross-reactivity with eosinophil peroxidase	
Form:	Liquid; Purity: >98%	
Packaging Size:	500 µg	
Storage and Stability:	Refrigerate. This product is stable for 1 year as supplied.	

## Anti-Lactoferrin

24311

*Mouse Anti-human Lf Monoclonal Antibody*

<b>Catalog Number</b>	24311	<b>For Research Use Only, Not Intended For Use in Diagnostic Procedures</b>
Description:	Physiology and pharmacology of human PMN cells (neutrophils/secondary granules).	
Source:	Mouse	
Specificity:	Human lactoferrin	
Form:	Liquid; Purity: >98%	
Packaging Size:	500 µg	
Storage and Stability:	Refrigerate. This product is stable for 1 year as supplied.	

## Anti-Nitrotyrosine

24312

*Sheep Anti-NT Polyclonal Antibody*

<b>Catalog Number:</b>	24312	<b>For Research Use Only, Not Intended For Use in Diagnostic Procedures</b>
Description:	Indication of peroxynitrite modification of proteins/oxidative stress	
Source:	Sheep	
Specificity:	Unmodified Human Serum Albumin: <ul style="list-style-type: none"> <li>• No significant binding with EIA</li> <li>• No signal in Western blot (at 1/5 dilution)</li> </ul> Nitrate Human Serum Albumin: <ul style="list-style-type: none"> <li>• 1/2 saturation at 1/30,000 dilution neat antibody in EIA (1 µg/mL antigen plated)</li> <li>• Binding of antibody completely blocked by 1 mM 3-nitrotyrosine</li> <li>• Optimal dilution in Western blotting of (1/100)</li> <li>• Western blot signal blocked by 1 mM 3-nitrotyrosine</li> </ul>	
Form:	Plasma diluted 1:1 with PBS/0.01% thimerasol and 0.2 µm filtered.	
Packaging Size:	1500 µL	
Storage and Stability:	Freeze. This product is stable for 2 years as supplied.	

## Anti-Catalase

24316

*Rabbit Anti-Human Catalase Polyclonal Antibody*

<b>Catalog Number</b>	24316	<b>For Research Use Only, Not Intended For Use in Diagnostic Procedures</b>
Description:	IgG fraction	
Buffer:	Phosphate buffered saline	
Source:	Rabbit	
Specificity:	Single arc by immunoelectrophoresis (IEP) vs. human red cell lysate. Single arc by IEP vs. 10 mg human erythrocyte catalase.	
Form:	Liquid Preservative: 0.1% NaN <sub>3</sub>	
Usage:	Radial immunodiffusion (RID). Working Concentration: 455 mg of antigen precipitated per mL of antiserum. Avoid freeze/thaw cycles.	
Packaging Size:	1000 µL	
Storage:	Freeze	

## Anti-Cytochrome P450 2E1

24317

*Rabbit Anti-Human P450 2E1 Polyclonal Antibody*

<b>Catalog Number</b>	24317	<b>For Research Use Only, Not Intended For Use in Diagnostic Procedures</b>
Source:	Rabbit. Immunogen: Recombinant human P450 2E1 expressed in E. coli.	
Specificity:	By Western blot, recognizes a band (P450 2E1) in human liver microsomes. Only slightly cross-reacts with recombinant human P450 1A1 after prolonged incubation. Does not cross-react with recombinant human cytochromes P450 1A2, 3A4 or 2C10.	
Form:	Liquid. Purified rabbit IgG. Preservative: 0.05% NaN <sub>3</sub>	
Usage:	A dilution of 1:1000 for Western blots will yield a positive result with ≥ 0.2 mg of recombinant human cytochrome P450 2E1 using alkaline phosphatase-conjugated anti-rabbit IgG as secondary antibody. Concentration: 5 mg/mL in 20 mM phosphate buffer, pH 8.0. Avoid freeze/thaw cycles.	
Packaging Size:	100 µL	
Storage:	Freeze	
Reference:	Gillarm, E.M., et al. 1994. Arch. Biochem. Biophys. 312, 59.	

## Anti-Cytochrome P450 3A4

24318

*Rabbit Anti-Human P450 3A4 Polyclonal Antibody*

<b>Catalog Number</b>	24318	<b>For Research Use Only, Not Intended For Use in Diagnostic Procedures</b>
Source:	Rabbit. Immunogen: Recombinant human P450 3A4 expressed in E. coli.	
Specificity:	By Western blot, reacts with human recombinant P450 3A4 and with P450 3A4 in human liver microsomes.	
Form:	Liquid. Purified rabbit IgG. Preservative: 0.05% NaN <sub>3</sub>	
Usage:	A dilution of 1:1000 for Western blots will yield a positive result with ≥ 0.2 mg of recombinant human cytochrome P450 3A4 using alkaline phosphatase-conjugated anti-rabbit IgG as secondary antibody. Avoid freeze/thaw cycles.	
Packaging Size:	100 µL	
Storage:	Freeze	



## Anti-Glutathione S-Transferase A1-1

24319

*Rabbit Anti-Human GST A1-1 Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures**

<b>Catalog Number</b>	24319
Description:	Polyclonal antibody produced by immunizing rabbits with recombinant human GST A1-1, expressed in <i>E. coli</i> (Cat. No. 25434).
Source:	Rabbit
Specificity:	Specific for the $\alpha$ -form of human and rat GST. Reacts with recombinant human GST A1-1. Does not react with $\mu$ or $\rho$ GST's.
Form:	Liquid. Undiluted serum. Preservative: 0.05% NaN <sub>3</sub>
Usage:	Immunoblotting. A dilution of 1/1000 will yield a positive result with $\geq 0.2$ mg of human A1-1 GST, using alkaline phosphatase-conjugated anti-rabbit IgG as secondary antibody. Avoid freeze/thaw cycles.
Packaging Size:	100 $\mu$ L
Storage:	Freeze
Reference:	Holmes, W., et al. 1990. <i>Br. J. Cancer</i> 62, 209.

## Anti-Glutathione S-Transferase M1-1

24320

*Rabbit Anti-Human GST M1-1 Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures**

<b>Catalog Number</b>	24320
Description:	Polyclonal antibody produced by immunizing rabbits with recombinant human GST M1-1, expressed in <i>E. coli</i> (Cat. No. 25435).
Source:	Rabbit
Specificity:	Specific for the $\mu$ -form of human and rat GST. Reacts with recombinant human GST M1-1. Does not react with $\alpha$ or $\rho$ GST's.
Form:	Liquid. Undiluted serum. Preservative: 0.05% NaN <sub>3</sub>
Usage:	Immunoblotting. A dilution of 1/1000 will yield a positive result with $\geq 0.2$ mg of human M1-1 GST, using alkaline phosphatase-conjugated anti-rabbit IgG as secondary antibody. Avoid freeze/thaw cycles.
Packaging Size:	100 $\mu$ L
Storage:	Freeze
Reference:	Holmes, W., et al. 1990. <i>Brit. J. Cancer</i> 62, 209.

## Anti-Glutathione S-Transferase P1-1

24321

*Rabbit Anti-Human GST P1-1 Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures**

<b>Catalog Number</b>	24321
Description:	Polyclonal antibody produced by immunizing rabbits with recombinant human GST P1-1, expressed in <i>E. coli</i> (Cat. No. 25436).
Source:	Rabbit
Specificity:	Specific for the $\rho$ -form of human and rat GST. Reacts with recombinant human GST P1-1. Does not react with $\alpha$ or $\mu$ GST's.
Form:	Liquid. Undiluted serum. Preservative: 0.05% NaN <sub>3</sub>
Usage:	Immunoblotting. A dilution of 1/1000 will yield a positive result with $\geq 0.2$ mg of human P1-1 GST, using alkaline phosphatase-conjugated anti-rabbit IgG as secondary antibody. Avoid freeze/thaw cycles.
Packaging Size:	100 $\mu$ L
Storage:	Freeze
Reference:	Holmes, W., et al. 1990. <i>Br. J. Cancer</i> 62, 209.

## Anti-Cu/Zn - Superoxide Dismutase

24322

*Sheep Anti-Human Cu/Zn-SOD Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 24322

Superoxide dismutase Cu/Zn enzyme (Cu/Zn SOD) is a homodimeric metalloenzyme of aerobic organisms, comprised of subunits of approximately 16 kDa. Found in a variety of cell types, including erythrocytes, liver, and brain, it serves to protect cells against the oxidative damage of free radicals by catalyzing the conversion of superoxide anions ( $O_2^-$ ) to molecular oxygen ( $O_2$ ) and hydrogen peroxide ( $H_2O_2$ ). This anti-Cu/Zn SOD polyclonal antibody is supplied as an IgG fraction obtained by DEAE ion-exchange chromatography.

Source: Sheep  
 Immunogen: Highly purified Cu/Zn SOD enzyme from human erythrocytes  
 Specificity: Monospecific as determined by immunoelectrophoresis against human red blood cell lysate. Confirmed by immunodiffusion vs. a human red blood cell lysate and a known anti-Cu/Zn SOD antibody. No cross-reactivity with Mn SOD by double diffusion or with human serum.  
 Form: Liquid, passed through a 0.2  $\mu$ m filter. Buffer: Glycine-buffered saline, pH 7.4. Preservatives: 0.1%  $NaN_3$ , 0.1%  $\epsilon$ -amino-n-caproic acid, 0.01% benzamidine, 1 mM EDTA.  
 Usage: Immunohistochemistry, 1:20 – 1:200; Immunoblotting: 1:1000. Avoid freeze/thaw cycles.  
 Packaging Size: 1000  $\mu$ L  
 Storage: Freeze

## Anti-Mn Superoxide Dismutase

24323

*Sheep Anti-Human Mn-SOD Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 24323

Superoxide dismutase enzyme Mn SOD is a homotetrameric metallo enzyme of aerobic organisms, comprised of subunits of approximately 25 kDa. Localized primarily in the mitochondria, it serves to protect cells against the oxidative damage of free radicals by catalyzing the conversion of superoxide anion ( $O_2^-$ ) to molecular oxygen ( $O_2$ ) and hydrogen peroxide ( $H_2O_2$ ). This anti-SOD (Mn) polyclonal antibody is supplied as an IgG fraction prepared by DEAE ion-exchange chromatography.

Source: Sheep  
 Immunogen: Highly purified Mn SOD enzyme from human liver  
 Specificity: Monospecific as determined by immunoelectrophoresis against a human liver cell lysate. No cross-reactivity with SOD (Cu/Zn) by double diffusion or with human serum.  
 Form: Liquid, passed through a 0.2  $\mu$ m filter. Preservatives: 0.1%  $NaN_3$ , 0.1%  $\epsilon$ -amino-n-caproic acid, 0.01% benzamidine, 1 mM EDTA. Buffer: Glycine buffered saline, pH 7.4  
 Usage: Immunohistochemistry, 1:20–1:200. Avoid freeze/thaw cycles.  
 Packaging Size: 1000  $\mu$ L  
 Storage: Freeze

## Anti-4-Hydroxy-2-nonenal

24325

*Mouse Anti-4-HNE Monoclonal Antibody*

24327

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

<b>Catalog Number</b>	24325 - 100 µg/vial (lyophilized) 24327 - 20 µg/vial (lyophilized)
Source:	Mouse (IgG1)
Immunogen:	4-HNE-modified keyhole limpet hemocyanine. Use Immunohistochemistry, Western blotting Dilution: 25 µg/mL for Immunohistochemistry, 15 µg/mL for Western blotting
Form:	Each vial contains 100µg (20 µg) of Anti-4-HNE monoclonal antibody. If dissolved in 1 mL (0.2mL) of distilled water, the composition of the solution will be 100 µg/mL with 0.05M TBS containing 1% BSA
Specificity:	This antibody shows negligible reactivity with proteins that were treated with other aldehydes such as: 2-nonenal, 2-hexenal, 1-hexenal, 4-hydroxy-2-hexenal, formaldehyde or glutaraldehyde.
Storage:	Store at <4°C
Stability:	3 years at -20°C

## Anti-8-Hydroxy-2'-deoxyguanosine

24326

*Mouse Anti-8-OHdG Monoclonal Antibody*

24328

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

<b>Catalog Number</b>	24326 - 100 µg/vial (lyophilized) 24328 - 20 µg/vial (lyophilized)
Source:	Mouse (IgG1)
Immunogen:	8-OHdG conjugated keyhole limpet hemocyanin. Use Immunohistochemistry Dilution 5-10 µg/mL
Form:	Each vial contains 100µg (20 µg) of Anti-8-OHdG monoclonal antibody. If dissolved in 1 mL (0.2mL) of distilled water, the composition of the solution will be 100 µg/mL with 0.05M TBS containing 1% BSA
Specificity:	21 Analogues of 8-OHdG were examined with only two exhibiting mild cross-reactivity. 8-sulfhydrylguanosine and 8-OHG demonstrate some cross-reactivity at high concentrations (at least 2 orders of magnitude were required for positive cross-reactivity).
Storage:	Store at <4° C
Stability:	3 years at -20° C

## Anti-Endothelial Nitric Oxide Synthase

25310

*Rabbit anti Bovine eNOS Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 25310

The antibody recognizes endothelial nitric oxide synthase (eNOS, 140 kDa) protein in human, rat and mouse. The immunogen used was a bovine eNOS synthetic peptide (599–613) plus C-terminal Cys (PYNSSPRPEQHKSYKC) conjugated to KLH. No cross-reactivity to iNOS or bNOS.

Source:	Rabbit
Form:	Undiluted serum. Preservative: 0.02% NaN <sub>3</sub>
Usage:	1:250 in immunocytochemistry; 1:1,000 in Western blotting. Avoid freeze/thaw cycles.
Packaging Size:	100 µL
Storage:	Freeze

## Anti-Inducible Nitric Oxide Synthase

25311

*Rabbit anti Mouse iNOS Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 25311

This antibody recognizes inducible nitric oxide synthase (iNOS, 130 kDa) protein in human, rat and mouse. The immunogen used was mouse macrophage NOS C-terminal peptide (1131-1144) plus N-terminal Cys (CKKGSALleepKATRL) conjugated to KLH. Exhibits no cross-reactivity with bNOS or eNOS.

Source: Rabbit  
 Form: Undiluted serum. Preservative: 0.02% NaN<sub>3</sub>  
 Usage: 1:500 in immunocytochemistry; 1:500 in immunohistochemistry; 1:2,000 in Western blotting. Avoid freeze/thaw cycles.  
 Packaging Size: 100 µL  
 Storage: Freeze

## Anti-Brain Nitric Oxide Synthase

25312

*Rabbit anti Rat bNOS Polyclonal Antibody***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number** 25312

The antibody recognizes brain nitric oxide synthase (bNOS, 155 kDa) protein in bovine, primates, mouse and rat. The immunogen used was a rat bNOS synthetic peptide (724-739) corresponding to the calmodulin-binding domain, plus C-terminal Cys (TKRRAIGFKKLAEAVKC) conjugated to KLH. Exhibits no cross-reactivity to iNOS or eNOS.

Source: Rabbit  
 Form: Undiluted serum. Preservative: 0.02% NaN<sub>3</sub>  
 Usage: 1:800 in immunohistochemistry; 1:1,000 in Western blotting. Avoid freeze/thaw cycles.  
 Packaging Size: 100 µL  
 Storage: Freeze

## Bovine Cu/Zn Superoxide Dismutase (bSOD)

25447 25414  
25415 25416

For Research Use Only, Not Intended For Use in Diagnostic Procedures

**Catalog Number** 25447 - 25 mg  
25414 - 250 mg  
25415 - 500 mg  
25416 - 1 g

Bovine Cu/Zn SOD can be used in research as a reference for quantification of SOD activity in biological samples. It can also be used as a reference for quality control studies of substances which contain SOD activity.

**Form:** Lyophilized solid  
**Source:** Bovine liver  
**Packaging Size:** Various  
**Formula Weight:** 32,500  
**Purity:** >99%  
**Specific Activity:** >3000 U/mg (McCord-Fridovitch)  
**Storage and Stability:** Freeze. This product is stable for 1 year as supplied.

Catalase, *Aspergillus niger*

25430

For Research Use Only, Not Intended For Use in Diagnostic Procedures

**Catalog Number** 25430

Long-acting, extremely stable form of catalase that is active from pH 2–12. Composed of four protein subunits, each containing a heme (Fe<sup>3+</sup>-protoporphyrin) group bound to its active site. Each subunit also contains one molecule of bound NADPH that helps to stabilize the enzyme. Catalyzes the reaction  $2\text{H}_2\text{O}_2 \rightarrow 2\text{H}_2\text{O} + \text{O}_2$ . Catalase, along with superoxide dismutase and glutathione peroxidase, controls the level of oxygen-derived free radicals in mammalian cells, and together may function as a somatic oxidant defense. Prevents apoptosis in cultured rat ovarian follicles.

**Form:** Light brown lyophilized solid  
**Packaging Size:** 100 KU  
**CAS Number:** 9001-05-2  
**RTECS:** FI4378000  
**EC Number:** 1.11.1.6  
**Formula Weight:** 250,000  
**Specific Activity:** 10 KU/mg protein  
**Activity:** >4 KU/mg dry weight  
**Contaminants:** Microbial contamination:  $\leq 50$  cfu/g dry weight  
**Unit Definition:** 1 KU = 1000 units. One unit is defined as the amount of enzyme that will decompose 1.0  $\mu\text{mol}$  of  $\text{H}_2\text{O}_2$  per minute at 30°C, pH 7.0.  
**Solubility:**  $\text{H}_2\text{O}$  (20 mg/ml) and physiological buffers  
**Storage:** Freeze. Following reconstitution, aliquot and refreeze for long term storage or store at room temperature for short term storage. This product is stable for 2 years as supplied. Aqueous stock solutions are stable for several weeks at room temperature or for several months frozen.  
**References:** Tilly, J.L., and Tilly, K.I. 1995. *Endocrinology* 136, 242.  
Kassim, E.A. 1982. *Microbiol. Immunol.* 26, 449.  
Kikuchi-Torii, K., et al. 1992. *J. Biochem.* 92, 1149.  
Beers, R.F., and Sizer, I.W. 1952. *J. Biol. Chem.* 195, 133.

## Catalase, Bovine Liver

25431

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 25431

Catalase, along with superoxide dismutase and glutathione peroxidase, controls the levels of oxygen-derived free radicals in mammalian cells. Together these molecules may function as a somatic oxidant defense. Prevents apoptosis in cultured rat ovarian follicles.

Form: Green lyophilized solid  
 Packaging Size: 5 MU  
 CAS Number: 9001-05-2  
 RTECS: F14378000  
 EC Number: 1.11.1.6  
 Formula Weight: 247,000  
 Specific Activity: 40,000 units/mg protein  
 Activity:  $\geq 5000$  units/mg dry weight  
 Unit Definition: 1 MU = 1,000,000 units. One unit is defined as the amount of enzyme that will decompose 1.0  $\mu\text{mol}$  of  $\text{H}_2\text{O}_2$  per minute at 25°C, pH 7.0.  
 Solubility:  $\text{H}_2\text{O}$  and aqueous buffers  
 Storage and Stability: Freeze. Protect from moisture. Following reconstitution, aliquot and refreeze. This product is stable for 2 years as supplied. Stock solutions are stable for at least 1 month when frozen.  
 References: Tilly, J.L., and Tilly, K.I. 1995. *Endocrinology* 136, 242.  
 Beers, R.F., and Sizer, I.W. 1952. *J. Biol. Chem.* 195, 133.

## Catalase, Human Erythrocyte

25432

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 25432

Catalase, along with superoxide dismutase and glutathione peroxidase, controls the levels of oxygen-derived free radicals in mammalian cells. Together these molecules may function as a somatic oxidant defense. Prevents apoptosis in cultured rat ovarian follicles.

Form: Liquid. Supplied in 50 mM Tris, pH 8.0.  
 Source: Prepared from blood that has been tested to be negative for HBsAg and for HIV and HCV antibodies. It is recommended that the product be handled using established good laboratory practices since no test method can guarantee that products derived from human blood will not transmit infection.  
 Packaging Size: 1 mg  
 CAS Number: 9001-05-2  
 RTECS: F14378000  
 EC Number: 1.11.1.6  
 Formula Weight: 256,000  
 Purity:  $\geq 95\%$  by SDS-Page  
 Specific Activity:  $>50,000$  units/mg  
 Unit Definition: One unit is defined as the amount of enzyme that will decompose 1.0  $\mu\text{mol}$  of  $\text{H}_2\text{O}_2$  per minute at 25°C, pH 7.0.  
 Storage and Stability: Freeze. Avoid freeze/thaw cycles. This product is stable for 2 years as supplied.  
 References: Tilly, J.L., and Tilly, K.I. 1995. *Endocrinology* 136, 242.  
 Bonaventura, J., et al. 1972. *Arch. Biochem. Biophys.* 150, 606.

## Glutathione Reductase, Yeast

25433

*GR*

Catalog Number 25433

For Research Use Only, Not Intended For Use in Diagnostic Procedures

A ubiquitous enzyme that catalyzes the oxidation of oxidized glutathione (GSSG) to glutathione (GSH).

Form:	Yellow crystalline suspension in 3.2 M (NH <sub>4</sub> ) <sub>2</sub> SO <sub>4</sub> , pH 6.0.
Packaging Size:	600 U
CAS Number:	9001-48-3
EC Number:	1.6.4.2
Formula Weight:	118,000
Contaminants:	NADPH oxidase and G-6-PDH: <0.01%; 6-PGDH: <0.1%
Specific Activity:	≥ 100 units/mg protein
Unit Definition:	One unit is defined as the amount of enzyme that will hydrolyze 1.0 mmol of oxidized glutathione per minute at 25°C, pH 7.6.
Solubility:	Aqueous buffers.
Storage and Stability:	Refrigerate. Do not freeze. This product is stable for 1 year as supplied.
Reference:	Smith, I.K., et al. 1988. <i>Anal. Biochem.</i> 175, 408.

## Glutathione-S-Transferase A1-1 Recombinant Human

25434

*rhGST A1-1*

Catalog Number 25434

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Affinity purified, biologically active, full-length recombinant GST A1-1 that possesses physical and catalytic properties indistinguishable from the native enzyme. GSTs catalyze the formation of thioester conjugates between glutathione and reactive xenobiotic compounds and are believed to be important for the cellular response to oxidative stress. Suitable for pro-carcinogen activation studies.

Form:	Liquid. Supplied in 50 mM Tris-HCl, pH 7.5, 50 mM NaCl, 1 mM DTT, 1 mM EDTA, and 50% glycerol.
Packaging Size:	100 µg
Formula Weight:	50,000
Purity:	>95% by SDS-PAGE
Specific Activity:	50 units/mg protein
Unit Definition:	One unit is defined as the amount of enzyme that will conjugate 1.0 mmol of 1-chloro-2,4-dinitrobenzene (CDNB) to reduced glutathione per minute at 20°C, pH 6.5.
Storage and Stability:	Freeze. This product is stable for 1 year as supplied.
References:	Kolm, R.H., et al. 1995. <i>Biochem. J.</i> 311, 453. Kolm, R.H., et al. 1995. <i>Protein Expr. Purif.</i> 6, 265. Stenberg, G., et al. 1992. <i>Protein Expr. Purif.</i> 3, 80. Widersten, M., et al. 1991. <i>Biochem. J.</i> 276, 519.

## Glutathione-S-Transferase M1-1 Recombinant Human

25435

*rhGST M1-1*

Catalog Number 25435

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

Affinity purified, biologically active, full length recombinant GST M1-1 that possesses physical and catalytic properties indistinguishable from the native enzyme. Corresponds to the type b allelic variant and GST Y purified from human liver. GSTs catalyze the formation of thioester conjugates between glutathione and reactive xenobiotic compounds and are believed to be important for the cellular response to oxidative stress. Suitable for procarcinogen activation studies.

Form: Liquid. Supplied in 50 mM Tris-HCl, pH 7.5, 50 mM NaCl, 1 mM DTT, 1 mM EDTA, and 50% glycerol.

Packaging Size: 100 µg

Formula Weight: 50,000

Purity: ≥ 95% by SDS-PAGE

Specific Activity: 100 units/mg protein

Unit Definition: One unit is defined as the amount of enzyme that will conjugate 1.0 mmol of 1-chloro-2,4-dinitrobenzene (CDNB) to reduced glutathione per minute at 20°C, pH 6.5.

Storage and Stability: Freeze. This product is stable for 1 year as supplied.

References: Kolm, R.H., et al. 1995. *Biochem. J.* 311, 453.  
Kolm, R.H., et al. 1995. *Protein Expr. Purif.* 6, 265.  
Stenberg, G., et al. 1992. *Protein Expr. Purif.* 3, 80.  
Widersten, M., et al. 1991. *Biochem. J.* 276, 519.

## Glutathione-S-Transferase P1-1 Recombinant Human

25436

*rhGST P1-1*

Catalog Number 25436

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

Affinity purified, biologically active, full length recombinant GST P1-1. GSTs catalyze the formation of the thioester conjugates between glutathione and reactive xenobiotic compounds and are believed to be important for the cellular response to oxidative stress. Suitable for procarcinogen studies. Higher levels of GST P1-1 are reported in many human tumors, including melanomas. GST P1-1 is widely distributed throughout the body, except for the liver.

Form: Liquid. Supplied in 50 mM Tris-HCl, pH 7.5, 50 mM NaCl, 1 mM DTT, 1 mM EDTA, and 50% glycerol.

Packaging Size: 100 µg

Formula Weight: 50,000

Purity: ≥ 95% by SDS-PAGE

Specific Activity: 75 units/mg protein

Unit Definition: One unit is defined as the amount of enzyme that will conjugate 1.0 mmol of 1-chloro-2,4-dinitrobenzene (CDNB) to reduced glutathione per minute at 20°C, pH 6.5.

Storage and Stability: Freeze. This product is stable for 1 year as supplied.

References: Kolm, R.H., et al. 1995. *Biochem. J.* 311, 453.  
Stenberg, G., et al. 1992. *Protein Expression Purif.* 3, 80.  
Widersten, M., et al. 1991. *Biochem. J.* 276, 519.



## Myeloperoxidase, Human Polymorphonuclear Leukocytes

25437

*MPO*

Catalog Number 25437

For Research Use Only, Not Intended For Use in Diagnostic Procedures

A lysosomal heme protein that catalyzes oxidations by hydrogen peroxide, including MPO-chloride-mediated killing of microbes and tumor cells, inactivation of chemotactic factors, cross-linking of proteins, and iodination of proteins.

Form: Solid lyophilized from 50 mM sodium acetate, pH 6.0, and 100 mM NaCl.  
 Packaging Size: 100 µg  
 Source: Prepared from blood that tested negative for HBsAg and for HIV and HCV antibodies. It is recommended that this product be handled using established good laboratory practices since no test method can guarantee that products derived from human blood will not transmit infection.  
 CAS Number: 9003-99-0  
 EC Number: 1.11.1.7  
 Formula Weight: 118,000  
 Purity: ≥ 95% by SDS-Page  
 Specific Activity: 50-200 units/mg protein  
 Unit Definition: One unit is defined as the amount of enzyme that will hydrolyze 1.0 mmol of H<sub>2</sub>O<sub>2</sub> per minute at 25°C, pH 6.0.  
 Solubility: Reconstitute with 1 mL of 50 mM sodium acetate, pH 6.0, and 100 mM NaCl. Enzyme may precipitate out of solution in low ionic strength buffers or in distilled H<sub>2</sub>O.  
 Storage and Stability: Freeze. Following reconstitution, aliquot and refreeze at -20°C or below. This product is stable for 2 years as supplied. Stock solutions are stable for 6-12 months at -20°C or below.  
 References: Panasenko, O.M., et al. 1994. Free Radical Biol. Med. 16, 143.  
 Mulligan, M.S., et al. 1992. Br. J. Pharmacol. 107, 1159.  
 Sharonov, B.P., and Churilova, I.V. 1992. Biochem. Biophys. Res. Commun. 189, 1129.

## Brain Nitric Oxide Synthase, Recombinant Rat

27510

*bNOS*

Catalog Number 27510

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Constitutive rat brain NOS isoform. Produced by a baculovirus expression system in the presence of hemin. Useful for studies involving NOS regulation, inhibitor screening and kinetic analysis.

Form: Liquid. Supplied as 0.1 mg/ml in 50 mM triethanolamine-HCl, 10 mM β-mercaptoethanol, 1 mM CHAPS, 500 mM EDTA, pH 7.0.  
 Packaging Size: 10 µg  
 CAS Number: 125978-95-2  
 EC Number: 1.14.13.39  
 Formula Weight: 150,000  
 Purity: ≥ 95% by SDS-PAGE  
 Specific Activity: ≥ 500 nmol L-citrulline/mg/min at 37°C.  
 Storage and Stability: Deep freeze. Avoid freeze/thaw cycles. After initial thawing, aliquot and refreeze (-70°C). This product is stable for 2 years as supplied.  
 References: Nakane, M., et al. 1995. Biochem. Biophys. Res. Commun. 206, 511.  
 Harteneck, C., et al. 1994. Biochem. J. 304, 683.  
 Richards, M.K., and Marletta, M.A. 1994. Biochemistry 33, 14723.  
 Charles, I.G., et al. 1993. Biochem. Biophys. Res. Commun. 196, 1481.

## Constitutive Nitric Oxide Synthase Recombinant Rat Neuronal 27511

cNOS

**Catalog Number** 27511

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

Constitutive rat brain NOS isoform cNOS. Recombinant enzyme from the 100,000 x g supernatant of a baculovirus expression system. Useful for studies involving NOS regulation, inhibitor screening and kinetic analysis.

**Form:** Liquid. Supplied in 50 mM HEPES, pH 7.4, 100 mM DTT, and 10% glycerol.  
**Packaging Size:** 50 U  
**Formula Weight:** 150,000  
**Specific Activity:** 35–80 units/mg protein  
**Unit Definition:** One unit is defined as the amount of enzyme that will produce 1.0 nmol of nitric oxide per minute in the presence of calmodulin at 37°C, pH 7.5.  
**Storage:** Deep freeze. Avoid freeze/thaw cycles. This product is stable for 1 year as supplied.  
**References:** Abu-Soud, H.M., et al. 1994. *J. Biol. Chem.* 269, 32047.  
 Abu-Soud, H.M., et al. 1994. *J. Biol. Chem.* 269, 32318.  
 Richards, M.K., et al. 1994. *Biochemistry* 33, 14723.  
 Sheta, E.A., et al. 1994. *J. Biol. Chem.* 269, 15147.

## Constitutive Nitric Oxide Synthase Recombinant Human, High Purity 27512

cNOS

**Catalog Number** 27512

**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

Constitutive rat brain NOS isoform. Catalyzes the production of nitric oxide (NO) by conversion of L-arginine to L-citrulline. Useful for studies involving NOS regulation, inhibitor screening and kinetic analysis.

**Form:** Liquid. Supplied in 20% glycerol, containing: 50 mM Hepes, 100 mM NaCl, and 10 mM tetrahydro-L-biopterin, at pH 7.4.  
**Source:** Recombinant enzyme purified from a baculovirus expression system.  
**Packaging Size:** 10 U  
**Formula Weight:** 150,000  
**Purity:** ≥ 95% by SDS-Page  
**Specific Activity:** 800-1300 units/mg  
**Unit Definition:** One unit is defined as the amount of enzyme that will produce 1.0 nmol of nitric oxide per minute in the presence of calmodulin at 37°C, pH 7.5.  
**Storage:** Deep freeze (-70°C). Avoid freeze/thaw cycles.  
**Stability:** This product is stable for 1 year as supplied.  
**References:** Abu-Soud, H.M., et al. 1994. *J. Biol. Chem.* 269, 32047.  
 Abu-Soud, H.M., et al. 1994. *J. Biol. Chem.* 269, 32318.  
 Richards, M.K., et al. 1994. *Biochemistry* 33, 14723.  
 Sheta, E.A., et al. 1994. *J. Biol. Chem.* 269, 15147.

## Inducible Nitric Oxide Synthase, Mouse Macrophage

27513

*iNOS***For Research Use Only, Not Intended For Use in Diagnostic Procedures****Catalog Number:** 27513

Inducible mouse brain NOS isoform iNOS. Catalyzes the production of nitric oxide by conversion of L-arginine to L-citrulline. Useful for studies involving NOS regulation, inhibitor screening and kinetic analysis.

**Form:** Liquid. Supplied in 50 mM HEPES and 10% glycerol, pH 7.4.  
**Source:** Recombinant enzyme from the 100,000 x g supernatant of immunostimulated murine macrophage (RAW 264.7 cells).  
**Packaging Size:** 50 U  
**Formula Weight:** 130,000  
**Specific Activity:** 7 units/mg  
**Unit Definition:** One unit is defined as the amount of enzyme that will produce 1.0 nmol of nitric oxide per minute at 37°C, pH 7.4.  
**Storage and Stability:** Deep freeze. Avoid freeze/thaw cycles. This product is stable for 2 years as supplied.  
**References:** Stevens-Truss, R., et al. 1995. *Biochemistry* 34, 15638.  
Marletta, M.A., 1994. *J. Med. Chem.* 37, 1899.  
White, K.A., et al. 1992. *Biochemistry* 31, 6627.  
Hevel, J.M. et al. 1991. *J. Biol. Chem.* 266, 22789.



## 4-Hydroxy-non-2-enal-diethylacetal

26515


4-HNE-DEA

26516

26595

**Catalog Number**26595 - 50 mg  
26515 - 100 mg  
26516 - 1.0 g**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

Description: A precursor to one of the major metabolites of lipid hydroperoxide degradation.  
 Form: Liquid  
 Source: Synthetic  
 Packaging Size: 10 mg, 100 mg, 1.0 g  
 CAS Number: 18445-69-7  
 Formula Weight: 230.3  
 Molecular Formula:  $C_{13}H_{26}O_3$   
 Purity:  $\geq 98\%$  (GC)  
 Storage and Stability: Freeze. This product is stable for 2 years when frozen.

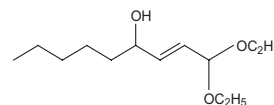

## 4-Hydroxy-hex-2-enal-diethylacetal

26518


4-HHE-DEA

26519

26594

**Catalog Number**26594 - 50 mg  
26518 - 100 mg  
26519 - 1.0 g**For Research Use Only, Not Intended For Use in Diagnostic Procedures**

Description: A precursor to one of the major metabolites of lipid hydroperoxide degradation.  
 Form: Liquid  
 Source: Synthetic  
 Packaging Size: 10 mg, 100 mg, 1.0 g  
 CAS Number: 18445-66-4  
 Formula Weight: 188.3  
 Molecular Formula:  $C_{10}H_{20}O_3$   
 Purity:  $\geq 96\%$  (GC)  
 Storage and Stability: Freeze. This product is stable for 2 years when frozen.

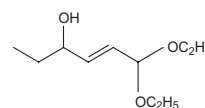

## 8-Hydroxyguanosine


26525

8-OHG

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 26525

Description: Product of oxidative damage to RNA.  
 Form: Crystalline solid  
 Packaging Size: 1 mg  
 CAS Number: 3868-31-3  
 Formula Weight: 299.2  
 Molecular Formula:  $C_{10}H_{13}N_5O_6$   
 Purity:  $\geq 98\%$  by TLC  
 Solubility:  $H_2O$   
 Storage and Stability: Refrigerate. This product is stable for 3 years as supplied.  
 References: Takeuchi, T., and Morimoto, K. 1993. *Carcinogenesis* 14, 1115.  
 Csillag, C., and Aldhous, P. 1992. *Science* 258, 1875.  
 Mo, Y., et al. 1992. *Proc. Natl. Acad. Sci. USA* 89, 11021.

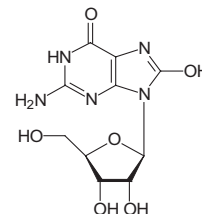

## 4-Hydroxy-2-nonenal

26526


4-HNE

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Catalog Number 26526

A major aldehydic product formed by peroxidation of  $\omega$ -6-unsaturated fatty acids that is regarded as a specific marker of lipid peroxidation. Inhibits proliferation and induces differentiation of HL-60 human leukemic cells. Also induces cell death in murine alveolar macrophages. An inhibitor of state 3 respiration at micromolar levels that causes a transient increase in cytosolic  $Ca^{2+}$ . Irreversibly inhibits  $Na^+-K^+$ -ATPase activity ( $IC_{50} = 120$  mM).

Form: Liquid. Supplied in ethanol.  
 Packaging Size: 1 mg  
 CAS Number: 75899-68-2  
 RTECS: RA8510000  
 Formula Weight: 156.2  
 Purity:  $\geq 98\%$  by HPLC  
 Molecular Formula:  $C_9H_{16}O_2$   
 Solubility: Ethanol  
 Storage and Stability: Deep freeze. This product is stable for 3 years as supplied.  
 References: Carini, R., et al. 1996. *Biochem. Biophys. Res. Commun.* 218, 772.  
 Li, L., et al. 1996. *Toxicol. Appl. Pharmacol.* 139, 135. Siems, W.G., et al. 1996. *Free Radic. Res.* 20, 215.  
 Ullrich, O., et al. 1996. *Free Radic. Res.* 24, 421. van Kuijk, F.J.G.M., et al. 1995. *Anal. Biochem.* 224, 420.  
 Esterbauer, H., et al. 1991. *Free Radic. Biol. Med.* 11, 81.

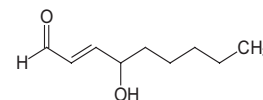

## 3-Nitro-L-Tyrosine


26527

Catalog Number 26527

For Research Use Only, Not Intended For Use in Diagnostic Procedures

A marker for peroxynitrite. A powerful oxidant and cytotoxic agent. Nitrotyrosine is the major product obtained by the nitration of tyrosine by peroxynitrite.

Form: Yellow-green solid  
 Packaging Size: 1 g  
 CAS Number: 621-44-3  
 Formula Weight: 226.2  
 Molecular Formula:  $C_9H_9N_2O_5$   
 Purity:  $\geq 98\%$  by TLC

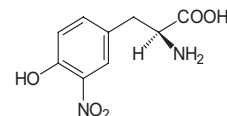

## 8-Hydroxy-2'-deoxyguanosine

26571


8-OHdG

Catalog Number 26571

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Product of oxidative damage to DNA.

Form: Crystalline solid  
 Packaging Size: 1 mg  
 CAS Number: 88847-89-6  
 Formula Weight: 283.2  
 Solubility:  $H_2O$   
 Molecular Formula:  $C_{10}H_{13}N_5O_5$   
 Purity:  $\geq 98\%$  by elemental analysis  
 Storage and Stability: Refrigerate. This product is stable for 3 years as supplied.  
 References: Spencer, J.P.E., et al. 1995. FEBS Lett. 374, 233.  
 Park, E.-M., et al. 1992. Proc. Natl. Acad. Sci. USA 89, 3375.  
 Lin, T.-S., et al. 1985. J. Med. Chem. 28, 1194.

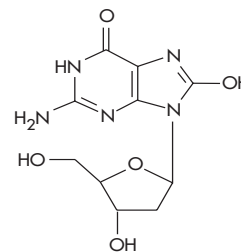

## 2,3-dimethoxy-1-naphthoquinone

26510


DMNQ

Catalog Number 26510

For Research Use Only, Not Intended For Use in Diagnostic Procedures

The use of naphthoquinoid compounds as free radical initiators is often compromised by their propensity to undergo nucleophilic addition, e.g. with sulfur-containing nucleophiles like glutathione. DMNQ is a novel, non-alkylating, non-thiol, adduct-forming compound which can be used to eliminate the previous mechanistic ambiguity involving redox-cycling quinoids as the source of reactive oxidant species/oxidative stress in biological studies.

Source: Synthetic Purity: >99% (TLC)  
 Form: Crystalline (solid)  
 Packaging Size: 200 mg  
 Storage and Stability: One year at 2° - 8° C  
 Molecular Weight: 218  
 Molecular Formula: C<sub>12</sub>H<sub>10</sub>O<sub>4</sub>

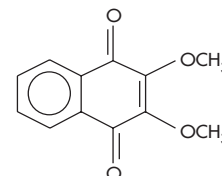

## 1-Methyl-4-vinyl-pyridinium Trifluoromethane Sulfonate

26513


M4VP

Catalog Number 26513

For Research Use Only, Not Intended For Use in Diagnostic Procedures

Specifically designed to trap mercaptans in aqueous samples. Applications include: elimination of interferences due to mercaptans in the measurement of biological parameters, inhibition of chemical or enzymatic reactions involving sulfhydryl groups and reversible protection of these groups in aqueous medium.

Source: Synthetic  
 Purity: >97% (GSH-DTNB titration)  
 Form: Crystalline white powder  
 Specificity: At neutral pH, only mercaptans are trapped. Other nucleophiles, such as amines or alcohols, do not react or react slowly.  
 Packaging Size: 500 mg  
 Storage and Stability: One year at 2° - 8° C  
 Molecular Weight: 269  
 Molecular Formula: C<sub>8</sub>H<sub>10</sub>N<sup>+</sup>(CF<sub>3</sub>SO<sub>3</sub><sup>-</sup>)

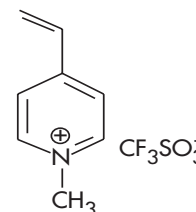

U.S. Patent 5543298



PRODUCT - CATALOG NUMBER .....	PAGE
1-Methyl-4-vinyl-pyridium Trifluoromethane Sulfonate - 26513 .....	95
2,3-dimethoxy-1-naphthoquinone - 26510 .....	95
3-Nitro-L-Tyrosine - 26527 .....	94
4-Hydroxy-2-nonenal - 26526 .....	93
4-Hydroxy-hex-2-enal-diethylacetal (100mg) - 26518 .....	92
4-Hydroxy-hex-2-enal-diethylacetal (1g) - 26519 .....	92
4-Hydroxy-hex-2-enal-diethylacetal (50mg) - 26594 .....	92
4-Hydroxy-non-enal-diethylacetal (100mg) - 26515 .....	92
4-Hydroxy-non-enal-diethylacetal (1g) - 26516 .....	92
4-Hydroxy-non-enal-diethylacetal (50mg) - 26595 .....	92
8-Hydroxy-2'-deoxyguanosine - 26571 .....	94
8-Hydroxyguanosine - 26525 .....	93
8-Isoprostane Assay - 21019 .....	40
8-OHdG Assay - 21026 .....	46
$\alpha_1$ -AP-410 Assay - 21047 .....	54
Aconitase-340 Assay - 21041 .....	48
Anti-4-Hydroxy-2-nonenal (100mg) - 24325 .....	82
Anti-4-Hydroxy-2-nonenal (20mg) - 24327 .....	82
Anti-8-Hydroxy-2'-deoxyguanosine (100mg) - 24326 .....	82
Anti-8-Hydroxy-2'-deoxyguanosine (20mg) - 24328 .....	82
Antibodies .....	78-83
Anti-Brain Nitric Oxide Synthase - 25312 .....	83
Anti-Catalase - 24316 .....	78
Anti-Cu/Zn - Superoxide Dismutase - 24322 .....	81
Anti-Cytochrome P450 2E1 - 24317 .....	79
Anti-Cytochrome P450 3A4 - 24318 .....	79
Anti-Endothelial Nitric Oxide Synthase - 25310 .....	82
Anti-Glutathione S-Transferase A1-1 - 24319 .....	80
Anti-Glutathione S-Transferase M1-1 - 24320 .....	80
Anti-Glutathione S-Transferase P1-1 - 24321 .....	80
Anti-Inducible Nitric Oxide Synthase - 25311 .....	83
Anti-Lactoferrin - 24311 .....	78
Anti-Mn Superoxide Dismutase - 24323 .....	81
Anti-Myeloperoxidase - 24310 .....	78
Anti-Nitrotyrosine - 24312 .....	78
Antioxidants .....	74
AOP-490 Assay - 21052 .....	36
Bioxytech® Antioxidant Biomarkers .....	16-37
Bioxytech® Inflammatory Biomarkers .....	60-63
Bioxytech® Nitrosative Biomarkers .....	64-73
Bioxytech® Oxidative Biomarkers .....	38-59
Bovine Cu/Zn Superoxide Dismutase (bSOD) (1g) - 25416 .....	84
Bovine Cu/Zn Superoxide Dismutase (bSOD) (250mg) - 25414 .....	84
Bovine Cu/Zn Superoxide Dismutase (bSOD) (25mg) - 25447 .....	84
Bovine Cu/Zn Superoxide Dismutase (bSOD) (500mg) - 25415 .....	84
Brain Nitric Oxide Synthase, Recombinant Rat - 27510 .....	88
Catalase Control ( <i>Aspergillus niger</i> ) - 27622 .....	77
Catalase, <i>Aspergillus niger</i> - 25430 .....	84
Catalase, Bovine Liver - 25431 .....	85
Catalase, Human Erythrocyte - 25432 .....	85
Catalase-520 Assay - 21042 .....	30
Cellular Glutathione Peroxidase Control - 27617 .....	76
cGPx-340 Assay - 21017 .....	22

PRODUCT - CATALOG NUMBER ..... PAGE

Constitutive Nitric Oxide Synthase Recombinant Human, High Purity - 27512 .....	89
Constitutive Nitric Oxide Synthase Recombinant Rat Neuronal - 27511 .....	89
Controls .....	76-77
Cu/Zn SOD Control - 27619 .....	76
Enzymes .....	84-90
G6PD   6PGD-340 Assay - 21045 .....	32
Glutathione Reductase Control - 27618 .....	76
Glutathione Reductase, Yeast - 25433 .....	86
Glutathione-S-Transferase A1-1 Recombinant Human - 25434 .....	86
Glutathione-S-Transferase M1-1 Recombinant Human - 25435 .....	87
Glutathione-S-Transferase P1-1 Recombinant Human - 25436 .....	87
GR-340 Assay - 21018 .....	24
GSH/GSSG-412 Assay - 21040 .....	28
GSH-400 Assay - 21011 .....	18
GSH-420 Assay - 21023 .....	26
GST-340 Assay - 21046 .....	34
H <sub>2</sub> O <sub>2</sub> -560 Assay - 21024 .....	42
HAE-586 Assay - 21043 .....	50
Inducible Nitric Oxide Synthase, Mouse Macrophage - 27513 .....	90
LactofEIA Assay - 21015 .....	62
L-Ergothioneine (100 mg) - 26511 .....	74
L-Ergothioneine (250 mg) - 26524 .....	74
LPO-560 Assay - 21025 .....	44
LPO-586 Assay - 21012 .....	38
Markers .....	92-94
MDA-586 Assay - 21044 .....	52
MPO-EIA Assay - 21013 .....	60
Myeloperoxidase, Human Polymorphonuclear Leukocytes - 25437 .....	88
Nitric Oxide Assay (enzymatic) - 22110 .....	66
Nitric Oxide Assay (non-enzymatic) - 22111 .....	68
Nitric Oxide Synthase Assay (colorimetric) - 22113 .....	72
Nitric Oxide Synthase Assay (radioactive arginine) - 22112 .....	70
Nitrotyrosine-EIA Assay - 21055 .....	64
Plasma GPx-EIA Assay - 21014 .....	20
SOD-525 Assay - 21010 .....	16
Specialty Chemicals .....	95
Urinary 8-EPI-PGF <sub>2α</sub> Assay - 21048 .....	56
Urinary F <sub>2</sub> -Isoprostane Metabolite Assay - 21049 .....	58

PRODUCT CATALOG NUMBER INDEX ON PAGE 98...

CATALOG NUMBER - PRODUCT .....	PAGE
21010 - SOD-525 Assay .....	16
21011 - GSH-400 Assay .....	18
21012 - LPO-586 Assay .....	38
21013 - MPO-EIA Assay .....	60
21014 - Plasma GPx-EIA Assay .....	20
21015 - Lactoferrin EIA Assay .....	62
21017 - cGPx-340 Assay .....	22
21018 - GR-340 Assay .....	24
21019 - 8-Isoprostane Assay .....	40
21023 - GSH-420 Assay .....	26
21024 - H <sub>2</sub> O <sub>2</sub> -560 Assay .....	42
21025 - LPO-560 Assay .....	44
21026 - 8-OHdG Assay .....	46
21040 - GSH/GSSG-412 Assay .....	28
21041 - Aconitase-340 Assay .....	48
21042 - Catalase-520 Assay .....	30
21043 - HAE-586 Assay .....	50
21044 - MDA-586 Assay .....	52
21045 - G6PD   6PGD-340 Assay .....	32
21046 - GST-340 Assay .....	34
21047 - α <sub>1</sub> -AP-410 Assay .....	54
21048 - Urinary 8-EPI-PGF <sub>2α</sub> Assay .....	56
21049 - Urinary F <sub>2</sub> -Isoprostane Metabolite Assay .....	58
21052 - AOP-490 Assay .....	36
21055 - Nitrotyrosine-EIA Assay .....	64
22110 - Nitric Oxide Assay (enzymatic) .....	66
22111 - Nitric Oxide Assay (non-enzymatic) .....	68
22112 - Nitric Oxide Synthase Assay (radioactive arginine) .....	70
22113 - Nitric Oxide Synthase Assay (colorimetric) .....	72
24310 - Anti-Myeloperoxidase .....	78
24311 - Anti-Lactoferrin .....	78
24312 - Anti-Nitrotyrosine .....	78
24316 - Anti-Catalase .....	78
24317 - Anti-Cytochrome P450 2E1 .....	79
24318 - Anti-Cytochrome P450 3A4 .....	79
24319 - Anti-Glutathione S-Transferase A1-1 .....	80
24320 - Anti-Glutathione S-Transferase M1-1 .....	80
24321 - Anti-Glutathione S-Transferase P1-1 .....	80
24322 - Anti-Cu/Zn - Superoxide Dismutase .....	81
24323 - Anti-Mn Superoxide Dismutase .....	81
24325 - Anti-4-Hydroxy-2-nonenal (100mg) .....	82
24326 - Anti-8-Hydroxy-2'-deoxyguanosine (100mg) .....	82
24327 - Anti-4-Hydroxy-2-nonenal (20mg) .....	82
24328 - Anti-8-Hydroxy-2'-deoxyguanosine (20mg) .....	82
25310 - Anti-Endothelial Nitric Oxide Synthase .....	82
25311 - Anti-Inducible Nitric Oxide Synthase .....	83
25312 - Anti-Brain Nitric Oxide Synthase .....	83
25414 - Bovine Cu/Zn Superoxide Dismutase (bSOD) (250mg) .....	84
25415 - Bovine Cu/Zn Superoxide Dismutase (bSOD) (500mg) .....	84
25416 - Bovine Cu/Zn Superoxide Dismutase (bSOD) (1g) .....	84
25430 - Catalase, <i>Aspergillus niger</i> .....	84
25431 - Catalase, Bovine Liver .....	85
25432 - Catalase, Human Erythrocyte .....	85

CATALOG NUMBER - PRODUCT .....	PAGE
25433 - Glutathione Reductase, Yeast .....	86
25434 - Glutathione-S-Transferase A1-1 Recombinant Human .....	86
25435 - Glutathione-S-Transferase M1-1 Recombinant Human .....	87
25436 - Glutathione-S-Transferase P1-1 Recombinant Human .....	87
25437 - Myeloperoxidase, Human Polymorphonuclear Leukocytes .....	88
25447 - Bovine Cu/Zn Superoxide Dismutase (bSOD) (25mg) .....	84
26510 - 2,3-dimethoxy-1-naphthoquinone .....	95
26511 - L-Ergothioneine (100 mg) .....	74
26513 - 1-Methyl-4-vinyl-pyridium Trifluoromethane Sulfonate .....	95
26515 - 4-Hydroxy-non-enal-diethylacetal (100mg) .....	92
26516 - 4-Hydroxy-non-enal-diethylacetal (1g) .....	92
26518 - 4-Hydroxy-hex-2-enal-diethylacetal (100mg) .....	92
26519 - 4-Hydroxy-hex-2-enal-diethylacetal (1g) .....	92
26524 - L-Ergothioneine (250 mg) .....	74
26525 - 8-Hydroxyguanasine .....	93
26526 - 4-Hydroxy-2-nonenal .....	93
26527 - 3-Nitro-L-Tyrosine .....	94
26571 - 8-Hydroxy-2'-deoxyguanasine .....	94
26594 - 4-Hydroxy-hex-2-enal-diethylacetal (50mg) .....	92
26595 - 4-Hydroxy-non-enal-diethylacetal (50mg) .....	92
27510 - Brain Nitric Oxide Synthase, Recombinant Rat .....	88
27511 - Constitutive Nitric Oxide Synthase Recombinant Rat Neuronal .....	89
27512 - Constitutive Nitric Oxide Synthase Recombinant Human, High Purity .....	89
27513 - Inducible Nitric Oxide Synthase, Mouse Macrophage .....	90
27617 - Cellular Glutathione Peroxidase Control .....	76
27618 - Glutathione Reductase Control .....	76
27619 - Cu/Zn SOD Control .....	76
27622 - Catalase Control (Aspergillus niger) .....	77

PRODUCT ALPHABETICAL INDEX ON PAGE 96...



